

Instructions for Use

M-100 Series (~~M1030~~, ~~M0830~~, M0820 and M0616)

- ① Screw the micro lens into the main unit.
- ② Hold the focusing ring and extend the lens barrel. The maximum magnification ratio is obtained when the lens barrel is fully extended. Use the fine focus ring to focus on a specimen (M1030 and M0830). The fine focus ring is not provided with the M0820 and M0616. Use only the lens barrel adjustment to focus.

M-200 Series (M820-S, S830-S, M820-AS and M820-SM)

- ① Screw the micro lens into the main unit.
- ② Adjust the focus of the built-in scale by the scale adjusting ring.
- ③ Focus on a specimen by using the fine focus ring of the micro lens.

