



User Manual

Liquid Withdrawal Device for Liquid Nitrogen 25, 35, and 50 LD/LDB Dewars

Item Number: 03773-11

Introduction

The liquid nitrogen withdrawal device (LWD) provides convenient pressure transfers of liquid nitrogen from 25, 35, and 50 LD/LDB Dewars. The device clamps directly to the special neck flange that is found under the plastic trim collar around the neck of these containers. Flow rates up to 8 liters per minute are possible at operating pressures up to 6.5 psig (0.45 bar/45 kPa). Liquid can be dispensed directly from the LWD via the spout provided, or through a standard cryogenic hose connected to the liquid outlet.

Safety

Before using this equipment, read the instructions carefully. Also, read and understand the user manual supplied with your Dewar. The user manual gives detailed health and safety information. Additional copies of these publications can be obtained from your supplier. Do not modify LWD or Dewar or use in a manner other than that described in accompanying instructions and documentation.

Contact with liquid nitrogen or cold gas may cause serious freezing (frostbite) injury. Do not touch liquid or cold metal surfaces with your bare skin. Always wear a face shield and gloves when transferring liquid nitrogen.

Nitrogen gas can cause suffocation without warning. In a confined area, nitrogen gas from this equipment may displace air required for breathing. Store and use this equipment only in well-ventilated areas.

Use only containers designed for low-temperature liquids. Cryogenic containers should be filled SLOWLY to minimize stresses which can damage the container.

Vent all pressure from a Dewar before attempting to remove the liquid withdrawal device. If the band clamp securing the LWD to a Dewar is loosened while there is pressure in the Dewar, cold gas or liquid could be expelled causing personal injury.

Liquid withdrawal device is for use with liquid nitrogen only. LWD is for use on specific IC Biomedical Dewars —models LD25, LD35 and LD50.

Pressure in the Dewar with LWD installed should never exceed 6.5 psig (0.45 bar/45 kPa). Fill Dewar from supply source at 22 psig (1.52 bar/152 kPa) or less. Do not pressurize using compressed air.

Installation

1. Remove the necktube cap/core from the Dewar
2. Remove 3 flat-head mounting screws in the base of the plastic trim collar (Figure 1) around the neck of the Dewar, remove the collar and discard or save for later. It is not used while the LWD is installed.

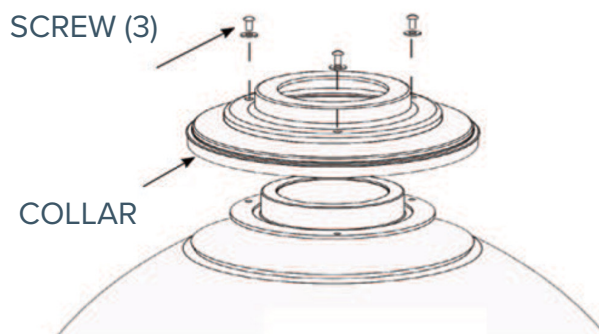


FIGURE 1

Installation (continued)

3. The liquid withdrawal tube (the longer of the two plastic tubes) has 3 scores (Figure 2)
 - The liquid withdrawal tube, without breaking off any of the sections, is for use with the older 50 LDB vessel.
 - Breaking off the tube at point 1, the first score from the bottom of the tube is for use with the newer 50 LD vessel.
 - Breaking off the tube at point 2, the second score from the bottom of the tube is for use with the 35 LD, 35 LDB and the older 25 LDB vessels.
 - Breaking off the tube at point 3, the third score from the bottom of the tube is for use with the newer 25 LD vessels.
 - Discard any removed pieces.

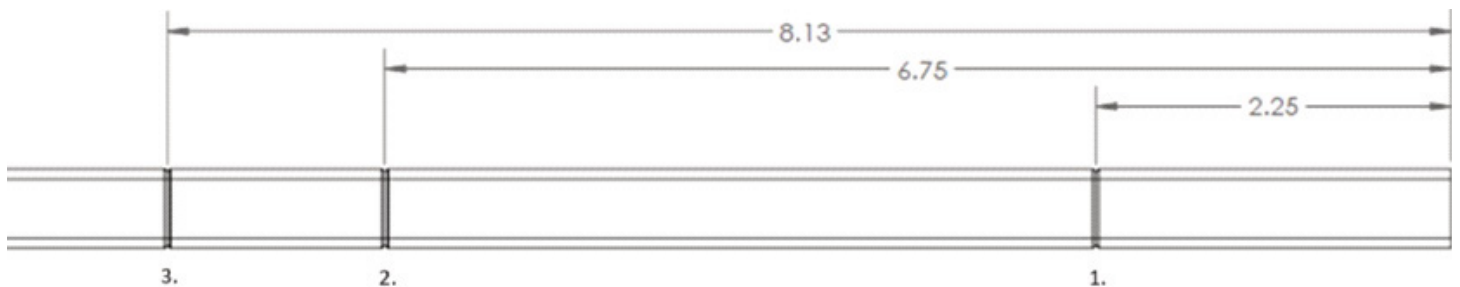


FIGURE 2

4. Insert the longer plastic tube into the underside of the flange assembly, beneath the LIQUID valve. Insert the shorter plastic tube into the threaded fitting on the underside of the flange assembly, beneath the VENT valve. Tighten the compression nuts with a wrench to secure.
5. Position the safety cable assembly around the top of the Dewar, underneath the flange. Loosen the V-band clamp until the clamp is large enough to remove from the LWD. Drop the band clamp around the flange on the Dewar (see Figure 3 on page 4).
6. Clean the flange O-ring and apply a thin layer of silicone grease before placing it on the grooved flange.
7. NOTE: Be sure flange surfaces are warm, clean and dry. The O-ring must be cleaned and lightly lubricated before installing the LWD.
8. Carefully lower the LWD unit into the Dewar. The LWD should rest firmly centered on the neck flange of the Dewar; the liquid withdrawal tube should not touch the bottom of the Dewar.
9. Place the band clamp around the flange of both the LWD unit and the Dewar. Be sure that the clamp is properly engaged on both flanges. Tighten the clamp—hand tighten only.
10. Attach the snap of the safety cable to the ring on the LWD (see Figure 3).
11. If the spout assembly is to be used, install it on the outlet fitting above the LIQUID valve. The phase separator will ensure a controlled flow of liquid.

Operation

Filling—It is not necessary to remove the LWD unit from the Dewar for filling. You may fill the Dewar by pressure—transfer of liquid nitrogen through the LIQUID valve of the LWD.

WARNING: Over pressurization of the Dewar could result in serious bodily injury or property damage. Never exceed the 6.5 psig (0.45 bar/45 kPa) relief valve pressure setting of the LWD.

To fill, remove the spout assembly and connect a 22 psig (1.52 bar/152 kPa) or lower pressure liquid supply cylinder to the withdrawal device using a liquid nitrogen transfer hose with female CGA 295 connections on both ends.

WARNING: The transfer hose must be used with a pressure relieve valve to prevent excessive pressure build-up in the hose when the valves are closed at both ends of the hose. If liquid nitrogen is trapped between the valves of the liquid cylinder and the LWD, it has no escape path as it warms and expands. The resulting pressure can cause the hose to rupture and may cause personal injury. The current model LWD is equipped with a pressure relief valve. Previous versions of the LWD must be used with a transfer hose equipped with a pressure relief valve.

Once the transfer hose is connected, open the LWD VENT valve first and then open the LWD LIQUID valve. Then slowly open the liquid valve on the supply cylinder until liquid flow is established. Adjust the VENT valve to maintain a back pressure of 6.5 psig (0.45 bar/45 kPa) or less in the Dewar. When liquid sputters from the VENT valve, quickly close the liquid valve on the supply cylinder, next the LWD LIQUID valve and then the LWD VENT valve.

WARNING: Closely monitor Dewar pressure throughout the fill process. Never let the pressure exceed 6.5 psig (0.45 bar/45 kPa). If the pressure approaches 6.5 psig (0.45 bar/45 kPa), immediately fully open the LWD VENT valve and/or close the LWD LIQUID valve as necessary to control the pressure in the Dewar.

Loosen the hose connection slowly and allow all pressure and any remaining liquid to dissipate before completely disconnecting the hose.

The primary relief valve, set at 6.5 psig (0.45 bar/45 kPa), maintains the normal operating pressure of the LWD. A secondary relief valve, set at 6.5 psig (0.45 bar/45 kPa), prevents hazardous pressure rise in the event of primary relief valve malfunction or if its capacity is exceeded. Never replace either of these parts with higher relief valves; increasing the pressure in the Dewar beyond 6.5 psig (0.45 bar/45 kPa) may burst the Dewar and cause personal injury/property damage.

LWD on the filled Dewar, leave the VENT valve open until the flanges are securely clamped.

NOTE: Be sure flanges are warm, clean and dry. The O-ring must be cleaned and lightly lubricated before replacing the LWD.

Liquid Withdrawal—when the liquid valve is opened, pressure in the Dewar forces liquid to go up through the withdrawal tube and out through the valve. The liquid nitrogen can be dispensed through the spout or through the transfer hose connected to the liquid outlet.

Pressure Building—If the Dewar is filled properly, liquid nitrogen evaporation will maintain a pressure sufficient to sustain liquid withdrawal.


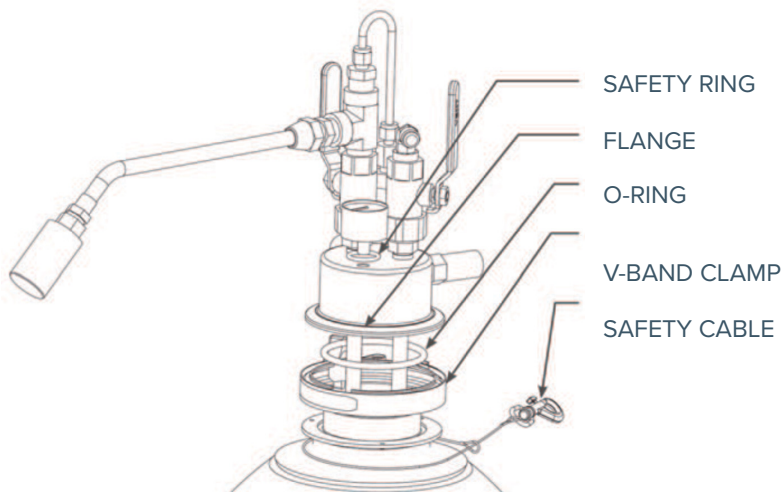


FIGURE 3

Maintenance

It is recommended to replace the relief valves at least every three years to ensure correct operation of the withdrawal device. Other parts should be replaced as required. See item number and information below.

Parts

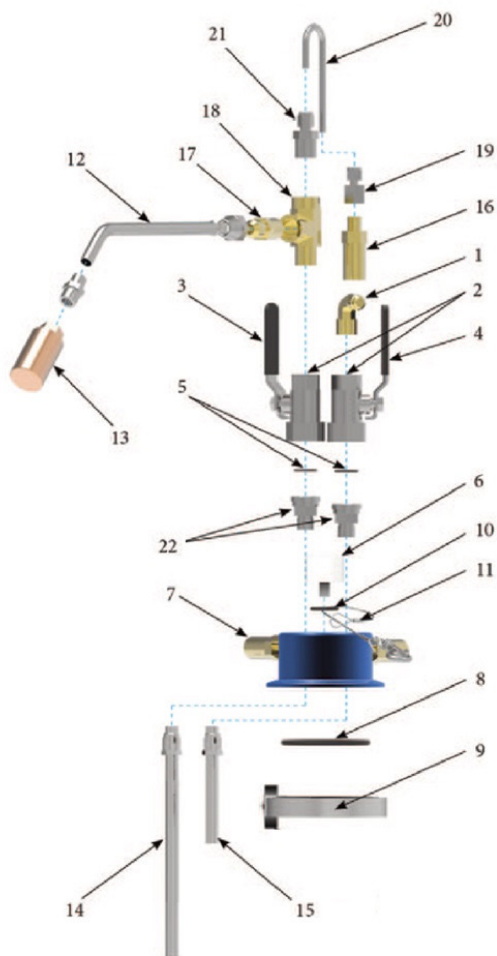

Item	Description	Item Number
1	Elbow, Vent	6814-9225
2	Valve (Liquid/Vent)	6986-9004
3	Valve Handle Cover, Liquid	373608
4	Valve Handle Cover, Vent	373607
5	O-Ring for Valve	8830-6142
6	Pressure Gauge 0—15 psig	7702-6095
7	Relief Valve, 6.5 psig (0.45 bar/45 kPa)	6913-9095
8	O-Ring for Flange	8830-6135
9	V-Band Clamp, with Warning Label	D050-8C12
10	Ring for Safety Cable	8830-6315
11	Safety Cable Assembly	D050-8C20
12	Spout Assembly	D050-8C14
13	Phase Separator	1193-8C80
14	Withdrawal Tube Assembly	D050-8C37
15	Vent Tube Assembly	D050-8C39
16	Relief Valve, 22 psig (2.4 bar/24 kPa)	6913-9069
17	Connector, CGA, 295 x ½ in. FTP	7355-4698
18	Street Tee, ½ in. NPT	6916-9211
19	Connector, ¼ in. ODT x ¼ in. FTP	45701930
20	Tube, Relief	D050-9C10
21	Connector, ¼ in. ODT x ½ in. MPT	45701929
22	Valve Adapter	D050-8C29

Ordering Information

Order all replacement parts and accessories from your distributor. Please include the part and model number of your Dewar, quantity, and description of each part requested. For more information, contact Cole-Parmer.

Cole-Parmer®

625 East Bunker Court
Vernon Hills, IL 60061
United States

United States

T: +1.800.323.4340 or
+1.847.549.7600
E: sales@antylia.com
W: coleparmer.com

Canada

T: +1.514.355.6100
E: info@antylia.ca
W: coleparmer.ca

China

T: 86.21.5109.9909
E: sales@antylia.com
W: coleparmer.cn

France

T: +33 (0) 1486 37800
E: fr.sales@antylia.com
W: coleparmer.fr

Germany

T: +49 (0) 9377 92030
E: de.sales@antylia.com
W: coleparmer.de

India

T: +9122 61394444
E: info@coleparmer.in
W: coleparmer.in

Italy

T: +39 (0) 284349215
E: it.sales@antylia.com
W: coleparmer.it

UK

T: +44 (0) 1480 272279
E: uk.sales@antylia.com
W: coleparmer.co.uk

Other

T: +1.847.549.7600