

08904-00

Instruction for Dry Keeper Desiccator Cabinet 42053-0000

The 42053-0000 Dry Keeper Desiccator Cabinet is very simple to use. Roll back the two roller latches to open the door. Remove the jar of Silica Gel Beads and remove the plastic bag with the beads from inside the jar. Carefully cut open one end of the plastic bag and empty the Silica Gel Beads into the jar. Place the jar with the cap removed back into the cabinet and close the door and replace the roller latches back to their locked position.

When the Silica Gel Beads become pink in color, they can be removed, emptied into a small heat resistant dish and regenerated in a vented oven at 150°C (300°F) for three hours or until they are blue in color.

Post-1125-5-11-2008 7671	
[REDACTED]	
Phone #	Phone #
Fax #	Fax #