# **polystat®**

 Cole-Parmer®

## **Cooling/Heating Circulating Baths**

Manual P/N U01206 Rev. 03/13/2017

### **Cooling/Heating Circulating Baths**

**1C6            2C15            3C6**

**1C6F 2        2C15++        3C6F**

**1C15                            3C15**

**1C15++                        3C15++**

### **User's Manual**

### **Heating Circulating Baths**

**1W7            2W7            3W7**

**1W11            2W11            3W11**

**1W24            2W24            3W24**

**1W5P**

**1W14P**

**1W21P**

# Table of Contents

<b>Preface .....</b>	<b>i</b>
Warranty .....	i
Unpacking .....	i
WEEE .....	ii
Contact Information .....	ii
<b>Section 1     Safety .....</b>	<b>1-1</b>
Safety Warnings .....	1-1
Element Assessment .....	1-3
Personal Protective Equipment .....	1-4
Training.....	1-4
<b>Section 2     General Information.....</b>	<b>2-1</b>
Description .....	2-1
Intended Use .....	2-1
Sample Nameplate.....	2-1
Equipment Ratings .....	2-2
Wetted Parts.....	2-2
Specifications .....	2-3
Approved Fluids.....	2-10
Additional Fluid Precautions .....	2-12
<b>Section 3     Installation .....</b>	<b>3-1</b>
Cooling Heating Circulating Bath .....	3-1
Ventilation.....	3-1
Electrical Requirements .....	3-1
External Circulation .....	3-3
Filling Requirements .....	3-4
Draining.....	3-4
<b>Section 4     Operation .....</b>	<b>4-1</b>
Circulating Bath.....	4-1
Setup .....	4-2
Start Up .....	4-2
Status Display .....	4-3
Standby Mode.....	4-3
Changing the Setpoint.....	4-4
Menu Displays.....	4-5
Menu .....	4-5
Menu Tree .....	4-6
Settings .....	4-7
High Temperature Cutout.....	4-15

High Temperature Cutout .....	4-16
Stopping the Circulating Bath.....	4-17
Power Down.....	4-17
Shut Down.....	4-17
Restarting .....	4-17
Decommissioning/Disposal .....	4-18
Storage .....	4-18
<b>Section 5      Preventive Maintenance.....</b>	<b>5-1</b>
Cleaning .....	5-1
Condenser Fins .....	5-1
Hoses .....	5-1
Testing the Safety Features.....	5-2
<b>Section 6      Troubleshooting.....</b>	<b>6-1</b>
Error Displays .....	6-1
Check List .....	6-3

## **Appendix**

# Preface

## Warranty

Cole-Parmer warrants to the direct purchaser that the Cole-Parmer Polystat Circulating Baths, distributed by Cole-Parmer, will be free from defects in material or workmanship for a period of two years from the date of delivery.

Cole-Parmer will repair or replace the product or provide credit, as its sole option, upon prompt notification and compliance with its instructions.

The Distributor warrants to Customer that upon prompt notification and compliance with Distributor's instructions, that the Distributor will repair or replace, at Distributor's sole option, any Product which is defective in material or workmanship.

Distributor expressly disclaims all other warranties, whether expressed, implied or statutory, including the warranties of merchantability, and fitness for a particular purpose. Distributor's sole responsibility and the Customer's exclusive remedy for any claim arising out of the purchase of any Product is repair or replacement, as described above. In no event shall Distributor's liability exceed the purchase price paid therefor; nor shall Distributor be liable for any claims, losses or damage of any third party or for lost profits or any special, indirect, incidental, consequential, or exemplary damages, howsoever arising, even if Distributor has been advised of the possibility of such damages.

## Unpacking

Retain all cartons and packing material until the circulating bath is operated and found to be in good condition. If the circulating bath shows external or internal damage contact the transportation company and file a damage claim. Under ICC regulations, this is your responsibility.

The circulating bath does not have handles. Take into account its weight when unpacking and transporting, see Section 2. We recommend a minimum of two people, or using a harness, when lifting. ▲

Cooling/heating circulating baths should be left in an upright position at room temperature for 24 hours before starting. This will ensure the lubrication oil has drained back into the compressor. ▲

## **WEEE**

This product is required to comply with the European Union's Waste Electrical & Electronic Equipment (WEEE) Directive 2002/96/EC. It is marked with the following symbol:

## **Contact Information**

Cole-Parmer Instrument Company  
625 East Bunker Court  
Vernon Hills, IL 60061  
1-800-323-4340

## Section 1 Safety

### Safety Warnings

Make sure you read and understand all instructions and safety precautions listed in this manual before installing or operating your circulating bath. If you have any questions concerning the operation or the information in this manual, please contact us.

**DANGER** indicates an imminently hazardous situation which, if not avoided, *will* result in death or serious injury.

**WARNING** indicates a potentially hazardous situation which, if not avoided, *could* result in death or serious injury.

**CAUTION** indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury. It is also be used to alert against unsafe practices.

The lightning flash with arrow symbol, within an equilateral triangle, is intended to alert the user to the presence of non-insulated "dangerous voltage" within the circulating bath's enclosure. The voltage magnitude is significant enough to constitute a risk of electrical shock.

This label indicates the presence of hot surfaces.

This label indicates read the manual.

**Observe all warning labels. ▲**

**Never remove warning labels. ▲**

**Cooling circulating baths should be left in an upright position for 24 hours before starting. This will ensure the lubrication oil has drained back into the compressor. ▲**

**The circulating bath's construction provides protection against the risk of electrical shock by grounding appropriate metal parts. The protection will not function unless the power cord is connected to a properly grounded outlet. It is the user's responsibility to assure a proper ground connection is provided. ▲**

**The circuit protector located on the rear of the circulating bath is not intended to act as a disconnecting means. ▲**

**Operate the circulating bath using only the supplied line cord. The circulating bath's power cord is used as the disconnecting device, it must be easily accessible at all times. ▲**

Never place the circulating bath in a location or atmosphere where excessive heat, moisture, or corrosive materials are present. ▲

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Ensure the tubing you select will meet your maximum temperature and pressure requirements. ▲

Ensure all communication and electrical connections are made prior to starting. ▲

Many refrigerants which may be undetectable by human senses are heavier than air and will replace the oxygen in an enclosed area causing loss of consciousness. Refer to the circulating bath's nameplate and the manufacturer's most current SDS for additional information. ▲

Never operate the circulating bath without fluid in the reservoir. ▲

Use only the approved fluids listed in this manual. ▲

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Other than water, before using any other approved fluid, or when performing maintenance where contact with the fluid is likely, refer to the manufacturer's SDS and EC Safety Data sheet for handling precautions. ▲

The user is responsible for decontamination if hazardous materials are spilled. Consult the manufacturer regarding decontamination and or cleaning agents compatibility. ▲

Ensure, that no toxic gases can be generated by the fluid. Flammable gases can build up over the fluid during usage. ▲

When using glycols and water, check the fluid concentration and pH on a regular basis. Changes in concentration and pH can impact system performance. ▲

Ensure the fluid is at a safe temperature (below 40°C) before handling or draining. ▲

Never operate damaged or leaking equipment, or with any damaged cords. ▲

Never operate the circulating bath or add fluid to the reservoir with panels removed. ▲

Do not clean the reservoir with solvents, only use a soft cloth and water. ▲

Drain the circulating bath before it is transported and/or stored in, near or below freezing temperatures.

Always turn the circulating bath off and disconnect the supply voltage from its power source before moving or before performing any service or maintenance procedures. ▲

Transport the circulating bath with care. Sudden jolts or drops can damage its components. ▲

Refer service and repairs to a qualified technician. ▲

Performance of installation, operation, or maintenance procedures other than those described in this manual may result in a hazardous situation and voids the manufacturer's warranty. ▲

# Element Assessment

## **Chemical Used —**

For cooling/heating circulating baths only:

R134A Refrigerant - CFC Free

C6/C6F 3.5 ounces

C15 6.2 ounces

C15++ (R404A) 9.3 ounces

Location: Refrigeration system

Lubricating Oil:

C6, C6F and C15 POE – 243ml

C15++ Ester Oil – 450 ml

Location: Compressor

## **Hazard Communication — None**

**Electrical** — All electrical energy sources for the circulating bath are provided by the user.

**Hazardous Energy Isolation** — Locate energy isolation devices in a location that is readily accessible.

**Electrical** Remove power from the circulating bath by turning the circuit protector off and disconnecting the line cord, see Section 3 **Electrical Requirements**.

**Fluid** Before using any fluid refer to the manufacturer's SDS and EC safety data sheets for handling precautions.

**Refrigeration** The cooling/heating circulating bath's refrigeration system is sealed and requires no lockout/tagout.

Remove any isolation devices before starting the circulating bath.

**Chemicals** — Prior to decontamination procedures should be developed and employed to minimize or mitigate opportunities for exposure to chemicals present in or around the circulating bath.

**Radioactive Material** — None

**Magnetic Fields** — None

## **Sampling and Analysis for External Contaminants from Other Sources**

**or from Spill Events** — It is the responsibility of the user to identify the presence of all external chemicals before the decontamination. This should be accomplished by collecting samples and analyzing them for any suspected chemicals used or stored in areas adjacent to the bath.

**Oxygen Deficiency** — Many refrigerants which may be undetectable by human senses are heavier than air and will replace the oxygen in an enclosed area causing loss of consciousness. Refer to the circulating bath's nameplate and the manufacturer's most current SDS for additional information. Test the atmosphere in a confined space area or area where gases are being used to purge the circulating bath. If necessary, use an air supplied respirator.

## **Personal Protective Equipment**

The are no special personal protective equipment requirements needed to perform normal operation. We do recommend wearing gloves and eye protection.

## **Training**

The user must review and understand all the sections in this manual before operating the circulating bath.

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## **Fluorinated Greenhouse Gases**

The following information is included to comply with REGULATION (EU) No 517/2014 OF THE EUROPEAN PARLIAMENT AND OF THE COUNCIL of 16 April 2014 on fluorinated greenhouse gases:

**This product contains fluorinated greenhouse gases in a hermetically sealed system.**

**If a leak in the sealed system is detected, the operator shall repair without undue delay.**

Refer to the F-Gas Declaration of Conformity for additional information.

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## F-GAS DECLARATION OF CONFORMITY

### Declaration of conformity with Article 14 of Regulation (EU) No 517/2014 of the European Parliament and of the Council

We, Thermo Fisher Scientific (Asheville) LLC, F-Gas Portal Registration Number 23643, declare under our sole responsibility that when placing on the market pre-charged equipment, which we import to or manufacture in the Union, the hydrofluorocarbons contained in that equipment are accounted for within the quota system referred to in Chapter IV of Regulation (EU) No 517/2014 as:

We hold authorisation(s) issued in accordance with Article 18(2) of Regulation (EU) No 517/2014 and registered in the registry referred to in Article 17 of that Regulation, at the time of release for free circulation to use the quota of a producer or importer of hydrofluorocarbons subject to Article 15 of Regulation (EU) No 517/2014 that cover(s) the quantity of hydrofluorocarbons contained in the equipment.

The hydrofluorocarbons contained in the equipment have been placed on the market in the Union, subsequently exported and charged into the equipment outside the Union, and the undertaking that placed the hydrofluorocarbons on the market made a declaration stating that the quantity of hydrofluorocarbons has been or will be reported as placed on the market in the Union and that it has not been and will not be reported as direct supply for export in the meaning of Article 15(2)(c) of Regulation (EU) No 517/2014 pursuant to Article 19 of Regulation (EU) No 517/2014 and Section 5C of the Annex to Commission Implementing Regulation (EU) No 1191/2014.

January 1, 2017

Mark Pearson, Director Global Regulatory Affairs

---

Date

[name and position of legal representative]

---

[signature of legal representative]

Manufacturer:

Thermo Fisher Scientific (Asheville) LLC  
275 Aiken Road  
Asheville, NC 28804  
U.S.A.

EU Only Representative:

Thermo Electron LED GmbH  
Robert-Bosch-Strasse 1  
D-63505 Langenselbold  
Germany  
VAT ID Number: DE 812 403 137

## Section 2 General Information

### Description

The Cole Parmer Polystat® Cooling/Heating Circulating Baths can pump to an external system. All have a digital display and easy-to-use touch pad, five programmable setpoint temperatures, acoustic and optical alarms, and offer adjustable high temperature protection.

### Intended Use

Intended use is to provide heating, cooling or constant temperature to applications placed into the reservoir and/or pumped to an external application in accordance with all the procedures and requirements stated in this manual.

Intended for use in a laboratory environment, on a bench top only.

Typical lab applications include:

Analytical Instrumentation	Lasers
General Laboratory Cooling	Condensers
Rotary Evaporators	Bio-reactors
Fermenters	Histology

### Sample Nameplate

Use the nameplate on the rear to identify your immersion circulator/circulating bath, its electrical requirements and, if applicable, the refrigeration data. Cole Parmer catalog part numbers are listed in the specification tables in this section.

## Equipment Ratings

Compliance

UL File Number: E473332

Ambient Temperature Range	10°C to 40°C (50°F to 104°F)
Maximum Relative Humidity (Non Condensing)	10% to 80% (non-condensing)
Operating Altitude	Sea Level to 2000 meters (6560 feet)
Oversupply Category	II
Pollution Degree	2
Degree of Protection	IP 20
Sound Power Level (Varies with pump speed)	55 to 60 dBA
Refrigerant	C6 and C6F R134A 99 grams (3.5 ounces) C15 R134A 176 grams (6.2 ounces) C15++ R404A 264 grams (9.3 ounces)
Storage Temperature Range	-25°C to 60°C (-13°F to 140°F)

## Wetted Materials

### Circulator

Viton  
EPDM  
Ryton  
Ultem  
Vectra  
Stainless Steel

### Stainless Steel Baths/Circulators

Stainless Steel 316  
Stainless Steel 304  
EPDM (drain fitting)  
Ryton  
Zotek-N (cover seal)

### Polyphenylene oxide (PPO) Baths/ Circulators

Polyphenylenoxid

## Cooling/Heating Circulating Baths Specifications

Stainless Steel Refrigerated/Heated Bath Circulators with Standard Circulator				
Catalog Part Number	12122-02	12122-16	12122-32	12122-46
Nameplate Identification	1C6	1C6F	1C15	1C15++
<b>Temperature Range</b> °C °F	-20 to 100 -4 to 212	-20 to 100 -4 to 212	-28 to 100 -18 to 212	-35 to 100 -31 to 212
<b>Heater Capacity</b> KWatts	1.0	1.0	1.0	1.0
<b>Stability</b> °C	0.05	0.05	0.05	0.05
<b>Bath Volume</b> liters gallons	5.4 - 6.5 1.4 - 1.7	5.4 - 6.5 1.4 - 1.7	6.8 - 8.6 1.8 - 2.3	6.8 - 8.6 1.8 - 2.3
<b>Cooling Capacity</b> watts @20°C	250	250	500	800
<b>Refrigerant</b>	R134a	R134a	R134a	R404a
<b>Dimensions (H x W x L)*</b> mm inches	622.3 x 203.2 x 416.6 24.5 x 8.0 x 16.4	426.7 x 467.4 x 416.6 16.8 x 18.4 x 16.4	650.2 x 259.1 x 490.2 25.6 x 10.2 x 19.3	688.3 x 370.8 x 528.3 27.1 x 14.6 x 20.8
<b>Electrical</b>				
Nominal VAC	115V	115V	115V	115V
Operating VAC Range	103V - 127V	103V - 127V	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz	60 Hz	60 Hz
Input Current Rating Amps	11.5	11.5	11.7	14.4
Total Wattage	1328	1328	1345	1662
Line Cord Plug (NEMA)	N5-15	N5-15	N5-15	N5-20
<b>Net Weight</b> kg lb	26.3 58.0	29.0 64.0	35.8 79.0	54.9 121.0

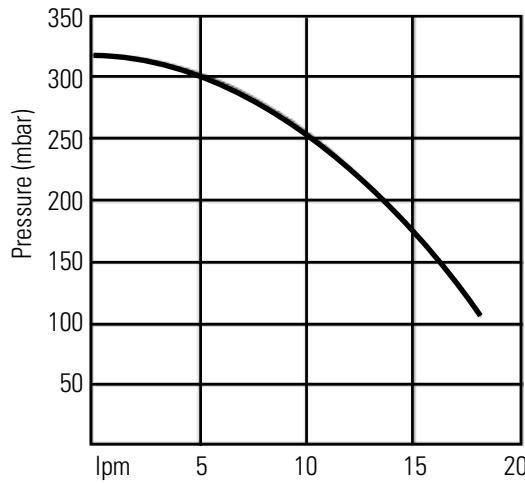
\*Overall dimensions. Add ~15 mm to L for drain fitting

- Stability is measured as follows:


Fluid (specific heat of 0.55 Btu/lb-F) at -10°C, work area cover is on. Baths run in factory ambient at nominal line voltage. Pump speed set to high with short insulated loop on pump lines.

Stability is defined as ½ the total span of measured data over approximately 30 minutes.

- Cole Parmer reserves the right to change specifications without notice.

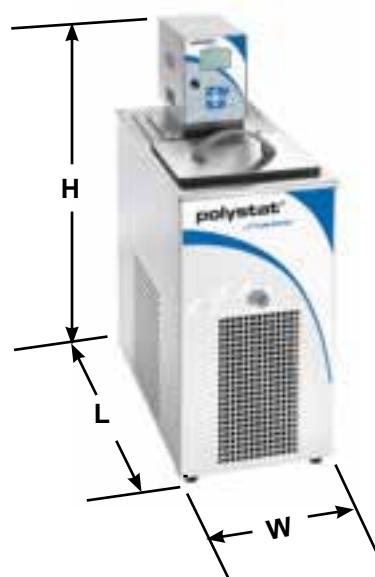- Pump testing is done with water at 20°C bath at nominal line voltage on high pump speed. Approximately one meter loop on pump with flow transducer and (2) pressure transducers with a ball valve to adjust the flow rate.

<b>Stainless Steel Cooling/Heating Circulating Baths with Standard Plus</b>		
<b>Catalog Part Number Nameplate Identification</b>	<b>12122-36 2C15</b>	<b>12122-52 2C15++</b>
<b>Temperature Range °C °F</b>	-28 to 150 -18 to 302	-35 to 150 -31 to 302
<b>Heater Capacity KWatts</b>	1.0	1.0
<b>Stability °C</b>	0.05	0.05
<b>Bath Volume liters gallons</b>	6.8 - 8.6 1.8 - 2.3	6.8 - 8.6 1.8 - 2.3
<b>Cooling Capacity watts @20°C</b>	500	800
<b>Refrigerant</b>	R134a	R404a
<b>Dimensions (H x W x L)* mm inches</b>	650.2 x 259.1 x 490.2 25.6 x 10.2 x 19.3	688.3 x 370.8 x 528.3 27.1 x 14.6 x 20.8
<b>Electrical</b>		
Nominal VAC	115V	115V
Operating VAC Range	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz
Input Current Rating Amps	11.7	14.4
Total Wattage	1345	1662
Line Cord Plug (NEMA)	N5-15	N5-20
<b>Net Weight kg lb</b>	35.4 78.0	54.9 121.0

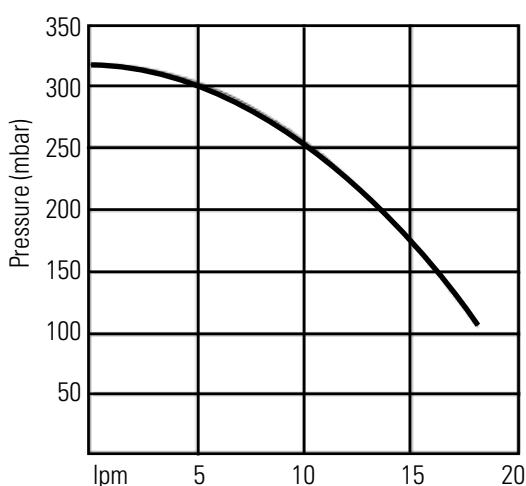
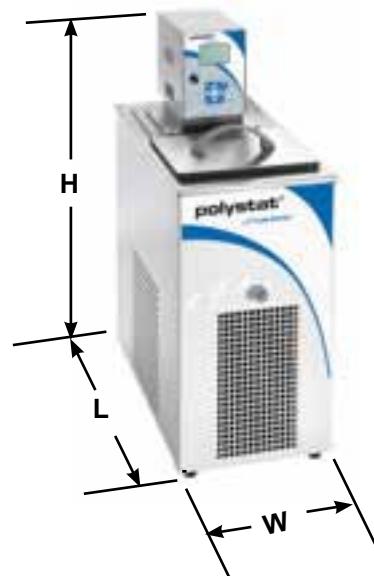
\*Overall dimensions. Add ~15 mm to L for drain fitting

- Stability is measured as follows:

Fluid (specific heat of 0.55 Btu/lb-F) at -10°C, work area cover is on. Baths run in factory ambient at nominal line voltage. Pump speed set to high with short insulated loop on pump lines.

Stability is defined as  $\frac{1}{2}$  the total span of measured data over approximately 30 minutes.

- Cole Parmer reserves the right to change specifications without notice.

- Pump testing is done with water at 20°C bath at nominal line voltage on high pump speed. Approximately one meter loop on pump with flow transducer and (2) pressure transducers with a ball valve to adjust the flow rate.

<b>Stainless Steel Cooling/Heating Circulating Baths with Advanced Circulator</b>				
<b>Catalog Part Number Nameplate Identification</b>	<b>12122-12 3C6</b>	<b>12122-26 3C6F</b>	<b>12122-42 3C15</b>	<b>12122-56 3C15++</b>
<b>Temperature Range °C °F</b>	-20 to 100 -4 to 212	-20 to 100 -4 to 212	-28 to 200 -18 to 392	-35 to 200 -31 to 392
<b>Heater Capacity KWatts</b>	1.2	1.2	1.2	1.2
<b>Stability °C</b>	0.025	0.025	0.025	0.025
<b>Bath Volume liters gallons</b>	5.4 - 6.5 1.4 - 1.7	5.4 - 6.5 1.4 - 1.7	6.8 - 8.6 1.8 - 2.3	6.8 - 8.6 1.8 - 2.3
<b>Cooling Capacity watts @20°C</b>	250	250	500	800
<b>Refrigerant</b>	R134a	R134a	R134a	R404a
<b>Dimensions<sup>1</sup> (H x W x L) mm inches</b>	640.0 x 203.2 x 416.6 25.2 x 8.0 x 16.4	444.5 x 467.4 x 416.6 17.5 x 18.4 x 16.4	655.3 x 259.1 x 490.2 25.8 x 10.2 x 19.3	706.1 x 370.8 x 528.3 27.8 x 14.6 x 20.8
<b>Electrical</b>				
Nominal VAC	115V	115V	115V	115V
Operating VAC Range	103V - 127V	103V - 127V	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz	60 Hz	60 Hz
Input Current Rating Amps	11.5	11.5	11.7	14.4
Total Wattage	1328	1328	1345	1662
Line Cord Plug (NEMA)	N5-15	N5-15	N5-15	N5-20
<b>Net Weight kg lb</b>	27.2 60.0	29.9 66.0	36.3 80.0	55.8 123.0

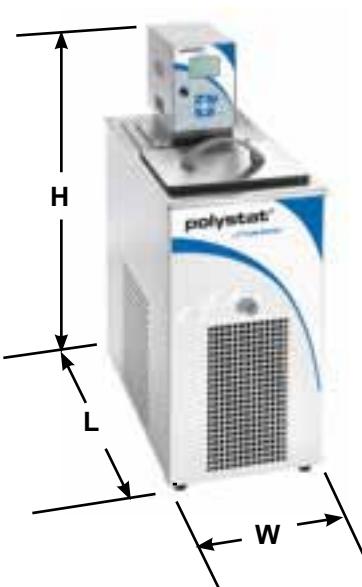
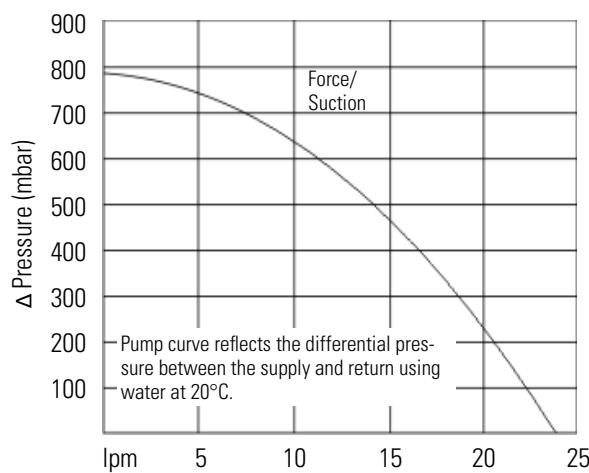
\*Overall dimensions. Add ~15 mm to L for drain fitting

- Stability is measured as follows:

Fluid (specific heat of 0.55 Btu/lb-F) at -10°C, work area cover is on. Baths run in factory ambient at nominal line voltage. Pump speed set to high with short insulated loop on pump lines.

Stability is defined as ½ the total span of measured data over approximately 30 minutes.

- Cole Parmer reserves the right to change specifications without notice.

- Pump testing is done with water at 20°C bath at nominal line voltage on high pump speed.  
Approximately one meter loop on pump with flow transducer and (2) pressure transducers with a ball valve to adjust the flow rate.

## Heating Circulating Baths Specifications

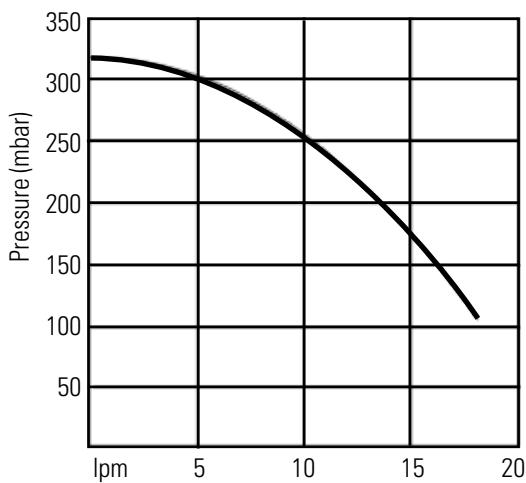
Stainless Steel Heating Circulating Baths with Standard Circulator			
Catalog Part Number Nameplate Identification	12134-60 1W7	12134-64 1W11	12134-68 1W24
<b>Temperature Range</b> °C* °F*	Ambient +10 to 100 Ambient +18 to 212	Ambient +10 to 100 Ambient +18 to 212	Ambient +10 to 100 Ambient +18 to 212
<b>Heater Capacity</b> KWatts	1.0	1.0	1.0
<b>Stability</b> °C	0.05	0.05	0.05
<b>Bath Volume</b> liters gallons	5.1 - 6.5 1.3 - 1.7	7.5 - 9.5 2.0 - 2.5	15.4 - 19.6 4.1 - 5.2
<b>Dimensions** (H x W x L)</b> mm inches	396.2 x 215.9 x 363.2 15.6 x 8.5 x 14.3	396.2 x 302.2 x 363.2 15.6 x 11.9 x 14.3	396.2 x 363.2 x 561.3 15.6 x 14.3 x 22.1
<b>Electrical</b>			
Nominal VAC	115V	115V	115V
Operating VAC Range	103V - 127V	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz	60 Hz
Input Current Rating Amps	9.6	9.6	9.6
Total Wattage	1007	1007	1007
Line Cord Plug (NEMA)	N5-15	N5-15	N5-15
<b>Net Weight</b> kg lb	10.0 22.0	11.3 25.0	15.9 35.0

\*Without work area cover installed.

Lower temperature ranges require supplemental cooling.

\*\*Overall dimensions. Add ~15 mm to L for drain fitting.

- Lower process fluid temperature ranges available with supplemental cooling.
- Stability is measured as follows:  
Fluid is water at 70°C, work area cover is on. Baths run in factory ambient at nominal line voltage. Pump speed set to high with no external pumping.  
Stability is defined as ½ the total span of measured data over approximately 30 minutes.
- Cole Parmer reserves the right to change specifications without notice.

- Pump testing is done with water at 20°C bath at nominal line voltage on high pump speed. Approximately one meter loop on pump with flow transducer and (2) pressure transducers with a ball valve to adjust the flow rate.


<b>Stainless Steel Heating Circulating Baths with Standard Plus Circulator</b>			
<b>Catalog Part Number Nameplate Identification</b>	<b>12121-02 2W7</b>	<b>12121-22 2W11</b>	<b>12121-42 2W24</b>
<b>Temperature Range</b> °C*	Ambient +10 to 150	Ambient +10 to 150	Ambient +10 to 150
°F*	Ambient +18 to 302	Ambient +18 to 302	Ambient +18 to 302
<b>Heater Capacity</b> KWatts	1.0	1.0	1.0
<b>Stability</b> °C	0.05	0.05	0.05
<b>Bath Volume</b> liters gallons	5.1 - 6.5 1.3 - 1.7	7.5 - 9.5 2.0 - 2.5	15.4 - 19.6 4.1 - 5.2
<b>Dimensions** (H x W x L)</b> mm inches	396.2 x 215.9 x 363.2 15.6 x 8.5 x 14.3	396.2 x 302.2 x 363.2 15.6 x 11.9 x 14.3	396.2 x 363.2 x 561.3 15.6 x 14.3 x 22.1
<b>Electrical</b>			
Nominal VAC	115V	115V	115V
Operating VAC Range	103V - 127V	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz	60 Hz
Input Current Rating Amps	9.6	9.6	9.6
Total Wattage	1007	1007	1007
Line Cord Plug (NEMA)	N5-15	N5-15	N5-15
<b>Net Weight</b> kg lb	10.0 22.0	11.3 25.0	15.9 35.0

\*Without work area cover installed.

Lower temperature ranges require supplemental cooling.

\*\*Overall dimensions. Add ~15 mm to L for drain fitting.

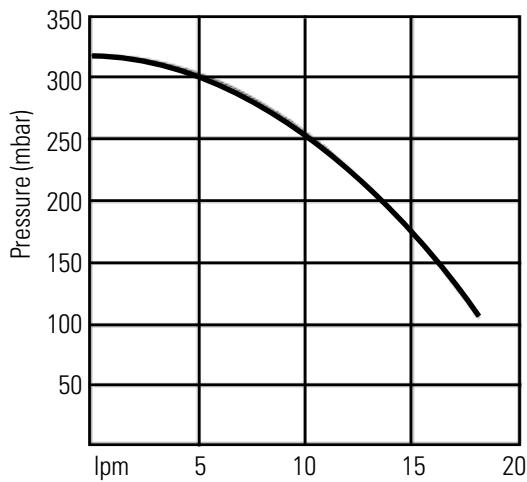- Lower process fluid temperature ranges available with supplemental cooling.
- Stability is measured as follows:  
Fluid is water at 70°C, work area cover is on. Baths run in factory ambient at nominal line voltage. Pump speed set to high with no external pumping.  
Stability is defined as ½ the total span of measured data over approximately 30 minutes.
- Cole Parmer reserves the right to change specifications without notice.

- Pump testing is done with water at 20°C bath at nominal line voltage on high pump speed. Approximately one meter loop on pump with flow transducer and (2) pressure transducers with a ball valve to adjust the flow rate.


<b>Stainless Steel Heating Circulating Baths with Advanced Circulator</b>			
<b>Catalog Part Number Nameplate Identification</b>	<b>12121-06 3W7</b>	<b>12121-26 3W11</b>	<b>12121-46 3W24</b>
<b>Temperature Range °C*</b> (Low pump speed) °F*	Ambient +10 to 200 Ambient +18 to 392	Ambient +10 to 200 Ambient +18 to 392	Ambient +10 to 200 Ambient +18 to 392
<b>Temperature Range °C*</b> (High pump speed) °F*	Ambient +25 to 200 Ambient +45 to 392	Ambient +20 to 200 Ambient +36 to 392	Ambient +10 to 200 Ambient +18 to 392
<b>Heater Capacity KWatts</b>	1.2	1.2	1.2
<b>Stability °C</b>	0.025	0.025	0.025
<b>Bath Volume liters gallons</b>	5.1 - 6.5 1.3 - 1.7	7.5 - 9.5 2.0 - 2.5	15.4 - 19.6 4.1 - 5.2
<b>Dimensions** (H x W x L) mm inches</b>	414.0 x 215.9 x 363.2 16.3 x 8.5 x 14.3	414.0 x 302.2 x 363.2 16.3 x 11.9 x 14.3	414.0 x 363.2 x 561.3 16.3 x 14.3 x 22.1
<b>Electrical</b>			
Nominal VAC	115V	115V	115V
Operating VAC Range	103V - 127V	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz	60 Hz
Input Current Rating Amps	11.3	11.3	11.3
Total Wattage	1304	1304	1304
Line Cord Plug (NEMA)	N5-15	N5-15	N5-15
<b>Net Weight kg lb</b>	10.9 24.0	12.2 27.0	16.8 37.0

\*Without work area cover installed.

Lower temperature ranges require supplemental cooling.

\*\*Overall dimensions. Add ~15 mm to L for drain fitting.

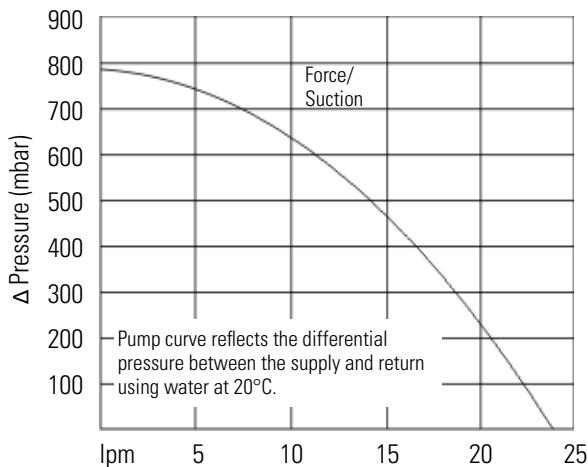- Lower process fluid temperature ranges available with supplemental cooling.
- Stability is measured as follows:  
Fluid is water at 70°C, work area cover is on. Baths run in factory ambient at nominal line voltage. Pump speed set to high with no external pumping.  
Stability is defined as ½ the total span of measured data over approximately 30 minutes.
- Cole Parmer reserves the right to change specifications without notice.

- Pump testing is done with water at 20°C bath at nominal line voltage on high pump speed.

Approximately one meter loop on pump with flow transducer and (2) pressure transducers with a ball valve to adjust the flow rate.

## Open Heating Circulating Baths Specifications

<b>Polyphenylene oxide (PPO) Heating Circulating Baths with Standard Circulator</b>			
<b>Catalog Part Number</b>	<b>12134-30 1W5P</b>	<b>12134-34 1W14P</b>	<b>12134-38 1W21P</b>
<b>Temperature Range</b> °C*	Ambient +10 to 100	Ambient +10 to 100	Ambient +10 to 100
°F*	Ambient +18 to 212	Ambient +18 to 212	Ambient +18 to 212
<b>Heater Capacity</b> KWatts	1.0	1.0	1.0
<b>Stability</b> °C	0.05	0.05	0.05
<b>Bath Volume</b> liters gallons	4.3 - 5.3 1.1 - 1.4	11.3 - 14.1 3.0 - 3.7	18.0 - 22.5 4.8 - 5.9
<b>Overall Dimensions</b> (H x W x L)			
mm	348.0 x 190.5 x 388.6	348.0 x 358.1 x 452.1	348.0 x 358.1 x 642.6
inches	13.7 x 7.5 x 15.3	13.7 x 14.1 x 17.8	13.7 x 14.1 x 25.3
<b>Electrical</b>			
Nominal VAC	115V	115V	115V
Operating VAC Range	103V - 127V	103V - 127V	103V - 127V
Operating Frequency	60 Hz	60 Hz	60 Hz
Input Current Rating Amps	9.6	9.6	9.6
Total Wattage	1007	1007	1007
Line Cord Plug (NEMA)	N5-15	N5-15	N5-15
<b>Net Weight</b> kg lb	5.3 11.6	6.8 14.9	7.5 16.6

\*Lower temperature ranges require supplemental cooling.

• Cole Parmer reserves the right to change specifications without notice.

## Approved Fluids

Cole Parmer takes no responsibility for damages caused by the selection of an unapproved fluids.

Only use the approved fluids listed on the chart below. ▲

<b>Approved Fluid</b>	Water	EG/Water	PG/Water	Sil 200 (50 cSt)	Sil 300	Synth 260
<b>Circulator</b>	All Models	All Models	All Models	Standard Plus/ Advanced only	Advanced only	Advanced only
<b>Flash Point*</b>	NA	NA	NA	>318°C	>300°C	>260°C
<b>Fire Point*</b>	NA	NA	NA	>350°C	>325°C	>275°C
<b>Operating Temp Range</b>						
200						
150						
100						
125						
100						
75	+5°C to +95°C	-30°C to +80°C	-30°C to +80°C	+20°C to +200°C	+80°C to +200°C	+45°C to +200°C
50						
25						
0						
-25						
-50						

\*The flash point and fire point refer to the fluid's flammability characteristics. The flash point is defined as the lowest temperature at which the vapor formed above a pool of the liquid ignites. The fire point is the lowest temperature at which, on further heating beyond the flash point, the sample will support combustion for 5 seconds. The bath liquid is limited to a temperature level 25°C below the fire point as defined by the EN 61010.

Handling and disposal of liquids other than water should be done in accordance with the fluid manufacturers specification and/or the SDS for the fluid used. ▲

When using water above 80°C closely monitor the fluid level, frequent top-offs will be required. It will also create steam. ▲

Water/glycol mixtures require top-offs with pure water, otherwise the percentage of glycol will increase resulting in high viscosity and poor performance. ▲

Polyphenylene oxide (PPO) Bath Circulators are used with water only. ▲

## Additional Fluid Information

### 5°C to 95°C — Distilled Water or Deionized Water (up to 3 MΩ-cm)

Normal tap water leads to calcareous deposits necessitating frequent decalcification, see table on next page.

Calcium tends to deposit itself on the heating element. The heating capacity is reduced and service life shortened.

### -30°C to 80°C — 50/50 Water with Laboratory Grade Ethylene or with Propylene Glycol

Below 5°C water has to be mixed with a glycol. The amount of glycol added should cover a temperature range 5°C lower than the operating temperature of the particular application. This will prevent the water/glycol from gelling (freezing) near the evaporating coil.

Excess glycol deteriorates the temperature accuracy due to its high viscosity.

- Other than water, do not use any approved fluid until you have read and understood the label and the material Safety Data Sheet (SDS).
- Ensure any fluid residue or any other material is thoroughly removed before filling the reservoir with a different fluid.
- Always wear protective clothing, especially a face shield and gloves.
- Avoid spattering on any of the circulating bath's components, always *slowly* add fluid. When adding, point the opening of a container away from yourself.
- Use fume hoods.
- Do not allow any ignition sources in the vicinity.

## Water Quality and Standards

<b>Process Fluid</b>	<b>Permissible (PPM)</b>	<b>Desirable (PPM)</b>
<b>Microbiologicals</b> (algae, bacteria, fungi)	0	0
<b>Inorganic Chemicals</b>		
Calcium	<25	<0.6
Chloride	<25	<10
Copper	<1.3	<1.0
		0.020 ppm if fluid in contact with aluminum
Iron	<0.3	<0.1
Lead	<0.015	0
Magnesium	<12	<0.1
Manganese	<0.05	<0.03
Nitrates\Nitrites	<10 as N	0
Potassium	<20	<0.3
Silicate	<25	<1.0
Sodium	<20	<0.3
Sulfate	<25	<1
Hardness	<17	<0.05
Total Dissolved Solids	<50	<10
<b>Other Parameters</b>		
pH	6.5-8.5	7-8
Resistivity	0.01*	0.05-0.1*

\* MΩ·cm (compensated to 25°C)

Unfavorably high total ionized solids (TIS) can accelerate the rate of galvanic corrosion. These contaminants can function as electrolytes which increase the potential for galvanic cell corrosion and lead to localized corrosion such as pitting. Eventually, the pitting will become so extensive that refrigerant will leak into the reservoir.

As an example, raw water in the United States averages 171 ppm (of NaCl). The recommended level for use in a water system is between 0.5 to 5.0 ppm (of NaCl).

Recommendation: Initially fill the reservoir with distilled or deionized water. Do not use untreated tap water as the total ionized solids level may be too high. This will reduce the electrolytic potential of the water and prevent or reduce the galvanic corrosion observed.

## Section 3 Installation

### Cooling/Heating Circulating Bath

**CAUTION**

**CAUTION**

### Ventilation

**CAUTION**

### Electrical Requirements

**DANGER**

The circulating bath is designed for continuous operation and for indoor use.

**Never place the circulating bath in a location where excessive heat, moisture, inadequate ventilation, or corrosive materials are present. ▲**

**Cooling/heating circulating baths should be left in an upright position for 24 hours at room temperature before starting. This will ensure the lubrication oil has drained back into the compressor. ▲**

Blocked ventilation will increase the circulating bath's temperature, reduce its cooling capacity and, on cooling/heating baths, eventually lead to premature compressor failure.

**When operating above 95°C ensure there is minimum clearance of 31 centimeters (12 inches) on all 4 sides. ▲**

The circulating bath construction provides protection against the risk of electrical shock by grounding appropriate metal parts. The protection will not function unless the power cord is connected to a properly grounded outlet. It is the user's responsibility to ensure a proper ground connection is provided. ▲

The circulating bath is intended for use on a dedicated outlet.

The circuit protector on the rear of the circulating bath is designed to protect the circulator's internal components.

**Note** If the circuit protector activates allow the temperature to cool before resetting. Restart the circulating bath. Contact us if it activates again. ▲

**The circulating bath's power cord is used as the disconnecting device, it must be easily accessible at all times. ▲**

Refer to the nameplate on the rear of the bath circulator for specific electrical requirements.

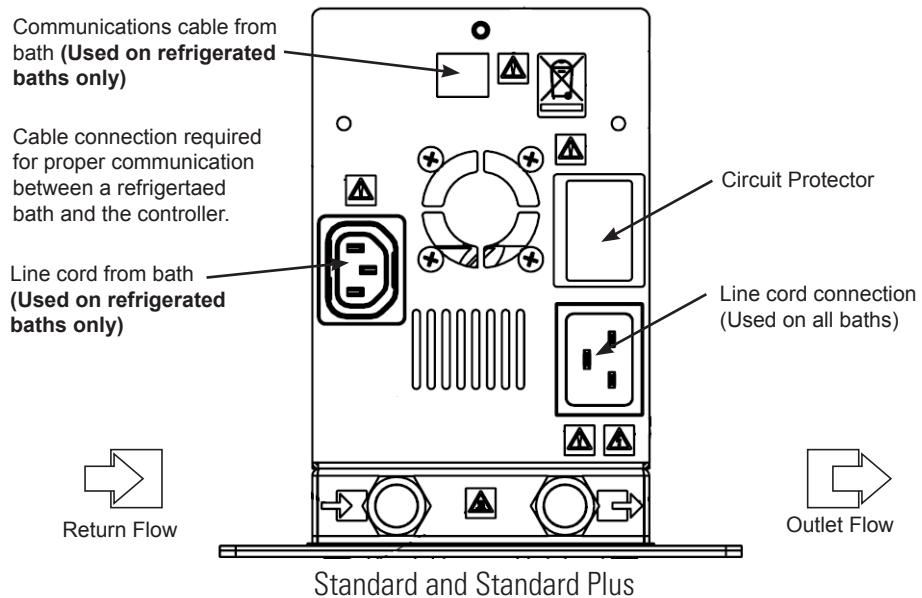
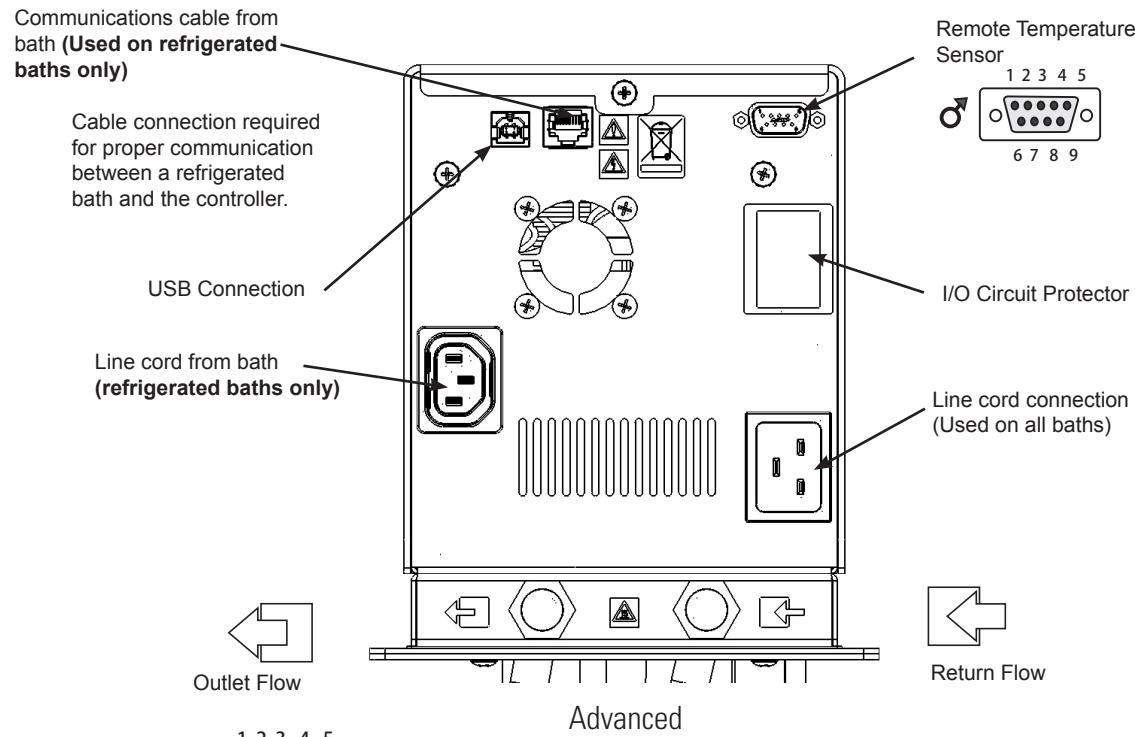
Ensure the electrical cords do not come in contact with any of the plumbing connections or tubing. ▲

Ensure all communication and electrical connections are made prior to starting the bath circulator and that the cords do not come in contact with any of the plumbing connections or tubing. ▲

- For refrigerated bath circulators, ensure the line cord from the bath is securely connected to the rear of the circulator.
- For refrigerated bath circulators, ensure the communications cable from the bath is connected to the RJ45 connectors (similar to Ethernet) on the rear of the circulator.
- For all circulators, connect the supplied line cord from the bath circulator to a properly grounded outlet.

## Remote Temperature Sensor (Advanced only)

**Never apply line voltage to the connection. ▲**

The optional remote temperature sensor is enabled using the bath circulator, see Section 4.

Pin		Pin	
1	White	7	Red
2, 3	NA	8	NA
4	White	9	Red (4th wire not connected to the control board)
5, 6	NA		

## External Circulation

The plumbing connections for external circulation are located on the rear of the bath circulator. is the return flow from the external application. is the outlet flow to the external application (supply side). The connections are 16 mm O.D. Remove the union nuts and plates to install the supplied 8 mm or 12 mm hose barbs and clamps.

**To prevent damage to the circulator's plumbing, use a 19 mm backing wrench when removing/installing the external connections. ▲**

## Filling Requirements

Ensure the reservoir drain port on the front of the circulating bath is *closed* and that all plumbing connections are secure. Also ensure any residue is thoroughly removed before filling.

**Before using any fluid refer to the manufacturer's SDS and EC safety data sheets for handling precautions. ▲**

To avoid spilling, place your containers into the circulating bath before filling.

With a low level WARNING the circulating bath continues to run, with a FAULT the circulating bath will shut the refrigeration, pump and heater will shut down, see Section 6. The low level warning is at approximately 4 cm (1 5/8") below the top of the bath, the low level fault is at approximately 5.5 cm (2 1/8") below it.

**Avoid overfilling, oil-based fluids expand when heated. ▲**

When pumping to an external system, keep extra fluid on hand to maintain the proper level in the circulating lines and the external system.

**Note Monitor the fluid level whenever heating the fluid. ▲**

**Before draining any fluid refer to the manufacturer's SDS and EC safety data sheets for handling precautions. ▲**

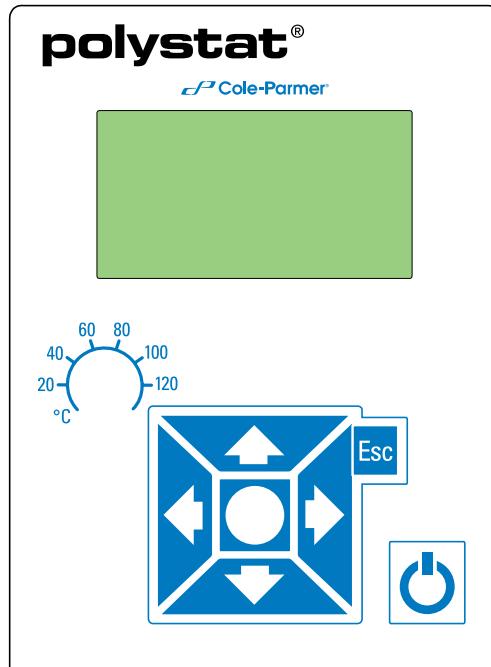
**Draining**  
**Ensure the fluid is at a safe handling temperature, below 40°C. Wear protective clothing and gloves. ▲**

- place a suitable receptacle underneath the drain. If desired, attach an 8 mm id tube on the drain.
- slowly turn the drain plug until flow is observed.


## Section 4 Operation

### Circulating Bath

The Cole Parmer Heating Cooling/Heating Circulating Baths have a digital display and easy-to-use touch pad, five programmable setpoint temperatures, acoustic and optical alarms and adjustable high temperature protection.

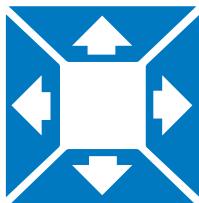Use this button to place the circulating bath in and out of standby mode.

Use these navigation arrows to move through the circulating bath's displays and to adjust values.

Use this button to start/stop the circulating bath. This button is also used to make and save changes on the display screen.

Use this button to cancel any changes and to return the circulating bath to its previous display. Canceling a change can only be made before the change is saved. In some cases, it is also used to save changes.

**Note** Holding this button for five seconds resets the display contrast to the default level and also brings up the language menu to change, if needed, the displayed language. See **Settings - Display Options** in this Section. ▲

Use the adjacent dial for adjusting the High Temperature Cutout. Details are explained in this Section.

## Setup

Leave cooling/heating circulating baths upright at room temperature (~25°C) for 24 hours before starting. This ensures the lubrication oil has drained back into the compressor. ▲

## Start Up

Before starting, double check all USB (optional), electrical and plumbing connections. ▲

Do not run the circulating bath until fluid is added. Have extra fluid on hand. If the circulating bath will not start refer to Section 5 Troubleshooting.

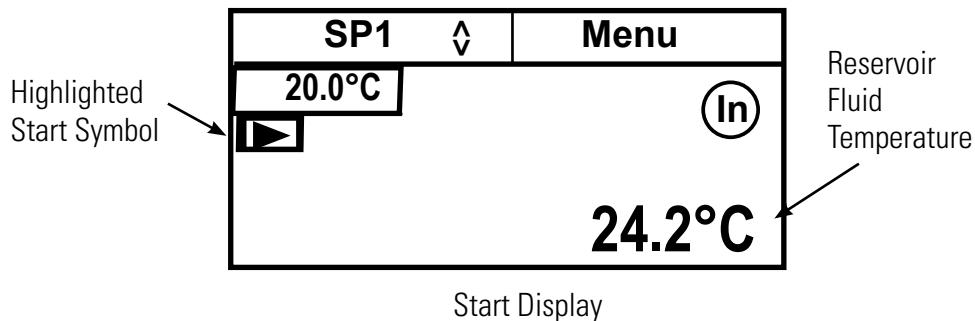
- Place the circuit protector to the **I** position.


- Press , the Start Display appears.

- Ensure the start symbol has a highlight box around it, if not use the arrow keys to navigate to the symbol.

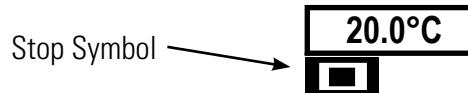- Press . The circulating bath starts and the start symbol turns into a stop symbol ( ).

**Note** It takes 30 seconds for the compressor to start. ▲


**Note** After start, check external plumbing connections for leaks. ▲

The **SP1** and **Menu** portions on the top of the display are used to view and/or change the circulating bath's settings. They are explained in detail later in this Section.

On Advanced circulating baths only, indicates the circulator is using its internal sensor for temperature control. indicates the circulator is using an external sensor for temperature control.

## Status Display

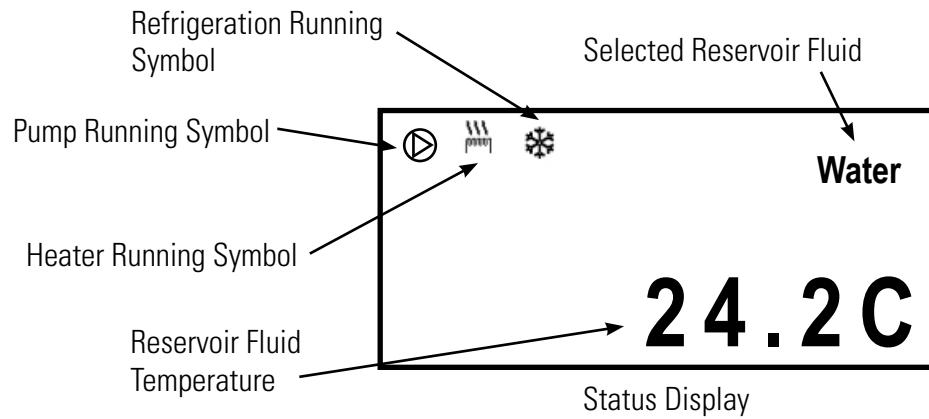If desired, press **Esc** to toggle between the Start/Status Displays.

## Standby Mode

Press , the display goes blank and the circulating bath enters the standby mode.

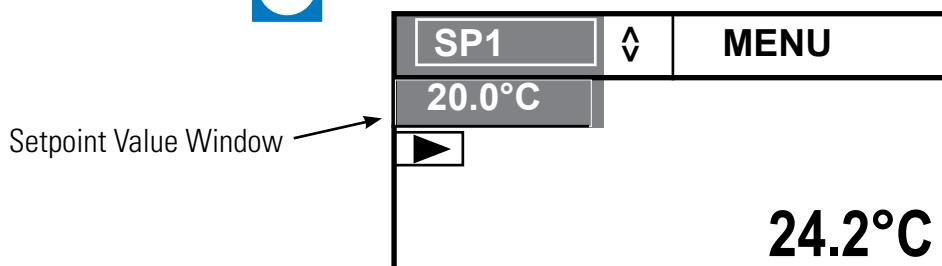
## Changing the Setpoint

**Note** You cannot adjust the setpoint closer than 0.1°C to either of the fluid's system limits, see Fluids Type in this Section, or beyond the circulating bath's temperature range. ▲

The setpoint can be changed with the circulating bath running or not.


The Setpoint is the desired fluid temperature. The circulating bath can store up to five setpoints, **SP1** through **SP5**. The procedure for changing the stored setpoint values is discussed later in this Section. Their factory default settings are 20°C, 37°C, 50°C, 70°C and 100°C.

Use the navigation arrows and move to the **SP1** window and then press  to highlight it as shown below.


Use the up and down navigation arrows to bring up the desired setpoint and then press .

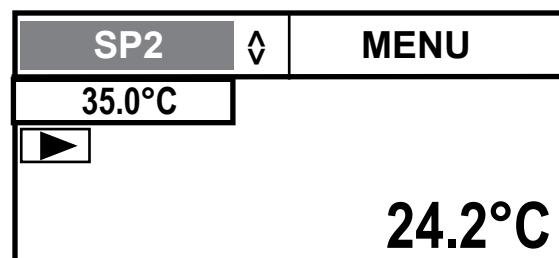The display on the Setpoint Value Window now indicates the corresponding setpoint's stored value.

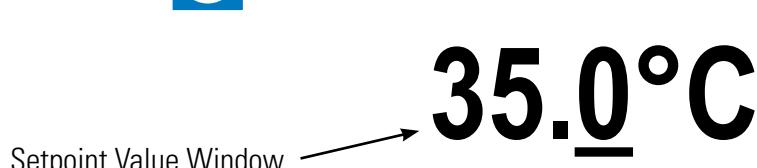If desired, you can change the displayed setpoint value by using the navigation arrows to highlight the Setpoint Value Window and then pressing .

Use the left and right arrows to move the cursor to the desired digit and then use the up and down arrows to change the value. Once all the desired changes are made, press  to save the change.

**Note** Using this procedure also changes that setpoint's stored value. ▲


## Menu Displays

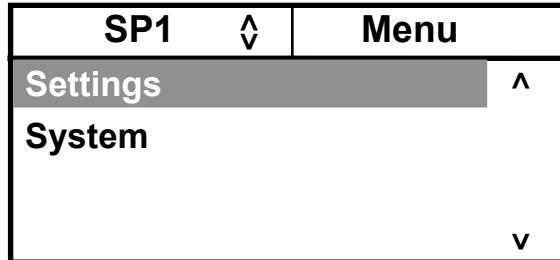The circulating bath uses menus to view/change its settings.


**Note** The circulating bath does not need to be running to view/change these settings. ▲

For all Menu displays, once  is pressed to change a display, you can press  to return to the previous screen.



1. Use the right arrow button to highlight **Menu** and the circulator brings up the Main Menu Display.

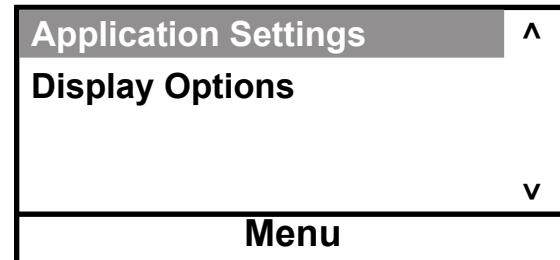2. Use the up and down arrow to highlight the desired setting and then press  to bring up additional submenus.

See page 4-7.

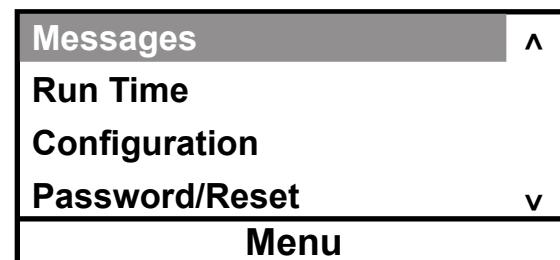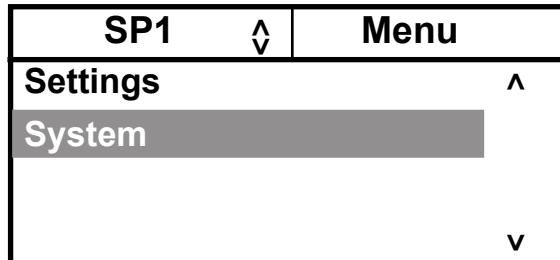
See page 4-12.


## Menu

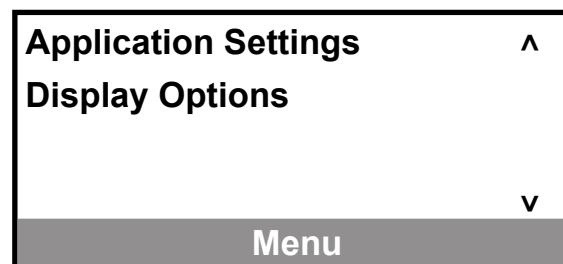The **Menu** line, at the bottom of all the submenu displays, is another way to return the circulating bath back to the Main Menu Display.

1. From any submenu display, use the down arrow button to highlight **Menu**.

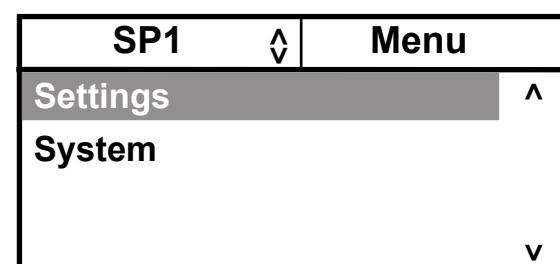
2. Press  to return to the Main Menu Display.

Pressing  from the Menu line returns the circulating bath to the previous screen.


## Menu Tree

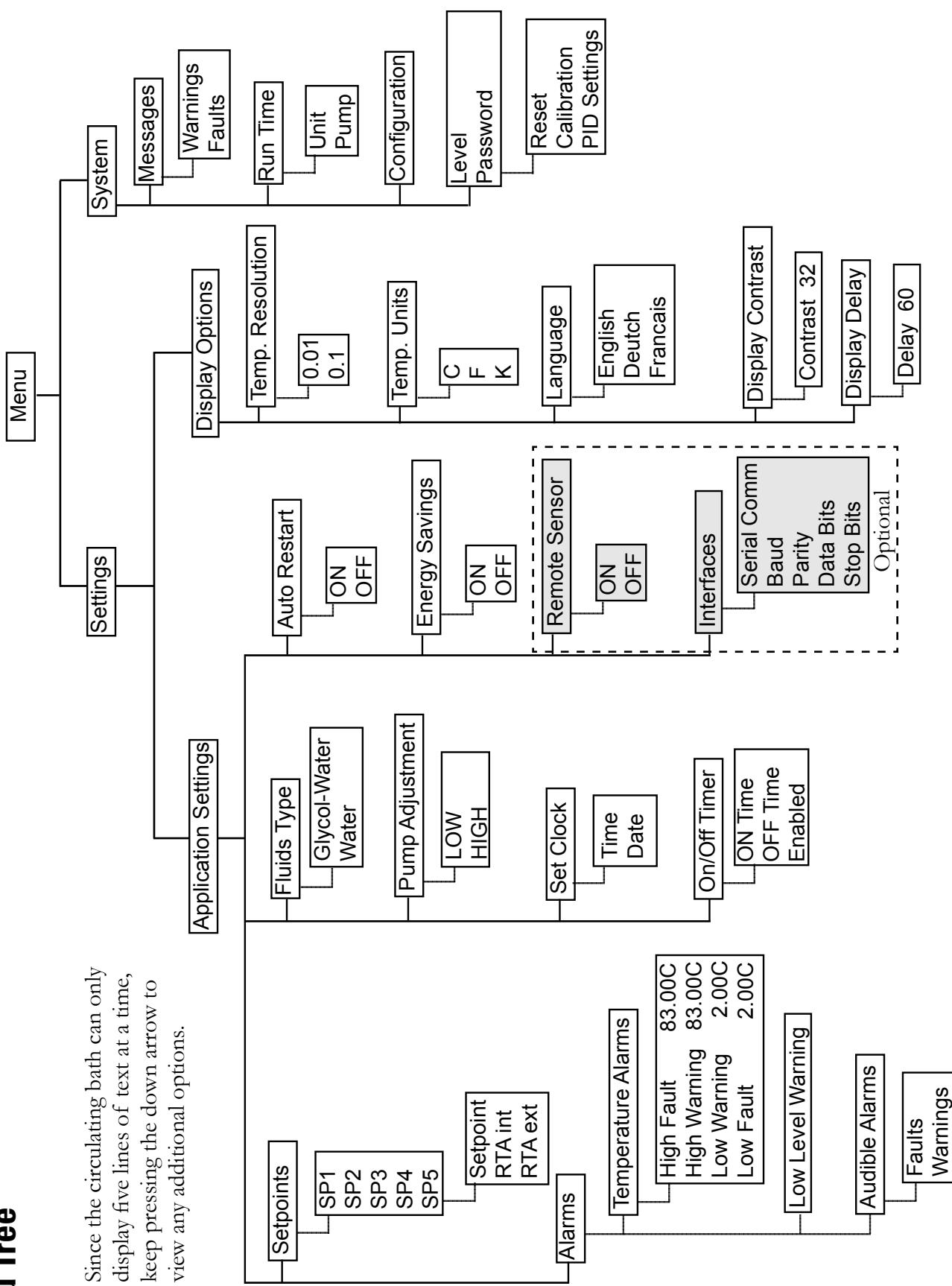Since the circulating bath can only display five lines of text at a time, keep pressing the down arrow to view any additional options.


**Settings - Application Settings** is used to view/adjust the circulating bath's five Setpoints and Real Temperature Adjustments (RTA), enable/disable the alarms, change the fluid type, set the pump speed, configure the interfaces (optional), set the clock, turn the timer on or off, and turn auto restart and energy savings on or off.

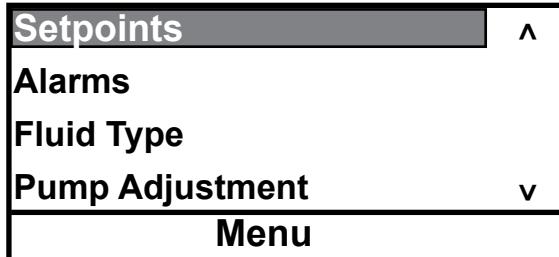
- With **Application Settings** highlighted press  to view:

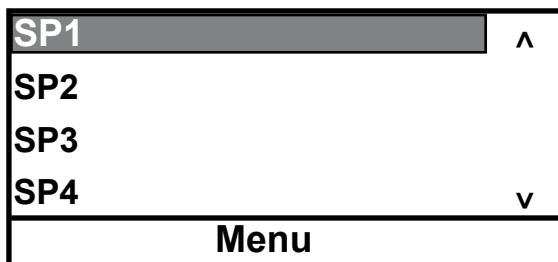- With **Setpoints** highlighted, press  to display the list. Use the up/down arrows to highlight the desired **SP**. **Note** Use the down arrow to display **SP5**. ▲


Scroll down for additional options:

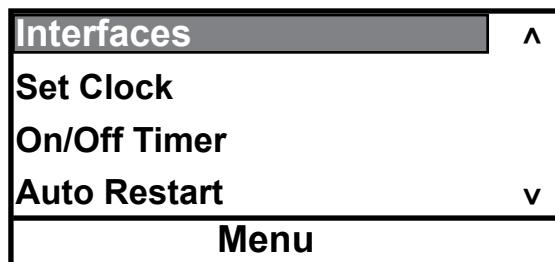- Press .

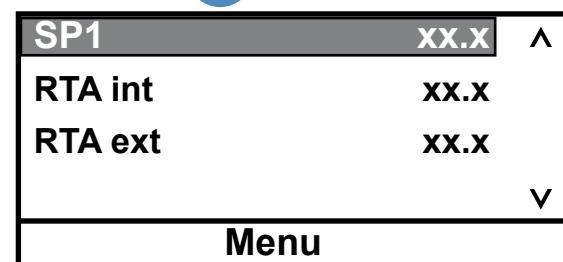The Setpoint and RTA are changed using the same procedure. With the desired setpoint highlighted press  to display the submenu.

If this temperature on the Start/Status Displays does not accurately reflect the actual temperature in the reservoir, an RTA can be applied. The RTA can be set  $\pm 10^{\circ}\text{C}$  ( $\pm 18^{\circ}\text{F}$ ).

As an example, if the temperature is stabilized and displaying  $20^{\circ}\text{C}$  but a calibrated reference thermometer reads  $20.5^{\circ}\text{C}$ , the RTA should be set to  $-0.5^{\circ}\text{C}$ . After you enter a RTA value allow display to stabilize before verifying the temperature. **Note** If display accuracy is required, we recommend repeating this procedure at various setpoint temperatures and on a regular basis. ▲

**Note** You cannot adjust the setpoint closer than  $0.1^{\circ}\text{C}$  to either of the fluid's system limits, see Fluids Type in this Section. ▲


- With the desired line highlighted press .

The right-most digit will have a cursor beneath it. Use the left and right arrows to move the cursor to the desired digit and then use the up and down arrows to change the value. Once all the desired changes are made, press  to save the change, or  to cancel it.

**35.0°C**

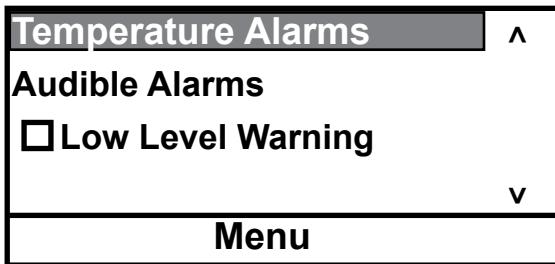**Alarms** is used to view/adjust the high and low temperature alarm limits, to enable/disable the audible alarms and to configure the low level warning reaction.

- With **Alarms** highlighted, press  to display:

- With **Temperature Alarms** highlighted, press  to display:

<b>High Fault</b>	<b>83.0°C</b>	
<b>High Warn</b>	<b>83.0°C</b>	
<b>Low Warn</b>	<b>2.0°C</b>	
<b>Low Fault</b>	<b>2.0°C</b>	

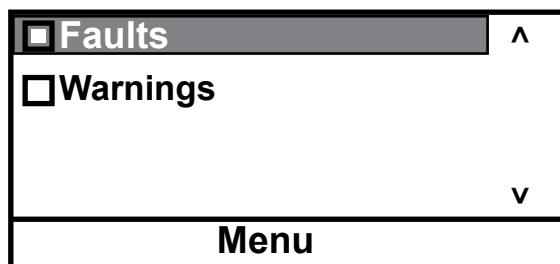
- Highlight the desired limit and press . Follow the same procedure used to change a setpoint. If the Fault temperature is exceeded the circulating bath shuts down and, if enabled, the audible alarm sounds. If the Warn temperature is exceeded the circulating bath continues to run and, if enabled, the audible alarm sounds. In both cases a message is displayed.

**High Fault** cannot be set below **High Warn**.  
**High Warn** cannot be set below **Low Warn**.  
**Low Fault** cannot be set above **High Warn**.

Press , or use the **Menu** window, to save and return to the previous display.

- With **Audible Alarms** highlighted, press  to display the alarms.


Highlight the desired alarm and press  to toggle between enable and disable mode.

If **Faults** is enabled the alarm sounds when a fault occurs. If **Warnings** is enabled the alarm sounds when a warning occurs.

Press , or use the **Menu** window, to save and return to the previous display.

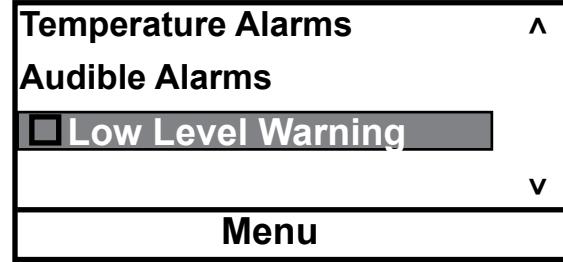- With **Low Level** highlighted, press  to toggle the low level warning alarm on/off:

Press , or use the **Menu** window, to save and return to the previous display.

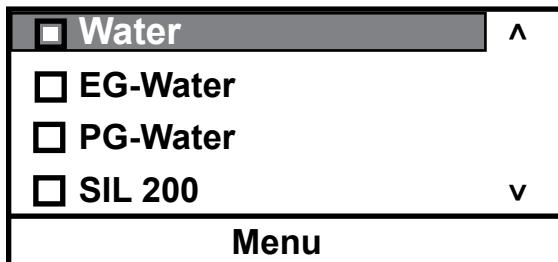
**Fluids Type** is used to identify the type of fluid used. The circulating bath uses the fluid type to automatically set certain operating parameters.

Polyphenylene oxide (PPO) Bath Circulators are used with water only. ▲

1. With **Fluid Type** highlighted, press to display the list of acceptable fluids.

Highlight the desired fluid and press to select it.


2. With the desired fluid selected press , or use the **Menu** window, to save and return to the previous display.

**Pump Adjustment** is used to review/set the desired pump speed.

1. With **Pump Adjustment** highlighted, press to display the speeds.

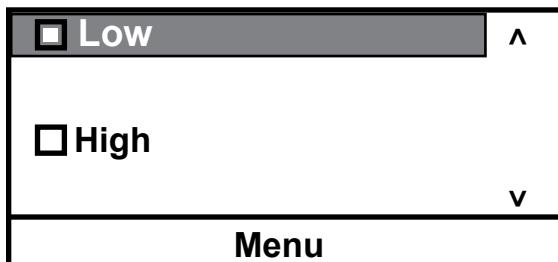Highlight the desired speed and press to select it.

#### Fluid temperature alarm limits

	High °C	Low °C
--	---------	--------

#### All circulating baths:

Distilled Water	+98	+2
DI Water (up to 3 Ω)	+98	+2
EG-Water	+103	-33
PG-Water	+103	-33

#### For Standard Plus and Advanced circulating baths only:

SIL 200	+200	+40
---------	------	-----

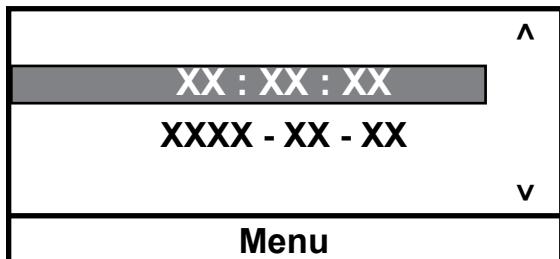
**Use only XIAMETER(r) PMX-200 SILICONE FLUID 50CS. ▲**

#### For Advanced circulating baths only:

SIL 300	+200	+80
SYNTH 260	+200	+40

**Note** The range is also limited by the bath circulator temperature range, see Section 2. ▲

**Set Clock** is used to set the time (hr : min : sec) and date (year - month - day).

**On/Off Timer** is used to enable and set the timer.

- With **On/Off Timer** highlighted, press

After setting the on and off times select **Enable** to activate the timer.

**Auto Restart** is used to enable the auto restart feature. When enabled, the circulating bath will automatically restart after a power failure or power interruption condition. The default setting is **OFF**. Note Consider any possible risks before enabling this mode of operation. ▲

- With **Auto Restart** highlighted, press

**Energy Saving** is used to enable the energy savings mode. The Energy Saving mode is primarily designed for applications running under a stable load. Enabling the mode saves energy by reducing the circulator's heater power and cooling requirements. This can result in substantial energy savings over the life of the circulator. The default setting is **ON**.

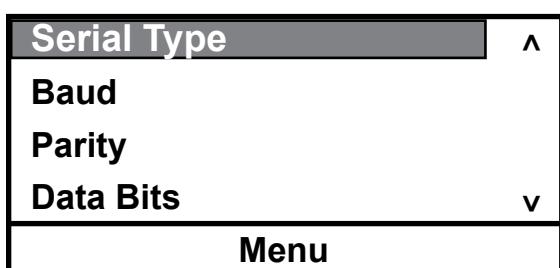
- With **Energy Saving** highlighted, press

- With optional **Remote Sensor** highlighted, press

**Interfaces** (optional) is used to enable/configure the optional serial communications feature.

- With **Interfaces** highlighted, press

Highlight the desired parameter and press

#### Available options:

**Serial Type** RS232, RS485, Off

**Protocol** Standard, NC

**Baud** 19200, 9600, 4800, 2400, 1200, 600, 300

**Parity** None, Odd, Even

**Data Bits** 8

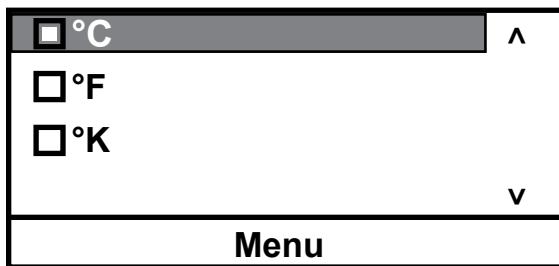
**Stop Bits** 1, 2

See the Appendix for additional information.

**Settings - Display Options** is used to view/adjust the circulating bath's Temperature Units, the Temperature Resolution, the displayed Language, and the Display Contrast and Display Delay.


1. With **Temp. Unit** highlighted press . Use the up/down arrows to highlight the desired temperature scale.

Press .

3. With **Language** highlighted press . Use the up/down arrows to highlight the desired language.

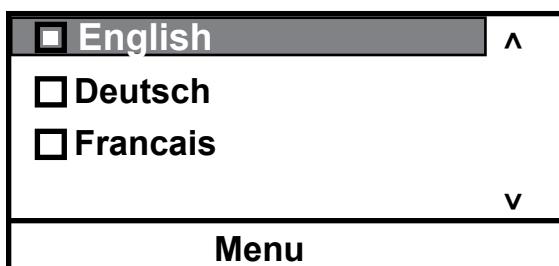Press .


2. With **Temp. Resolution** highlighted press .

Use the up/down arrows to highlight the desired resolution. Press .

4. With **Display Contrast** highlighted press .

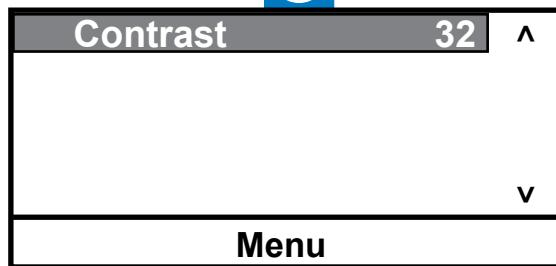Press  again and use the up/down arrows keys to change the contrast. With the desired contrast showing, press  again.


**Note** Holding **Esc** for five seconds resets the display contrast to the default level and also brings up the language menu to change, if needed, the displayed language. ▲

5. With **Display Delay** highlighted press  to enable/disable it.

Use the up/down arrows to highlight the time and press  again.


Use the up/down arrows to change the value. Once the desired delay is displayed press .

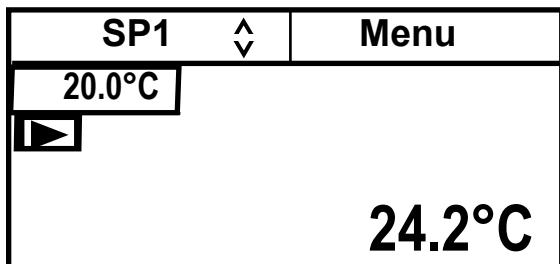With **Delay** enabled and the Start Display showing, if no arrows keys on the circulator are pressed the Start Display will change to the Status Display after the delay expires, in this example after 60 seconds.

Start Display

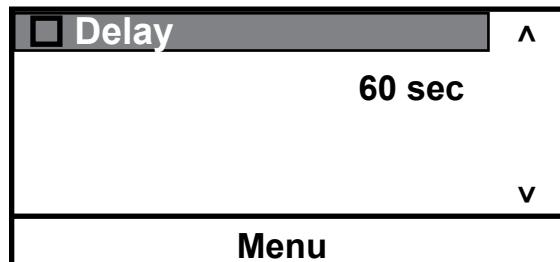If desired, press  to toggle between the Start/Status Displays.


With **Display Delay** enabled and the Start Display showing, if no arrows keys on the circulator are pressed the Start Display will change to the Status Display after the delay expires, see pages 4-2 and 4-3.

Status Display


**System Messages** is used to view any Warning or Fault messages.

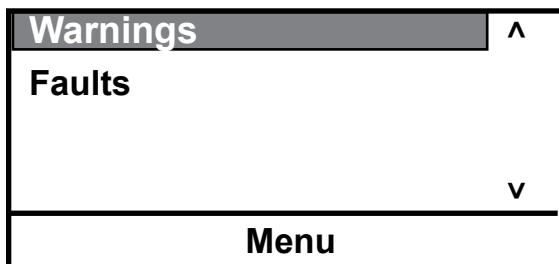- With **Messages** highlighted, press  to display the options.


**System Run Time** is used to view the circulating bath (**Unit**) and pump operating hours.

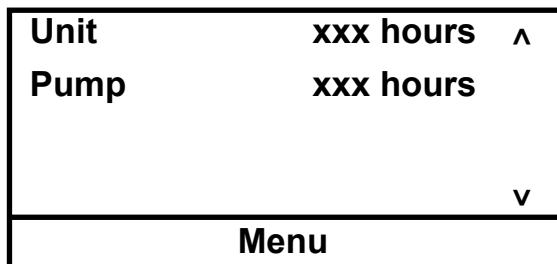- With **Run Time** highlighted, press  to display the times.

**System Configuration** is used to view the circulating bath's configuration.

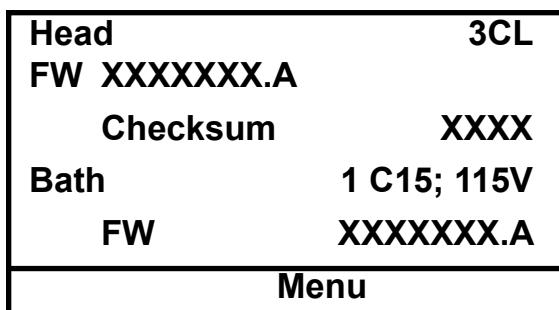- With **Configuration** highlighted, press  to display the settings.

Note Heating baths display  
**Bath Unknown.** ▲

**System - Password/Reset** is used only by a qualified technician. Changing the password enables circulator reset options, the temperature sensor calibration procedure and displays PID values.

1. With **Password/Reset** highlighted, press  to display:

Level	User	^
<b>Password</b>	0	
		v
Menu		

2. Press  and change the number to **1**.

Level	User	^
<b>Password</b>	<b>1</b>	
		v
Menu		

3. Press  to display:

Level	Operator	^
<b>Password</b>	<b>1</b>	
<b>Reset</b>		
<b>Calibration</b>		v
Menu		

Scroll down to display **PID Tuning**, see page 4-15.

1. If desired, highlight **Reset** and press  to display:

<b>Reset user settings</b>	^
<b>Reset PID settings</b>	
<b>Reset both</b>	
	v
Menu	

Highlight the desired reset option and press .

A confirmation message appears, press  again.

**Note** The circulating bath resets to the **User** mode whenever it is turned off. It also resets to the **User** mode whenever the Start/Status Display is displayed continuously for 10 minutes. ▲

**Note** Ensure the RTA is set to 0 before doing a calibration. ▲

- To calibrate the temperature sensor highlight **Calibration** and press  to display:

<b>Internal RTD</b>	^
<b>External RTD</b>	v
<hr/>	

- With **Calibrate** highlighted, press  to display:

<b>Internal RTD</b>	<b>xx.x</b> ^
<b>High</b>	<b>xx.x</b>
<b>Low</b>	<b>xx.x</b>
<b>Calibrate</b>	<b>SP xx.x</b> v
<hr/>	
<b>Menu</b>	

- Press  and depending on which end you are calibrating highlight **High** or **Low**. Once the **Internal RTD** temperature stabilizes enter the temperature displayed on your reference thermometer and press .

<b>Internal RTD</b>	<b>xx.x</b> ^
<b>High</b>	<b>xx.x</b>
<b>Low</b>	<b>xx.x</b>
<b>Calibrate</b>	<b>SP xx.x</b> v
<hr/>	
<b>Menu</b>	

- Repeat for the other calibration point.

<b>Internal RTD</b>	<b>xx.x</b> ^
<b>High</b>	<b>xx.x</b>
<b>Low</b>	<b>xx.x</b>
<b>Calibrate</b>	<b>SP xx.x</b> v
<hr/>	
<b>Menu</b>	

- With the desired sensor highlighted, press  to display:

<b>Calibrate</b>	^
<b>Restore User Cal</b>	
<b>Save User Cal</b>	
<b>Restore Factory Cal</b>	v
<hr/>	
<b>Menu</b>	

- Highlight the **SP** temperature box and enter either the desired high or low setpoint value.

<b>Internal RTD</b>	<b>xx.x</b> ^
<b>High</b>	<b>xx.x</b>
<b>Low</b>	<b>xx.x</b>
<b>Calibrate</b>	<b>SP xx.x</b> v
<hr/>	
<b>Menu</b>	

- Highlight **Calibrate** and the press  to complete the procedure.


<b>Internal RTD</b>	<b>xx.x</b> ^
<b>High</b>	<b>xx.x</b>
<b>Low</b>	<b>xx.x</b>
<b>Calibrate</b>	<b>SP xx.x</b> v
<hr/>	
<b>Menu</b>	

Once the calibration is complete you can store it into memory by selecting **Save User Cal** and pressing .


You can later restore the same calibration by highlighting **Restore User Cal** and pressing .

Another option is to restore the factory calibration values by highlighting **Restore Factory Cal** and pressing .

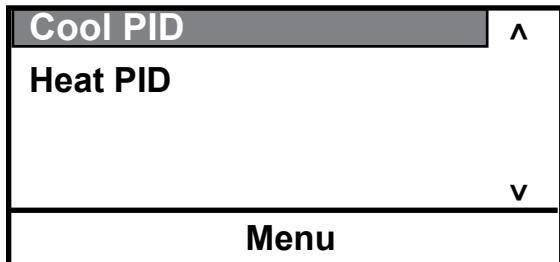
1. With **PID Tuning** highlighted, press  to display:

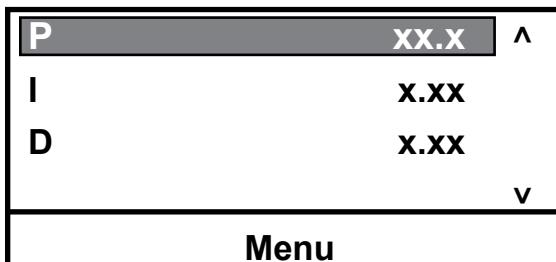2. Highlight the desired PID and press  to display:

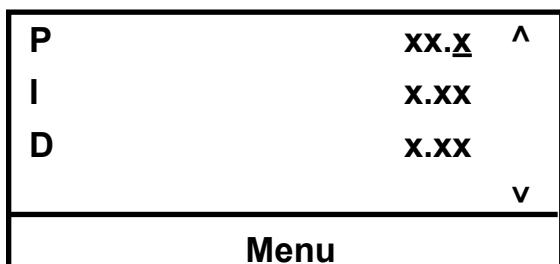3. If required, press  to change the value.


Factory values are:

P = 05.0

I = 0.10

D = 0.00

## High Temperature Cutout

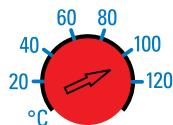HTC (Temperature range varies with type of recirculating bath.)

To protect your application, the adjustable High Temperature Cutout (HTC) ensures the heater will not exceed temperatures that can cause serious damage. A temperature sensor is located in the bath fluid. A High Temperature fault occurs when the temperature of the sensor exceeds the set temperature limit.

In the event of a fault the circulating bath shuts down and displays a fault message, see Section 6. The cause of the fault must be identified and corrected before the circulating bath can be restarted.

The cutout is not preset and must be adjusted. To set the cutout, with the circulating bath off use a flat head screwdriver to turn the red dial fully clockwise. Start the circulating bath. Adjust the setpoint for a few degrees higher than the highest desired fluid temperature and allow the bath to stabilize at the temperature setpoint. Then slowly turn the dial counterclockwise until the circulating bath shuts down and the message appears. Press  to clear the message.

Before you can restart the circulating bath it has to cool down a few degrees. To restart the circulating bath press the black ring surrounding the red dial and then press  again.


If Auto Restart is enabled the circulating bath restarts, if disabled use the Start Up procedure.

**Note** We recommend periodically rechecking operation or if the circulating bath is moved. ▲

## Stopping the Circulating Bath

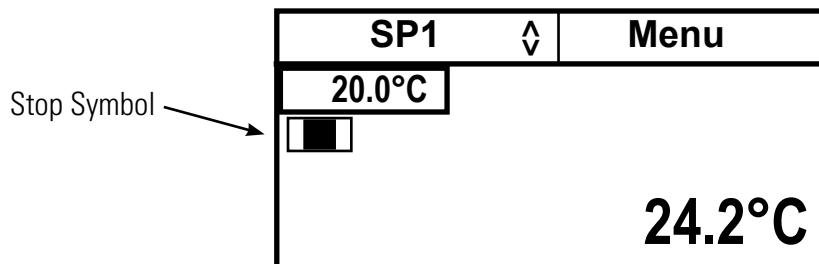Ensure the stop symbol is highlighted, if not use the arrow keys to navigate to the symbol.

Press . The circulating bath stops and the stop symbol turns into a start symbol ().

## Power Down

Press . The screen will go blank.

## Shut Down

Place the circuit protector to the  position.

Using any other means to shut the circulating bath down can reduce the life of the compressor. ▲

Always turn the circulating bath off and disconnect it from its supply voltage before moving. ▲

The circuit protector is not intended to act as a disconnecting means. ▲

## Restarting

**Note** When quickly restarting, the compressor may take up to 10 minutes before it starts to operate. ▲

## Decommissioning/ Disposal

**WARNING**

Decommissioning prepares equipment for safe and secure transportation.

**Laboratory Grade Ethylene glycol (EG) is poisonous and flammable. Before disposing refer to the manufacturer's most current SDS for handling precautions. ▲**

**CAUTION**

**Decommissioning must be performed only by qualified dealer using certified equipment. All prevailing regulations must be followed. ▲**

Consider decommissioning the circulating bath when:

- It fails to maintain desired specifications
- It no longer meets safety standards
- It is beyond repair for its age and worth

**Note** Keep in mind any impact your application may have had on the circulating bath. ▲

Direct questions about decommissioning or disposal to our Sales, Service and Customer Support.

## Storage

**If the bath circulator is to be transported and/or stored in cold temperatures it needs to be drained and then flushed with a 50/50 laboratory grade glycol/water mixture. ▲**

The bath circulator can be stored for up to 90 days inside the temperature range of -25°C to +60°C (-13°F to +140°F).

If necessary, when removing the bath circulator from storage allow it time to warm up and dry out in order to prevent any condensation issues.

## Section 5 Preventive Maintenance

### **WARNING**

Laboratory Grade Ethylene glycol (EG) is poisonous and flammable. Before performing any preventive maintenance refer to the manufacturer's most current SDS for handling precautions. ▲

### **CAUTION**

Disconnect the power cord prior to performing any maintenance. ▲

Handle the circulating bath with care, sudden jolts or drops can damage its components. ▲

There are no user serviceable components within the equipment panels.

### **Cleaning**

After time, the circulating bath's stainless steel surfaces may show spots and become tarnished. Normal stainless steel cleaners can be used.

Clean the reservoir and built-in components at least every time the bath liquid is changed. Use water and a soft cloth.

### **CAUTION**

The inside of the circulating bath must be kept clean in order to ensure a long service life. Quickly remove substances containing acidic or alkaline substances and metal shavings as they could harm the surfaces causing corrosion. If corrosion (e.g., small rust marks) occur in spite of this, cleaning with stainless steel caustic agents has proved to be suitable. Apply these substances according to the manufacturer's recommendations.

Before cleaning, to protect labels, the nameplate, electrical connections, painted and plastic surfaces and to prevent the cleaning agent from entering through any vent openings, mask off all areas except the reservoir.

### **CAUTION**

### **Condenser Fins**

Wear protective clothing and take appropriate measures when handling the cleaning agent.

For cooling/heating circulating baths, in order to maintain cooling capacity, clean the fins two to four times per year, depending on the operating environment.

Switch off the circulating bath and unplug the power cord.

Clean the fins with compressed air.

For extreme soiling a qualified technician will need to remove the cooling compressor casing.

## Hoses

Inspect and tighten the external hoses and clamps daily.

## Testing the Safety Features

### High temperature protection

Set a cut-off temperature that is lower than the desired setpoint temperature.

Switch on the circulating bath and ensure it shuts down at the set cut-off temperature.

If the circulating bath does not shut down have it checked by a qualified technician.

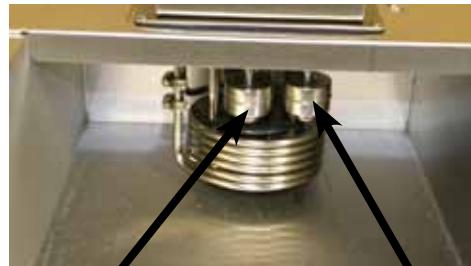
Reset the safety to the desired temperature.

### Low liquid level protection

With the circulating bath on, use a screwdriver and slowly push down on each sensor until the error message indicated below appears.

See Section 6 for details on error messages.

If either message does not appear, have the circulating bath checked by a qualified technician.

FAULT:

HTC/Low Level Fault

PRESS ENTER


FAULT:  
LLC

Low Level Cutout  
PRESS ENTER

## Section 6 Troubleshooting
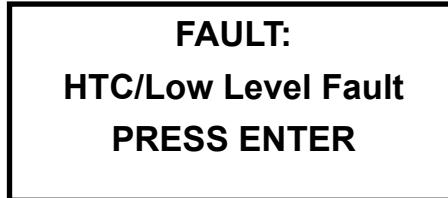
### Error Displays

Error messages are cleared by pressing the enter key (  ). Once the cause of the error message is identified and corrected, to restart the circulating bath - on circulating bath with an HTC error press the black ring surrounding the red dial - and then for all circulating bath press  again. If **Auto start** is enabled the circulating bath will restart, if disabled use the Start Up procedure.

Error Message (Typical)

#### Fault Displays

The heating element, pump and, if applicable, refrigeration shut down with a fault. A fault also sounds the alarm, if enabled.

Message	Action
<b>High Fixed Temp.</b>	<ul style="list-style-type: none"><li>• circulating bath's nonadjustable high temperature protection limit exceeded</li><li>• check fluid selection</li><li>• check environmental conditions</li></ul>
<b>High Temperature</b>	<ul style="list-style-type: none"><li>• adjustable high temperature protection limit exceeded</li><li>• check limit setting</li><li>• check fluid selection</li><li>• ensure circulating bath has adequate ventilation</li></ul>
<b>High Temperature Refrigeration</b>	<ul style="list-style-type: none"><li>• check voltage supply</li><li>• the refrigeration may need servicing</li></ul>
<b>HPC High Press. Cutout</b>	<ul style="list-style-type: none"><li>• check for obstructions to air flow</li><li>• the refrigeration may need servicing</li></ul>
<b>HTC/Low Level Fault</b>	<ul style="list-style-type: none"><li>• if the fluid level is greater than approximately 5.5 cm below the reservoir top it is a low level fault. Fill fluid to proper level.</li><li>• if not, high temperature protection limit exceeded</li><li>• check limit setting</li><li>• allow circulating bath to cool down to at least 10°C cooler than the HTC setting</li><li>• reset HTC by pressing black ring</li><li>• set HTC to desired setting</li></ul>

<b>LLC Low Level Cutout</b>	<ul style="list-style-type: none"> <li>• reservoir fluid level too low for safe operation</li> <li>• check fluid level</li> <li>• check for leaks</li> </ul>
<b>Low Fixed Temp.</b>	<ul style="list-style-type: none"> <li>• circulating bath's nonadjustable low temperature protection limit exceeded</li> <li>• check fluid selection</li> </ul>
<b>Low Temperature</b>	<ul style="list-style-type: none"> <li>• adjustable high temperature protection limit exceeded</li> <li>• check limit setting</li> <li>• check fluid selection</li> </ul>
<b>Motor Fault</b>	<ul style="list-style-type: none"> <li>• it can take over 10 minutes for the motor temperature to get low enough before the circulating bath can be restarted</li> </ul>
<b>MOL Motor Overload</b>	<ul style="list-style-type: none"> <li>• allow circulating bath to cool down</li> </ul>
<b>Open RTD1 Internal</b>	<ul style="list-style-type: none"> <li>• open internal temperature sensor</li> </ul>
<b>Open RTD2 External (Advanced only)</b>	<ul style="list-style-type: none"> <li>• open external temperature sensor</li> </ul>
<b>Shorted RTD1 Internal</b>	<ul style="list-style-type: none"> <li>• shorted internal temperature sensor</li> </ul>
<b>Shorted RTD2 External (Advanced only)</b>	<ul style="list-style-type: none"> <li>• shorted external temperature sensor</li> </ul>

### Warning Displays

The circulating bath will continue to run with a warning. A warning also sounds the alarm, if enabled.

<b>Bad Calibration</b>	<ul style="list-style-type: none"> <li>• redo calibration</li> </ul>
<b>High Temperature</b>	<ul style="list-style-type: none"> <li>• adjustable high temperature protection limit exceeded</li> <li>• check limit setting</li> <li>• check fluid selection</li> </ul>
<b>Low Level</b>	<ul style="list-style-type: none"> <li>• reservoir fluid level too low for safe operation</li> <li>• check fluid level</li> </ul>
<b>Low Temperature</b>	<ul style="list-style-type: none"> <li>• adjustable low temperature protection limit exceeded</li> <li>• check limit setting</li> <li>• check fluid selection</li> </ul>

## Checklist

### Circulating bath will not start

Check for error codes, see Error Codes in this section.

Ensure the circuit protector is in the on (I) position.

Make sure supply voltage is connected and matches the nameplate rating  $\pm 10\%$

### No display

Recycle the circulating bath's circuit protector.

### Circulating bath will not circulate process fluid

Check the reservoir level. Fill, if necessary.

Check the application for restrictions in the cooling lines.

The pump motor overloaded. The pump's internal overtemperature overcurrent device will shut off the pump causing the flow to stop. This can be caused by low fluid, debris in system, operating circulating bath in a high ambient temperature condition or excessively confined space. Allow time for the motor to cool down.

Make sure supply voltage matches the nameplate rating  $\pm 10\%$ .

### Inadequate temperature control

Verify the setpoint.

On cooling/heating circulating baths, make sure the condenser is free of dust and debris.

Check the fluid concentration.

Ensure installation complies with the site requirements in Section 3.

Make sure supply voltage matches nameplate rating  $\pm 10\%$ .

If the temperature continues to rise, make sure your application's heat load does not exceed the rated specifications.

Check for high thermal gradients (e.g., the application load is being turned on and off or rapidly changing).

### Circulating bath shuts down

Ensure  button wasn't accidentally pressed.

Ensure the circuit protector is in the on (I) position.

Check for error codes.

Make sure supply voltage is connected and matches the nameplate rating  $\pm 10\%$ .

Restart the circulating bath.

### **USB Driver Not Recognized**

If your operating system does not automatically recognize the optional driver log on to:

<http://www.ftdichip.com/FTDrivers.htm>

for instructions.

# Appendix AC Serial Communications Protocol

Serial communication is accomplished through the USB port on the chiller. If your operating system does not automatically recognize the optional driver log on to: <http://www.ftdichip.com/FTDrivers.htm> for instructions.

**Note** This appendix assumes you have a basic understanding of communications protocols. Information on the NC, STANDARD and NAMUR protocols is available upon request. ▲

**Note** Keypad operation is still available with serial communications enabled. ▲

All commands must be entered in the exact format shown in the tables on the following pages. The tables show all commands available, their format and responses. Chiller responses are either the requested data or an error message. The chiller response must be received before the host sends the next command.

The host sends a command embedded in a single communications packet, then waits for the chiller's response. If the command is not understood, the chiller responds with an error command. Otherwise, the chiller responds with the requested data.

Commands are not case sensitive. Upper or lower case letters may be used. Commands are listed in the Commands Table, error responses are given in the Errors Table, and symbols are shown in the Key Table.

Key	
Symbol	Meaming
[B]	A binary value 0 or 1 (0 = Off, FALSE or Disable(d); 1 = On, TRUE or Enable(d)).
[CR]	Carriage return – used as the termination character.
[U]	Text representing the units associated with a value.
[V]	A value that can be requested in a read command or sent as part of a set command.
[V <sub>MAX</sub> ]	Maximum allowed value. Part of error message when set value is too high.
[V <sub>MIN</sub> ]	Minimum allowed value. Part of error message when set value is too low.

**Value:** Read commands return analog [V] or bit [B] values or settings, while set commands send analog or bit settings. Read commands return values with the same displayed precision. Set command messages missing the space character between the command and the setting will be rejected, as the user's intent is unclear.

**Units:** A read command returning an analog [V] value or setting, will include the units [U] associated with that value or setting. A set command sending an analog value will not include the units. The units returned by the complementary read command are assumed.

**Termination character:** A carriage return [CR] is used to terminate command and response messages. (Typically the "Enter" key on the keyboard.)

**Note** The inter-character timeout (time between transmitted characters) is set to 30 seconds. Exceeding the timeout will clear the receiver buffer and require the message to be retransmitted.

**Note** Special characters (backspace, delete, insert, etc.) are not recognized by the protocol and generate error responses.

## Commands Table:

Commands		<i>All messages from master and slave are terminated with a carriage return [CR]</i>		
Command Description	Notes	Master Sends	Sample Slave Response <i>(echo off)</i>	<i>Alternate units</i>
Read Temperature	<i>Internal</i>	RT	[V]C	F K
Read Temperature 2	<i>External</i>	RT2	[V]C	F K
Read Displayed Setpoint		RS	[V]C	F K
Read Setpoint X (X = 1 to 5)		RSX	[V]C	F K
Read High Temperature Fault		RHTF	[V]C	F K
Read High Temperature Warn		RHTW	[V]C	F K
Read Low Temperature Fault		RLTF	[V]C	F K
Read Low Temperature Warn		RLTW	[V]C	F K
Read Proportional Heat Band Setting		RPH	[V]%	
Read Proportional Cool Band Setting		RPC	[V]%	
Read Integral Heat Band Setting		RIH	[V]Repeats per minute	
Read Integral Cool Band Setting		RIC	[V]Repeats per minute	
Read Derivative Heat Band Setting		RDH	[V]Minutes	
Read Derivative Cool Band Setting		RDC	[V]Minutes	
Read Temperature Precision		RTP	[V]	
Read Temperature Units		RTU	[V}	C,F,K
Read Unit On		RO	[B]	
Read External Probe Enabled		RE	[B]	
Read Auto Restart Enabled		RAR	[B]	
Read Energy Saving Mode		REN	[B]	
Read Time		RCK	hh:mm:ss	
Read Date		RDT	mm/dd/yyyy or dd/mm/yyyy	
Read Date Format		RDF	mm/dd/yyyy or dd/mm/yyyy	

<b>Commands</b>		<i>All messages from master and slave are terminated with a carriage return [CR]</i>	
Command Description	Notes	Master Sends	Sample Slave Response <i>(echo off)</i> <i>Alternate units</i>
Read Firmware Version		RVER	[V]
Read Firmware Checksum		RSUM	[V]
Set Displayed Setpoint		SS [V]	OK
Set Setpoint X (X = 1 to 5)		SSX [V]	OK
Set High Temperature Fault		SHTF [V]	OK
Set High Temperature Warning		SHTW [V]	OK
Set Low Temperature Fault		SLTF [V]	OK
Set Low Temperature Warning		SLTW [V]	OK
Set Proportional Heat Band Setting		SPH [V]	OK
Set Proportional Cool Band Setting		SPC [V]	OK
Set Integral Heat Band Setting		SIH [V]	OK
Set Integral Cool Band Setting		SIC [V]	OK
Set Derivative Heat Band Setting		SDH [V]	OK
Set Derivative Cool Band Setting		SDC [V]	OK
Set Temperature Resolution		STR [V]	OK
Set Temperature Units		STU [V] C,F,K	OK
Set Unit On Status		SO [B]	OK
Set External Probe On Status		SE [B]	OK
Set Auto Restart Enabled		SAR [B]	OK
Set Energy Saving Mode		SEN [V]	OK
Set Pump Speed		SPS [V] L, H	OK

Errors Table:

Errors		
Error Description	Notes	Slave Responds
<i>Not defined, not implemented or incorrectly formatted</i>		? Unsupported command
<i>Extra characters...</i>		? Format error
<i>Set value too high</i>		? Maximum allowed is $[V_{\text{MAX}}]$
<i>Set value too low</i>		? Minimum allowed is $[V_{\text{MIN}}]$
<i>Argument to binary set command not 0 or 1</i>		? Value must be 0 or 1
<i>Set command attempted while in read only mode</i>		? Mode is read only
<i>Set command failed (e.g. SO 1 with low level)</i>		? Failed

## RUFS Read Unit Fault Status

This command returns 5 values. These are decimal representations of hexadecimal values. Each individual bit of the value represents a different warning, fault or status.

decimal	hex	B7	B6	B5	B4	B3	B2	B1	B0
1	1	0	0	0	0	0	0	0	1
2	2	0	0	0	0	0	0	1	0
4	4	0	0	0	0	0	1	0	0
8	8	0	0	0	0	1	0	0	0
16	10	0	0	0	1	0	0	0	0
32	20	0	0	1	0	0	0	0	0
64	40	0	1	0	0	0	0	0	0
128	80	1	0	0	0	0	0	0	0

Value	Description of bits
$V_1$	B0 – B5 unused B6 rtd1 shorted B7 rtd1 open
$V_2$	B0 HTC fault B1 high RA temperature fault B2 – B7 unused
$V_3$	B0 low level warn B1 lo temperature warn B2 high temperature warn B3 low level fault B4 low temperature fault B5 high temperature fault B6 low temperature fixed fault B7 high temperature fixed fault
$V_4$	B0 PWM heat duty cycle > 0 B1 compressor On/Off B2 Pump On status B3 Controller On status B4 Controller Stopping B5 Controller fault status B6 unused B7 Beeper On status
$V_5$	B0 Pump speed fault B1 MOL fault B2 HPC fault B3 Cool Icon On steady (controller is cooling at max capacity) B4 Cool Icon flashing B5 Heat Icon On steady (controller is heating at max capacity) B6 Heat Icon flashing B7 External sensor controlling

Refer to Key table on page 1 for explanation of symbols and their meanings.

Examples:

Read Temperature:

Host

R	T		<b>CR</b>
Command		[CR]	

Controller:

2	0	.	0	C	<b>CR</b>
	[V]			[U]	[CR]

Set Setpoint:

Host

S	S		2	0	<b>CR</b>
Command			[V]	[CR]	

Controller:

O	K	<b>CR</b>
Command		[CR]
Accepted		

Read Temperature 2:

Host:	R	T	2	<b>CR</b>		
Controller:	2	0	.	0	C	[CR]

Set Setpoint to -22°C when minimum allowed is -20°C: Minimum allowed is [V<sub>MIN</sub>]

Host:	S	S		-	2	2	<b>CR</b>															
Controller:	?		M	i	n	i	m	u	m	a	1	1	o	w	e	d	i	s	-	2	0	<b>CR</b>

## Optional Serial Communications Adapter

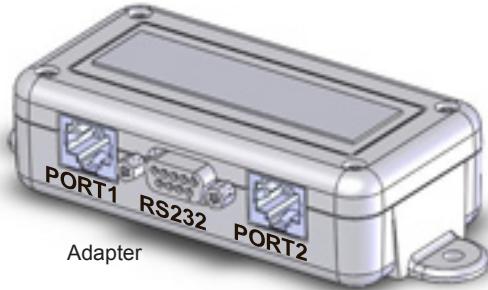
### Tools required:

- None

### Procedure:


Turn off the circulator before installing the adapter. ▲

Pin #	Function	Pin #	Function
1	No connection	1-7	No connection
2	TX	8	T+
3	RX	9	T-
4	No connection		
5	GND = Signal ground		
6 - 9	No connection		

TX = Transmitted data from circulator

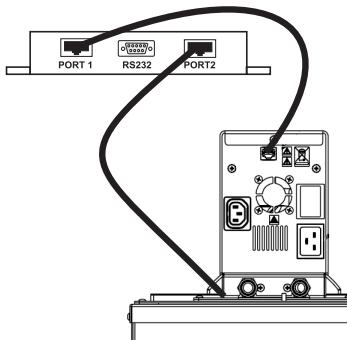RX = Received data to circulator

1. If the circulator already has a communication cable installed, remove the cable from the rear of the circulator and plug that cable into PORT 2 on the adapter.
2. Plug the supplied cable into PORT 1 on the adapter and the other end into the circulator. Regardless of the configuration, the supplied cable *always* goes from the circulator to PORT 1.
3. Plug the supplied serial communications cable into the communication port on the adapter and then the other end into your computer.
4. If desired, use the supplied Velcro® tape to attach the adapter to a convenient location on the circulator.