

Model 365510 Digital Stopwatch

OPERATION

1. NORMAL MODE

Time: Hours, minutes, and seconds displayed

Alarm: Press & hold the SPLIT/RESET button to display alarm time.

Alarm ON/OFF: Bell symbol ON (right-hand side of the LCD) when alarm is set

Calendar: Month/Date displayed when START/STOP button is pressed and held.

2. STOPWATCH MODE (Press MODE once to access this mode).

Accumulated Elapsed Timer:

1. Press Start/Stop to Start
2. Press Start/Stop to Stop
3. Press Start/Stop Restart
4. Press Start/Stop to Stop
5. Press Split/Reset to reset the display

Split Time Measurement Steps

1. Press Start/Stop to Start
2. Press Split/Reset to Split
3. Press Split/Reset to release Split
4. Press Start/Stop to Stop
5. Press Split/Reset to reset the display

Two Competitor Measurement

1. Press Start/Stop to Start
2. Press Split/Reset to Split
3. Press Start/Stop to Stop
4. Press Split/Reset to release Split
5. Press Split/Reset to reset the display

3. SET THE CLOCK/CALENDAR (Press MODE three times to access)

Press SPLIT/RESET to select the digit for editing (the selected digit will flash). Press START/STOP to reset the digit to zero or keep the button depressed for quick scrolling to other values. Press MODE to return to normal operation.

4. SET THE ALARM TIME (Press MODE twice to access this mode)

Press STOP/START to advance the hour then press SPLIT/RESET to select minutes then press STOP/START to advance the minutes. Press MODE to return to normal operation. Press and hold the SPLIT/RESET and START/STOP buttons to arm or dis-arm the alarm.

5. SET THE SNOOZE FUNCTION

When the alarm sounds press START/STOP to go to the 5-minute snooze mode. To silence the alarm without the snooze mode, press SPLIT/RESET after the alarm sounds.

6. SET THE HOUR CHIME

Press and hold SPLIT/RESET then press START/STOP while continuing to hold the SPLIT/RESET button to toggle between hourly chime ON and OFF.

7. BATTERY REPLACEMENT

This Stopwatch uses an LR-44 / A-76 alkaline button battery. Qualified personnel must replace the battery. Battery life is typically one year.

WARRANTY

EXTECH INSTRUMENTS CORPORATION warrants this instrument to be free of defects in parts and workmanship for one year from date of shipment (a six month limited warranty applies on sensors and cables). If it should become necessary to return the instrument for service during or beyond the warranty period, contact the Customer Service Department at (781) 890-7440 ext. 210 for authorization. A Return Authorization (RA) number must be issued before any product is returned to Extech. The sender is responsible for shipping charges, freight, insurance and proper packaging to prevent damage in transit. This warranty does not apply to defects resulting from action of the user such as misuse, improper wiring, operation outside of specification, improper maintenance or repair, or unauthorized modification. Extech specifically disclaims any implied warranties or merchantability or fitness for a specific purpose and will not be liable for any direct, indirect, incidental or consequential damages. Extech's total liability is limited to repair or replacement of the product. The warranty set forth above is inclusive and no other warranty, whether written or oral, is expressed or implied.

Copyright © 2001 Extech Instruments Corporation. All rights reserved
including the right of reproduction in whole or in part in any form.



Tech Support Hotlines

781-890-7440 ext. 200

extech@extech.com

www.extech.com