Cole-Parmer®

Digital Reversing Mixer

Language Index

English	1
French	13
Spanish.....	25
German.....	37

Contents

Packing List	2
Introduction	2
Safety Label Legend	2
Safety Instructions	3
General Instructions	4
Approvals	4
Installation	5
Environmental Operating Conditions	5
Auto Start.....	6
Torque Capabilities	6
Dimensions	7
Technical Data.....	8
General Operating Instructions.....	8
Cleaning and Service	11
Warranty and Product Liability	11

Packing List

Compare the contents of the shipment to the items on the following list to ensure all parts are received with the mixer. Do not discard the container and packing materials until all parts are accounted for. It is recommended that the original packing materials be kept for transportation and storage.

50004-00

- Digital Reversing Mixer with chuck
- Power supply
- Support Rod (including cap screw and lock washer)
- Chuck Key
- Instruction manual

50004-10 Contents of Digital Reversing Mixer System

- 50004-00 Digital Reversing Mixer
- 04561-24 Clamp
- 50801-50 Stand
- 50801-62 Pitched Blade Propeller

Introduction

This Mixer is specifically designed for professional use in the laboratory and for use in industry to mix liquids with various properties and applications.

This Mixer uses a high efficiency brushless DC motor. The speed is kept constant with continuous feedback to the motor. When torque demands are outside the Mixer's range, integrated electronic overload protection ensures no damage can occur to the electronics, motor or transmission.

The mixer is designed for reliability and durability. There are no wear parts that require user service during the normal life of the mixer. There are no user serviceable parts.


Safety Label Legend

CAUTION: Read operating instructions carefully and in full.

RECYCLE: Recycle in accordance with the waste electronic and electrical equipment (WEEE) directive.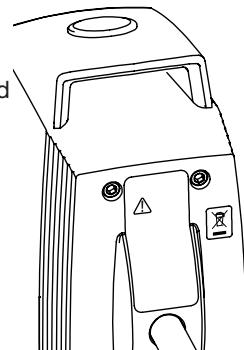

Safety Instructions

-  1. All operators must be familiar with the mixer and read this entire manual. If the mixer is used in a manner not specified by the manufacturer, the protection provided by the equipment may be impaired.
-  2. The mixer must be securely fixed to a stable support. Mount to a stand (50801-50) using a heavy duty clamp (04561-24). If other stands or clamps are used, the mixer must be secured such that it will not fall if tipped.
-  3. **CAUTION:** This is not an explosion proof mixer. Do not use in a highly flammable or explosive environment.
-  4. Spinning paddles or impellers can cause severe personal injuries. Operators must use extreme care and good judgment when mixing at any speed.
-  5. All mixing paddles and impellers must be in good condition with straight shafts. If the mixer vibrates at high speeds check the paddle shaft for damage and repair or replace if needed.
-  6. Extreme care must be taken when mixing chemicals to ensure that no chemicals are splashed outside the mixing vessel. Care must be taken when changing to faster mixing speeds.
-  7. Ensure that the mixing impeller does not contact the containment vessel.
-  8. Do not operate while standing in water. Keep the mixer dry and do not immerse any part, except the mixing paddle into any liquids. Protect from splashing.
-  9. Ensure that no loose clothing, jewelry, or hair can become entangled in any rotating parts. A fast spinning chuck can cause injury to operator.
-  10. Power can be interrupted to the mixer by pressing the power button. In an emergency, disconnect the mains cord.
-  11. Ensure the power button and mains power cord of the mixer can be accessed immediately, directly and without risk at any time, and above all, in case of fault.
-  12. The mixer must be used with supplied power cord. The cord and grounded plug type are important safety elements. Do not use an inadequately rated cord or ungrounded plug. Always connect the mixer to a grounded outlet. Do not alter the plug.

General Instructions

Read Safety Instructions: Unpack the mixer carefully. Ensure the mixer is not damaged and verify the contents against the packing list. If the mixer is damaged or the contents do not correspond with the packing list, notify your supplier immediately.

Read the manual carefully! Make sure that all users read this manual thoroughly before operating the mixer. For further questions or inquiries, please contact:

Cole-Parmer

625 E. Bunker Ct.
Vernon Hills, IL 60061
United States

Tel: 1-800-323-4340
Website: ColeParmer.com
E-mail: sales@antylia.com

Approvals

Standards		
North America	 	CAN/CSA C22.2 No. 61010-1:2012/R:2019-07 CSA-C22.2 No. 61010-2-051:2019 UL 61010-1:2022/R:2019-07 UL 61010-2-051:2019 FCC Part 15, Subpart B:2017 ICES-003:2017
Europe, United Kingdom, Australia, New Zealand	  	IEC 61010-1:2010/A1:2019 IEC 61010-2-051:2015 IEC 61326-1:2012 IEC 61326-1:2013

CSA	Canadian Standards Association
CE	Communauté Européenne (European Community)
UL	Underwriters Laboratories
FCC	Federal Communications Commission
IEC	The International Electrotechnical Commission
ICES	Interference Causing Equipment Standards
EMC	Electromagnetic Compatibility
EMI	Electromagnetic Interference
TÜV	TÜV SÜD
UKCA	UK Conformity Assessed

Installation

Installing the support rod

Remove the hex screw and lock washer from the support rod. Insert the support rod into the back of the mixer so that the holes line up for the hex screw. Secure with hex screw and lock washer and tighten with hex key.

Clamp the mixer to a Cole-Parmer stand on a stable, level surface (see Safety Instructions).

Connect mixer to supplied power supply. Connect power supply to a grounded electrical outlet.

Impeller attachment

Insert the impeller shaft into the open chuck and adjust the impeller to the desired height. The hollow shaft allows the impeller shaft to come through the top of the mixer if necessary. Tighten the chuck's outer ring by hand and then secure with the chuck key.

CAUTION:

 Operating a freely rotating impeller shaft through the top of the mixer is not safe. Do not allow the impeller shaft to touch or extend through the rubber gasket while the motor is powered on. The gasket is for sealing purposes only.

Environmental Operating Conditions

The Reversing Digital mixer must operate in the following conditions:

Indoors

1. Altitudes up to 2000 m (6500 ft).
2. Temperatures from 5°C to 40°C (41°F to 104°F).
3. Maximum relative humidity 80% for temperatures up to 31°C (88°F) decreasing linearly to 50% relative humidity at 40°C (104°F).
4. Pollution degree 2 in accordance with IEC 664.

Auto Start


This mixer includes an optional Auto Start feature. When set, the mixer will resume mixing after a power failure, when power has been removed by control equipment (i.e., PLC) or when the power has been cut (i.e., mixer unplugged). When power is restored, the mixer will automatically ramp up to the last set speed and resume mixing. The mixer ships from the factory with Auto Start disabled.

To set the Auto Start feature first ensure the mixer is powered off. Press and hold the Increase Button then press the Power Button. When active, Auto Start is indicated by flashing LED dots along the bottom of the display.


To disable Auto Start first ensure mixer is powered off. Press and hold the Decrease Button then press the Power Button.

Torque Capabilities

Torque Curve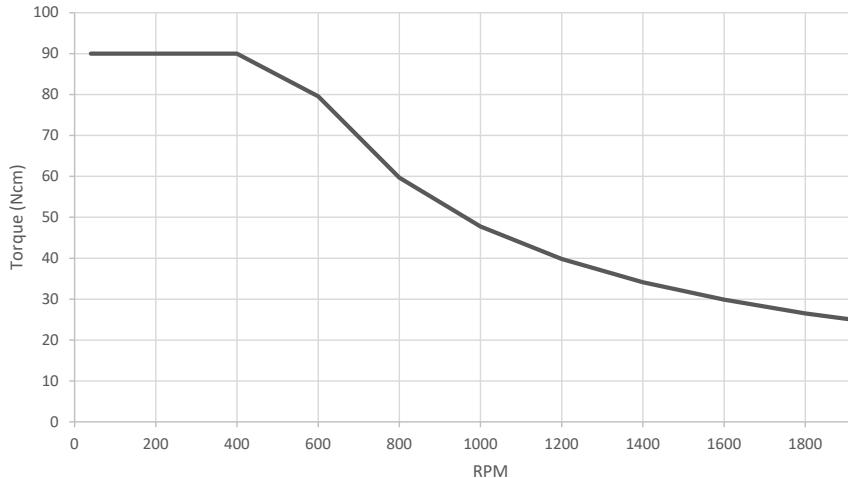

Dimensions

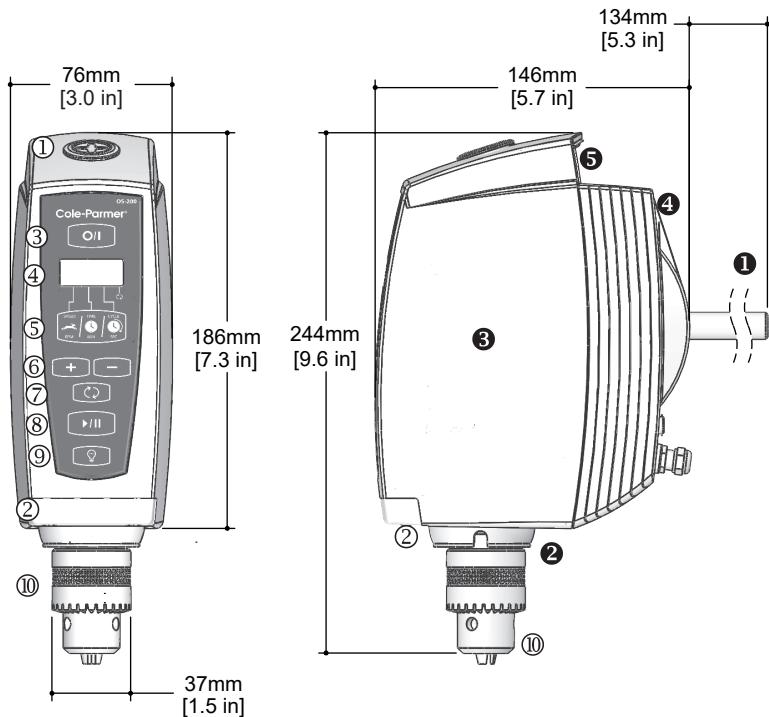- ① Rubber Gasket
 - ② Stir Light
 - ③ Button - Power on/off
 - ④ Four Digit LED Display
 - ⑤ Multifunction Button
 - ⑥ Button - Increase/Decrease rpm
 - ⑦ Button - Reverse Direction
 - ⑧ Button - Start/Pause
 - ⑨ Button - Light on/off
 - ⑩ Chuck
- ① 13mm (1/2 in) dia Support Rod
 - ② Splash Guard
 - ③ Housing - Molded Front
 - ④ Housing - Cast Back
 - ⑤ Ergonomic Carry Handle

Technical Data

Model Number	50004-00
Speed Range	40 - 2010 rpm
Functions	Countdown Timer Clockwise Counter Clockwise xRx Agitation
Speed Readout Accuracy	± 1 rpm or ± 3% of reading, whichever is greater, 4 digits
Maximum Torque	90 Ncm, (127 in-oz)
Input To Power Supply	AC 100-240 volts, 50/60 Hz, 2.5 amps (Must use supplied power supply.)
Electrical Output from Power Supply to Mixer	DC 24 volts, 3.75 amps
Mechanical Output Power	1/15 hp, 50 watts, (Brushless DC Motor)
Weight - Mixer Only	3.0 kg (7 lbs)
Weight - Power Supply	0.6 kg (1.2 lbs)
Maximum Volume	25 litres (6.6 US gallons)
Maximum Viscosity	20,000 cps - e.g. chocolate syrup
Molded Front Housing	Flame retardant, chemical resistant, high impact nylon. Recyclable (PA).
Cast Back Housing	Epoxy coated zinc
Combined Housing	Splash Proof (IP42)
Keypad	Waterproof, chemical resistant polyester.
Stir Light	5 Ultra Bright LEDs, 100,000 hour life.
Power Cord Length	183 cm (6 ft) from mixer to power supply-DC. 183 cm (6 ft) from power supply to wall-AC.
Adjustable Chuck	Fits onto 13 mm (33/64 in) dia output shaft of mixer. Adjusts up to 10 mm (3/8 in).

General Operating Instructions

Make sure the mixer is securely mounted as specified in the installation instructions. Ensure the mixing paddle or impeller is securely attached to the desired position.

The mixer is now ready for use to mix liquids and liquid/solid solutions. Appropriate impeller and container specifications must be determined by the operator, ensuring that all safety instructions are followed.

Keypad Functions

Power Button - Turns the mixer on and off.

Multifunction Button - Sets three parameters - speed (rpm), countdown timer (minutes) and xRx agitation (seconds).

Increase Button - Increases the speed by one rpm or the countdown timer by one minute or the xRx agitation by one second.

Decrease Button - Decreases the speed by one rpm or countdown timer by one minute or the xRx agitation by one second.

Reverse - Reverses the direction that the impeller is turning. The impeller turns clockwise by default.

Start/Pause Button - Allows the mixer to be paused and started again. Use this button to start the mixer when ready.

Light Power Button - Turns the light on and off. The mixer must be plugged in for the light to function.

LED Display - Numeric values reflect speed (rpm), countdown timer (minutes) or xRx agitation (seconds). Solid LED dots indicate which function is selected. Flashing LED dots indicate the Auto Start feature is active (see Auto Start). The LED dots do not act as numeric decimal points.

- ① LED dot (solid) when speed is displayed.
- ② LED dot (solid) when time is displayed.
- ③ LED dot (solid) when cycle time is displayed.
- ④ LED dot (solid) when stirrer is turning counter clockwise.

Clockwise Mixing

Turn the mixer on by pressing the power button. The display will read 40 rpm (or last set speed when Auto Start is enabled) and the speed LED dot will be illuminated. Increase to the desired speed and press the start/pause button. The mixer will start turning in a clockwise direction. The speed LED dot will be illuminated (solid) and the speed value in rpm will be displayed.

Counter Clockwise Mixing

Set by following the clockwise mixing instructions above. Press the reverse button. The reverse button can be pressed while the mixer is running to mix in a counter clockwise direction. The LED dot above the reversing icon will be illuminated (solid) when mixing in a counter clockwise direction.

Countdown Timer

Turn the mixer on by pressing the power button. Set the desired speed by pressing the “+” button. Press the multifunction button until the LED dot for time is illuminated. Press the “+” to set the countdown timer in minutes. Press the start/pause button to start mixing. Press the multifunction button to toggle the display between speed, time and cycle. Mixing stops when time is completed. Three beeps will sound. Note: Countdown Timer value cannot be changed during operation.

xRx Agitation Cycle

Turn the mixer on by pressing the power button. Set speed as above and countdown timer if desired. Press the multifunction button once more until the indicator light for cycle is illuminated. Press the “+” to set the number of seconds the impeller will rotate in one direction before changing directions. Press the start/pause button. Note: xRx Agitation Cycle value cannot be changed during operation.

For example: If set for 5 seconds the mixer impeller will turn clockwise for 5 seconds at the set speed then stop and change to turn counterclockwise for 5 seconds. This mixing pattern will continue until the pause button is pressed or until the time is completed (if set).

The display will show the cycle seconds counting down during each direction unless speed or the countdown timer is selected.

When the mixer is powered on, the default speed is 40 rpm, countdown timer is 0 minutes, and the xRx agitation cycle is 0 seconds. Previously set values will be displayed if Auto Start is enabled.

The mixer will remember the original values set for speed, the countdown timer and the xRx agitation cycle so that when the start/pause button is pressed again the operations will repeat.

Speed can be changed (increased or decreased) while the mixer is running. Press the multifunction button so the speed is displayed. Increase or decrease speed by pressing the “+” or “-” buttons.

The countdown timer and the xRx agitation cycle can be changed only when the mixer is paused and the indicator lights are solid.

The multifunction button can be pressed when the mixer is running to view the status of each function.

For safety, a “soft start” has been incorporated into the mixer. The mixer will ramp up to reach the set speed to minimize splashing.

Display Codes

- OL** Indicates an overload condition. The load is too great for the mixer and must be reduced. Press the power button and enter in new parameters and start.
- OH** Indicates an overheat condition. The operating conditions have caused the mixer to overheat. Reduce mixer loads and/or environmental temperature to continue mixing. To reset mixer turn power off, allow mixer to cool and then turn back on.
-

Cleaning and Service

Cleaning

Unplug mixer prior to cleaning. Using a mild soap and water solution, wipe the housing and control panel of the mixer with a damp cloth. Do not use chlorine bleach, chlorine-based cleanser, abrasives, ammonia, steel wool, or scouring pads with metal content or similar harsh solvents or abrasives. These may damage the surface of the mixer

Service

Do not attempt to service or repair mixer. If the mixer housing is opened, the warranty becomes void. Contact Supplier for return authorization and return instructions. Clean the mixer so that it is free from any materials which may be harmful. Provide an SDS or other safety data sheets for all chemicals used in conjunction with this mixer where appropriate. Place the mixer and its parts into the original packing to protect and prevent damage during transport. Follow shipping directions provided when return is authorized and RMA issued.

CAUTION:

 Failure to clean, service, and handle the mixer as outlined above can cause it to fail which could impair the safety of the user.

Warranty and Product Liability

Warranty

Cole-Parmer warrants the herein described product of its manufacture for three years from receipt of the equipment by the purchaser, against defects in material and workmanship.

This limited warranty covers parts and labor. If the mixer housing is opened, the warranty becomes void.

This warranty may only be altered by a specifically published amendment. No individual has authorization to alter the provisions of this warranty policy or its amendments. Expendable items are not covered by this warranty. Damages due to corrosion or accidental breakage are not covered.

To register the warranty online, go to the Cole-Parmer website address below:

www.coleparmer.com/product-warranty-registration

Product Liability

Under no circumstances shall Cole-Parmer Instrument Company be liable for indirect, consequential, or special damages of any kind.

Cole-Parmer Instrument Company reserves the right to make technical changes without prior notice.

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