
## SETTING A TIMER

1. Press CH/CLOCK button until the desired channel is solid (not flashing) on the screen.
2. Use the H/M/S buttons to configure timer.

**Note:** To clear and start over, press CLEAR button.

## STARTING A TIMER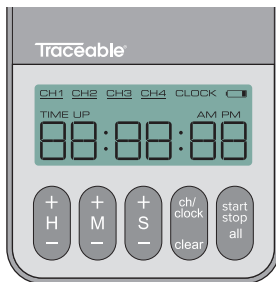

- When starting/stopping an individual timer, navigate to that timing channel using the CH/CLOCK button and then press START/STOP ALL.
- To start/stop all timers, press the START/STOP ALL button from the clock screen.
- Timers which are counting down will flash quickly.
- Timers which are counting up will flash slowly.
- Timers which have expired will flash rapidly.

## WARRANTY, SERVICE, OR RECALIBRATION

**Traceable**<sup>®</sup>  
an antylia scientific company

12554 Old Galveston Rd. Suite B230  
Webster, Texas 77598 USA  
Ph. 281-482-1714 • Fax 281-482-9448  
support@traceable.com  
traceable.com


Traceable<sup>®</sup> Products are ISO 9001:2015 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA.

Item No. 56000-16

Model No. 5667

©2022 1065T9\_M\_92-5667-00 Rev. 1 07172025

# Traceable<sup>®</sup> 4-Channel Timer Quick-Start Guide

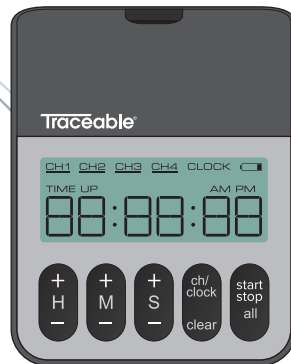## Traceable® 4-Channel Timer

Start/Stop All button


Channel/Clock Clear timer setting button

Seconds button

Minutes button


Hours button

Battery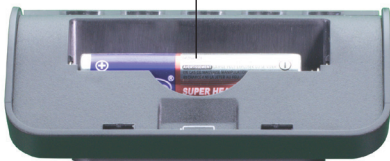

## INITIAL SETUP

- 1 Remove battery cover and insert batteries.

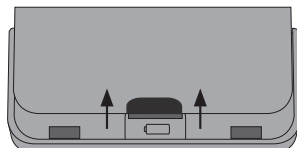
- 2 Replace battery cover.


- 3 Press CH/CLOCK button until clock appears on top row of display.


- 4 Hold CH/CLOCK until there is an audible beep and the time flashes.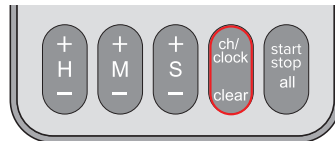

- 5 To adjust 12/24 hour clock, press and hold START/STOP ALL button.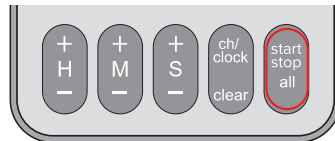


- 6 Use H/M/S buttons to adjust the time.

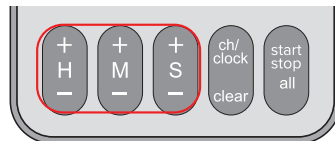- 7 Pressing Seconds button will round to the nearest minute.

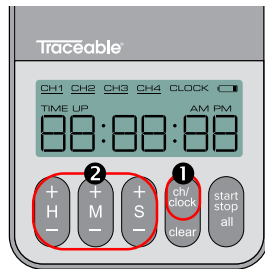- 8 Press the CH/CLOCK button to exit setup.

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