

SPECIFICATIONS

Accuracy: $\pm 1^{\circ}\text{C}$ / $\pm 2^{\circ}\text{F}$ between -4 and 248°F
(-20 and 120°C otherwise $\pm 2^{\circ}\text{C}$ / $\pm 4^{\circ}\text{F}$)

Resolution: 0.1°

Sampling: 1 second

Battery life: 5000 hours

Automatic

power off: 10 mins

OPERATION

1. Press [ON/OFF] to switch on.
2. Press [HOLD] to retain reading.
3. Press [C/F] to choose temperature unit.
4. Press [ON/OFF] again to power off.

ALL OPERATIONAL DIFFICULTIES

If this timer does not function properly for any reason, replace the battery with a new high quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

An incorrect display, no display or operational difficulties indicate that the battery should be replaced. Slide open the battery cover on the back of the thermometer. Insert the new battery with the positive side facing you. Replace the battery cover.

WARRANTY, SERVICE, OR CALIBRATION

For warranty, service, or calibration contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com

www.traceable.com

Traceable® Products are ISO 9001:2015 Quality-
Certified by DNV and ISO/IEC 17025:2017
accredited as a Calibration Laboratory by A2LA.

©2020 Traceable® Products. 92-7874-00 Rev. 1 072320

**TRACEABLE®
DIGITAL STEM
THERMOMETER
INSTRUCTIONS**