

OPERATION

Pull the Long-Stem Thermometer from the plastic sleeve. Slide the F/C switch to F to display the temperature in Celsius. Insert the stainless steel stem into the material to be measured. Slide the ON/OFF switch to OFF to turn the unit off. To conserve battery life, always turn the unit off when not in use.

BATTERY REPLACEMENT

A faint display, no display, erratic readings, or operational difficulties indicate that the battery should be replaced. Remove the battery cap, located at the top of the unit and marked "OPEN", by turning it in the direction of the arrow (counter-clockwise).

Replace the battery with a new, 1.5V, G-13 battery. Equivalent battery replacements are: RAYOVAC RW42, DURACELL D357, and EVEREADY 357. Insert the battery with the positive side facing you. Replace the battery cap.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230

Webster, Texas 77598 USA

Ph. 281 482-1714 • Fax 281 482-9448

E-mail support@traceable.com • www.traceable.com

Traceable® Products are ISO 9001:2015 Quality-Certified by DNV and ISO/IEC 17025:2017 accredited as a Calibration Laboratory by A2LA.

Traceable® is a registered trademark of Cole-Parmer.

©2020 Traceable® Products. 92-8922-00 Rev. 1 042420

**LONG-STEM
THERMOMETER
INSTRUCTIONS**