

SPECIFICATIONS

Timing Capacity: 19 hours 59 minutes
Timing Increments: 1 minute
Timing Accuracy: 0.01%

COUNTDOWN TIMING

The first press of the Hr (hour) or Min (minute) button turns the timer on (0:00 flashing).

Set the countdown time by pressing Hr (hour) or Min (minute) buttons until the desired time is on the display. (Press and hold the Hr or Min button to advance the display rapidly.)

Four (4) seconds after setting the countdown time, the timer will automatically begin counting down, this is indicated by a flashing colon (:).

TIME-OUT

Press the Hr and Min buttons simultaneously to stop the timer. Press the Min button to resume timing.

CLEARING THE DISPLAY

If the timer is running (indicated by a flashing colon), press Hr and Min buttons simultaneously to stop the timer. Press the Hr button to clear the display and turn the timer off.

ALARM

The alarm will automatically silence after 20 seconds and the timer will turn off to preserve the battery. The alarm may be silenced and the timer turned off manually by pressing the Hr button.

REMINDER ALARM

While the alarm is sounding, pressing the Min button once will silence the alarm and start the timer counting up. Once the timer reaches 10 minutes, a reminder alarm will sound.

MEMORY RECALL

While the alarm is sounding, pressing the Min button twice quickly will activate the memory feature. The display will return to the previously programmed time and begin counting down.

ALL OPERATIONAL DIFFICULTIES

If this timer does not function properly for any reason, replace the battery with a new high quality battery (see the “Battery Replacement” section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

An incorrect display, faint display or operational difficulties indicate that the battery should be replaced. To replace the battery, remove the four screws on the back of the unit. Remove the back. Equivalent battery replacements are: RAYOVAC 392 and TOSHIBA LR41. Insert the new battery with the positive side facing you. Replace the back of the unit. Replace the four screws and tighten securely.

TRACEABLE[®]
TOUCH TIMER
INSTRUCTIONS