Display:	%"-High LCD showing	Make sure tape has been removed from	pressing any button.	If this timer does not function properly for
, ,	minutes, and seconds	BATTERY COMPARTMENT. (See Battery	5. When alarm is turned off manually, press	any reason, please replace battery with
Accuracy:	0.01%	Replacement below). 00M 00S will appear	START/STOP to return to last programmed	a new, high-quality battery (see "Battery
Timing Channels:	one	on display.	setting.	Replacement" section). Low battery power can occasionally cause any number of
Timing Capacity:	99 minutes, 59 seconds	COUNTDOWN ALARM TIMING 1. Set MIN and SEC to desired time.	6. To return display to 00M 00S, press MIN	"apparent" operational difficulties. Replacing
Alarm:	Audible and visual	Note: Press and hold MIN or SEC to	and SEC simultaneously. Timer may be cleared to zero either when the count is	battery with a new, fresh battery will solve

Memory: Automatic return to last Function: Countdown timing

SPECIFICATIONS

Size:

Weight:

programmed time

3½" x 1¾ " x ½"

1¼ ounces

3. When time is up, alarm will sound for 30 seconds, turn off automatically, and return

to last programmed setting.

2. Press START/STOP button to begin timing.

BEFORE USING TIMER

advance setting rapidly.

TIME-OUT Timer may be stopped at any time by

STOP button again to resume timing.

running or when it is stopped.

4. Alarm may be turned off manually by

pressing START/STOP button. Press START/

most difficulties.

ALL OPERATIONAL DIFFICULTIES

BATTERY REPLACEMENT

An incorrect display, no display or operational difficulties indicate that battery should be replaced. Use a coin to open battery cover on back of timer (turn cover approximately 1/8 of a turn counter-clockwise). Insert battery with positive (+) side facing up. Replace battery cover. Catalog No. 1039.

