



DESCRIPTION

"HR" button: increase Hours value

"MIN" button: increase Minutes value

"SEC" button: increase Seconds value

"CLEAR" button: reset a timer setting

"CLOCK/TIMER" button: switch between clock and timer

"START/STOP" button: Start, Pause or Stop count up or count down timer. Memory recall last setting of TIMER 1
TIMER 2
TIMER 3.

Timing Channels

"TIMER 1" button: start or stop channel 1.

"TIMER 2" button: start or stop channel 2.

"TIMER 3" button: start or stop channel 3.

Press corresponding channel for channel memory recall.

SPECIFICATIONS

Clock: Displays Hr/Min/Sec (AM/PM)

Count Up Timer: 99 hours, 59 minutes and 59 seconds

Count Down Timer: 99 hours, 59 minutes and 59 seconds. Timer will beep for 1 minute once set time is reached.

Batteries: 2 x 1.5V lithium batteries

Material: ABS plastic

Dimensions: 8.82 x 6.62 x 1.32 cm, Weight: 90g.

OPERATION

To switch between "Time/Count Up" and "Count Down" mode, press the "CLOCK/TIMER" button.

1. In Time/Count Up mode, press the "CLOCK/TIMER" button for 2 seconds until screen flashes. This indicates time setting menu has been entered.
2. Press the "HR" button to increase HOURS; press "MIN" button to increase Minutes and press "SEC" button to increase Seconds. Continue pressing button to increase value.
3. Press "CLOCK/TIMER" button to confirm new setting.

COUNT UP TIMER

1. In Time/Count Up mode, press START/STOP button. Timer will begin counting up. Press START/STOP button again to pause and press again to continue counting up.
2. To clear count up time, press START/STOP to stop timer and then press "CLEAR" button.

COUNTDOWN TIMER

Press "CLOCK/TIMER" button to enter "Count Down" mode.

COUNTDOWN SETTINGS

Use three independent channels to time three different events; Channels: TIMER 1, TIMER 2, and TIMER 3.

1. Press "TIMER 1" button until TIMER 1 appears on display.
2. Press "HR" button to increase Hours value, press "MIN" button to increase Minutes value and press "SEC" button to increase Seconds. Continue pressing button to increase value.
3. Press "CLEAR" to reset the value to zero.
4. Press the TIMER 1 button to confirm new setting.
5. If multi-channel timing is desired, repeat steps above to set channels TIMER 2 and TIMER 3.

TO START A COUNTDOWN

1. Press the TIMER 1 button to start the channel 1 count down, the TIMER 2 button to start the channel 2 count down and the TIMER 3 button to start the channel 3 countdown.
2. Press the TIMER 1, TIMER 2 or TIMER 3 button again to pause timer. Press again to continue countdown.

NOTE: TIMER1 or TIMER2 or TIMER3 message will appear in the Time/Count Up mode while it is counting down.

3. Press the CLEAR button to reset a timer setting or to reset all 3 timer settings back to zero when the timer is in the pause status.
4. Press "START/STOP" button to recall or bring back last setting. When the timer is reset to zero, a "MEMORY" message will appear on the screen.

NOTE: Press "TIMER 1", "TIMER 2" or "TIMER 3" to recall or set the TIMER1, TIMER 2 or TIMER 3 timer back to zero respectively. At the same time, a "MEMORY" message will appear on its corresponding screen.

TO STOP A COUNTDOWN

1. Once it reaches set time, a "TIME'S UP" message will appear. The screen will flash, continue to count up and will beep for 1 minute.
2. When it reaches set time, press "TIMER 1" or "TIMER 2" or "TIMER 3" or "START/STOP" to stop the beep sound and the timer. Press "CLEAR" to reset it to zero.
3. You can set the timer only if the timer's value is zero.

NOTE: "TIME'S UP" and "TIMER 1" (TIMER 2 or TIMER 3) message will flash on the screen when you are in the "CLOCK/TIMER" mode.

ALL OPERATIONAL DIFFICULTIES

If this timer does not function properly for any reason, please replace battery with a new, high-quality battery (see Battery Replacement section). Low battery power can occasionally cause any number of “apparent” operational difficulties. Replacing battery with a fresh new battery will solve most difficulties.

BATTERY REPLACEMENT

An incorrect display, no display or operational difficulties indicate that the battery should be replaced. Slide open the battery cover on the back of the timer. Insert a new AAA alkaline battery. Make certain to note the proper polarity as indicated by the diagram in the battery compartment. Replace the battery cover.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

TRACEABLE® PRODUCTS

12554 Old Galveston Rd. Suite B230
Webster, Texas 77598 USA
Ph. 281 482-1714 • Fax 281 482-9448
E-mail support@traceable.com
www.traceable.com

Traceable® Products are ISO 9001:2015 Quality-
Certified by DNV and ISO/IEC 17025:2017
accredited as a Calibration Laboratory by A2LA.

Cat. No. 5008

Traceable® is a registered trademark of Cole-Parmer.

©2020 Traceable® Products. 92-5008-00 Rev 5 092524

TRACEABLE® THREE-LINE ALARM TIMER INSTRUCTIONS