

MOUNTING INSTRUCTIONS

Assembled by: **MASTERFLEX**[®]

77200-02


Mounting Instructions for two MASTERFLEX[®] L/S[®] EASY-LOAD[®] and L/S[®] EASY-LOAD[®] II PUMP HEADS.

Use only MASTERFLEX precision tubing with MASTERFLEX Pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

<u>Mounting Kit Catalog No.</u>	<u>Number of Pump Heads</u>	<u>Description</u>
77200-02	2	Stainless steel mounting hardware

Assembly Instructions:

1. Install studs into pump drive (short threaded end into drive).
2. Slide the pumps into place on the studs.
3. Install washers and wing nuts onto studs and hand-tighten.

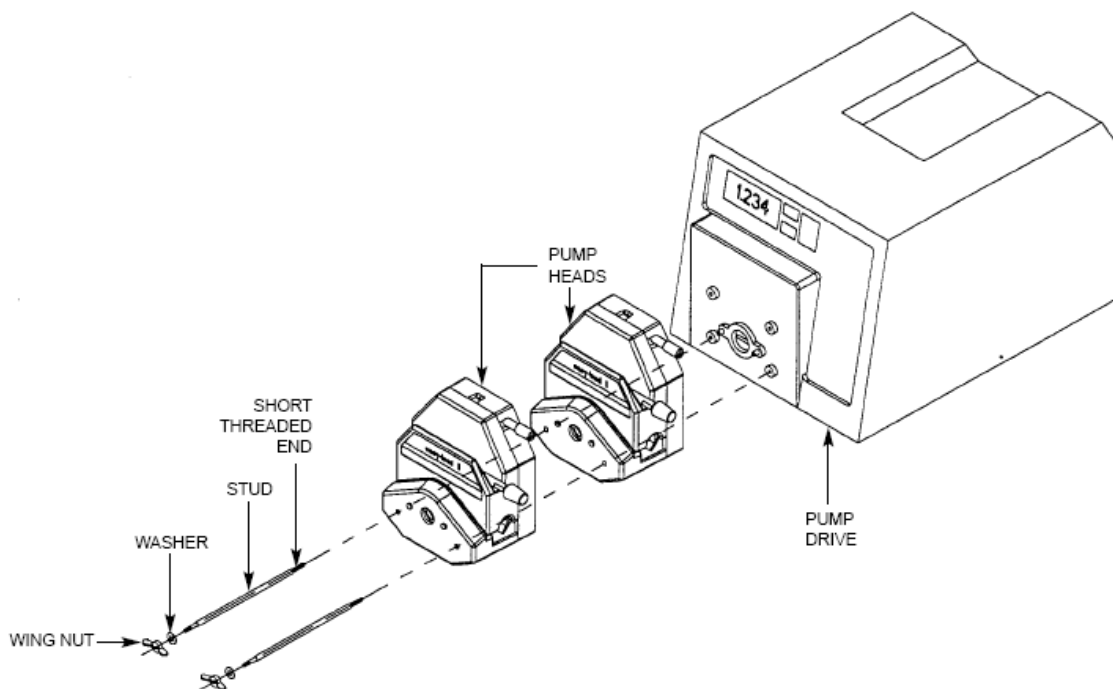

Trademarks bearing the ® symbol in this publication are registered in the U.S. and in other countries.