7523-40

OPERATING MANUAL:  
**PUMP DRIVES**

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**WARNING:** Tubing breakage may result in fluid being sprayed from pump. Use appropriate measures to protect operator and equipment.

**WARNING:** Turn drive off before removing or installing tubing. Fingers or loose clothing could get caught in drive mechanism.

### **WARNING: PRODUCT USE LIMITATION**

This product is not designed for, nor intended for use in patient connected applications; including, but not limited to, medical and dental use, and accordingly has not been submitted for FDA approval.

## Introduction

The Console Drive controls the speed of MASTERFLEX® L/S® (Laboratory/Standard) pumps to provide flow rates from 0.10 to 3400 mL/min.

Mount up to 2 (600 rpm) or 4 (100 rpm) MASTERFLEX® L/S® pump heads and all MASTERFLEX-compatible pump heads.

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**D) FLOW CONTROL**—Set flow rate for selected tubing size. To change flow rate, press ▼ or ▲ arrows. (If pump is running, its speed will change with new settings.)

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**E) CAL CONTROL**—Refine built-in calibration, using a measured volume.

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**G) PRIME**—Run pump at full speed to fill or clear lines.

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**H) DIRECTION**—To change motor direction.

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**J) POWER SWITCH**—All settings are retained in memory.

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**K) MODE SELECT**—INT for internal control; mA for remote current control; V for remote voltage control.

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## Setup and Drive Operation

1. Mount pump head and load tubing. (See pump head manual.)
2. Turn pump on and select TUBING SIZE.

**NOTE:** If CAL LED is lit, that tubing size has been previously field calibrated. If LED is not lit, the drive is operating with the built-in factory calibration. To clear a field calibration, press and hold the CAL switch until the CAL light goes out. This will take about 3 seconds. To recalibrate for better accuracy, see Calibration section.

3. MODE selection (INT, mA, V).
4. Select MOTOR DIRECTION (CW or CCW).
5. PRIME and CALibrate the pump (if required).
6. Press FLOW button and watch display to set the flow rate with UP/DOWN keys.
7. Press STOP/START button to begin pumping.

**NOTE:** Pump will restart automatically after a brownout or powerout condition.

## Calibration

*Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.*

1. Select correct tubing size and flow rate.
2. Press CAL, calibration volume appears.
3. Press STOP/START, the pump will use its stored memory to dispense the specified calibration sample quantity. The pump will stop automatically.
4. Weigh/measure the sample.
5. Use UP/DOWN arrow keys to correct the flashing display.

**NOTE:** If the adjusted calibration is too great, "Err" will appear in the display. If this occurs, press the CAL control and repeat the calibration procedure. The microprocessor will retain one special calibration value per tubing size, even when power is turned off. The next calibration will replace the existing value.

6. Press SIZE to exit the calibration cycle.

## Maximum Flow Rate ("OTH" Tubing)

1. To set the maximum flow rate for non-standard pumps or tubing sizes "OTH", press CAL, then FLOW. The maximum flow rate will then flash on the display.
2. Use UP/DOWN arrow keys to set desired flow rate.
3. Press SIZE to exit.

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## SETUP AND DRIVE OPERATION CALIBRATION

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## DISPense/copy

A first press of the DISP key results in the last entered dispense volume being displayed. The "mL" annunciator will illuminate and flash. The INC/DEC keys are used to change the dispense volume, if desired. The STOP/START key then initiates delivery of the set volume. The amount remaining to be dispensed will be displayed during countdown. The dispense function is exited by pressing any key except Increment, Decrement, DISP, or STOP/START.

A second press of the DISP key causes the COPY annunciator to illuminate and flash. The STOP/START key is then used to deliver the desired volume without the need to know the volume in specific units. A third press of the DISP key enters the volume dispensed. The COPY annunciator stops flashing. The STOP/START key is then used to initiate delivery of the copied volume. The number of copies dispensed will be displayed after each dispense. The maximum number of copies is 9999. The STOP/START key is used to pause the copy dispense during dispensing; copy dispense can then be continued using the STOP/START key.

A fourth press of the DISP key results in the last entered dispense interval being displayed. The SEC annunciator will illuminate and flash. The INC/DEC keys are used to change the dispense time, if desired, from 1 to 9999 seconds. The STOP/START key then initiates delivery for the set time interval. The remaining time will be displayed during countdown. Pressing the DISP key a fifth time exits this mode.

## Keypad Lockout Enable/Disable

Press and hold FLOW. After five (5) seconds, display will change to all dashes. Then, while holding FLOW, press PRIME five (5) times.

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## REMOTE CONTROL

Selectable input (0–20 mA, 4–20 mA, 0–10V DC )  
 ±0.5% linearity control  
 2300V isolation potential  
 START/STOP; CW/CCW; PRIME via contact closure

### Remote Control Setup

1. Place the power switch in the off position.

**CAUTION:** Power must be turned off before connecting the external remote control cable to prevent damage to the drive.

- Connect the cable from the external remote control to the DB15-pin receptacle on the rear panel.
- Select type of remote control input and output required as follows:
  - Press and hold the MODE control while turning the power switch to the on position. After two seconds, release the MODE control. The initial display will show: "inP". After two seconds the display will show either 0–20 or 4–20.

**NOTE:** Press the up (increment) or down (decrement) arrows to select between 4–20 and 0–20 for current loop control.

- Press the MODE control again. The initial display will show: "out". After two seconds the display will show either 0–20 or 4–20.

**NOTE:** Press the up (increment) or down (decrement) arrows to select between 4–20 and 0–20 for current loop output.

- Press the MODE control to select mode of operation. The LED's indicate the selected mode. Select either mA or V.

**NOTE:** If only remote STOP/START, PRIME and/or CW/CCW is to be used, the MODE control can be set to any of the three positions.

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## TROUBLESHOOTING AND MAINTENANCE

### Fuse Replacement

1. Place the power switch in the off position.
2. Disconnect the AC power input line cord from the AC receptacle.
3. Remove and check the fuse and replace if defective.
4. Reconnect the AC power input line cord to the AC receptacle.

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### A) 6-600 RPM GEAR ASSEMBLY (included in service kit 07553-06)

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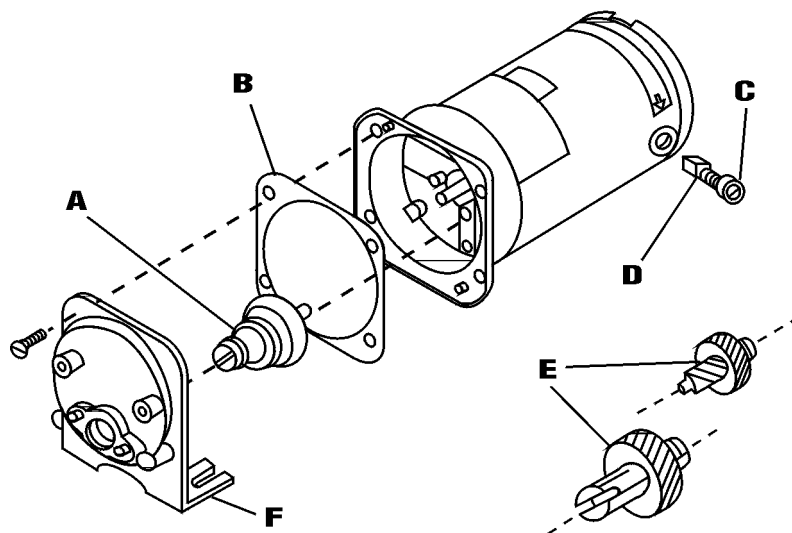
## MOTOR BRUSH CHECK/REPLACEMENT

**NOTE:** Brushes should be checked every 6 months or 2000 operating hours or if erratic operation occurs.

**DANGER:** High voltages exist and are accessible in the Console Drive. Use extreme caution when servicing internal components.

1. Place the POWER switch in the off position.
2. Disconnect the AC power input line cord from the AC receptacle.
3. Remove the three screws from each side of the housing and lift off the housing.
4. Carefully unscrew each brush cap on opposite sides of the motor. Withdraw the brush, and examine it for wear.

**NOTE:** One brush is accessible through a hole in the PC board.

**NOTE:** Replace both brushes, if either brush is less than 7.6 mm (0.300 in) long from base to point.

5. Insert brushes and install brush cap on each side of motor.
6. Install housing and secure with the three screws on each side.
7. Connect the AC power input line cord to the AC receptacle.

Contact your dealer if you have service needs.

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## Troubleshooting (cont.)

If an **error message** is displayed, refer to the following list for possible corrective action you can take. If these do not correct the problem, contact your dealer.

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## Cleaning

Keep the drive enclosure clean with mild detergents. Never immerse nor use excessive fluid.

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**REPLACEMENT PARTS  
AND ACCESSORIES;  
CLEANING;  
SPECIFICATIONS**

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<b>SPECIFICATIONS</b>	
<b>Output:</b>	
Speed:	
07523-40, -47	10 to 600 r/min
07523-50, -57	1.6 to 100 r/min
Torque output, Maximum:	
07523-40, -47	180 oz-in (13 kg•cm)
07523-50, -57	360 oz-in (26 kg•cm)
Speed regulation:	
Line	±0.25 % F.S.
Load	±0.25 % F.S.
Drift	±0.25 % F.S.
Display:	Four-digit, seven segment LED
Remote outputs:	Voltage speed output (0-10V DC) Current speed output (0-20 mA or 4-20 mA) Tach output (TTL, 128 to 7680 Hz) Motor running output (N.O. & N.C. contact closure)
<b>Input:</b>	
Supply voltage limits:	
07523-40, -50	90 to 130 Vrms @ 50/60 Hz
07523-47, -57	200 to 260 Vrms @ 50 Hz
Current, max.:	
07523-40, -50	2.3A
07523-47, -57	1.2A
Remote Inputs:	START/STOP, CW/CCW, PRIME (Contact closure) Voltage input (0-10V DC) Current input (0-20 mA or 4-20 mA)
<b>Construction:</b>	
Dimensions (L × W × H):	11.5 in × 7 in × 7 in (292 × 178 × 178 mm)
Weight:	15 lb (6.8 kg)
Enclosure Rating:	IP 23 per IEC 529
<b>Environment:</b>	
Temperature, Operating:	0° to 40°C (32° to 104°F)
Temperature, Storage:	-45° to 65°C (-49° to 149°F)
Humidity (non-condensing):	10% to 90%
Altitude:	Less than 2000 m
Pollution Degree:	Pollution Degree 2 (Indoor Usage — lab, office)
Chemical Resistance:	Exposed material is painted aluminum, plastic and vinyl
Compliance:	115V: UL508C, CSA C22.2, No 14 230V (For CE Mark): EN61010-1 (EU Low Voltage Directive) and EN61326 (EU EMC Directive)

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## WARRANTY

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is noted on the Warranty Card.

## PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

## TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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Vernon Hills, Illinois U.S.A. 60061-1844
Barrington, Illinois U.S.A. 60010-2392