



# Operation Manual

## Instruction Notes for 3622561, 3622562, 3622563, 3622564

Digital Control Module

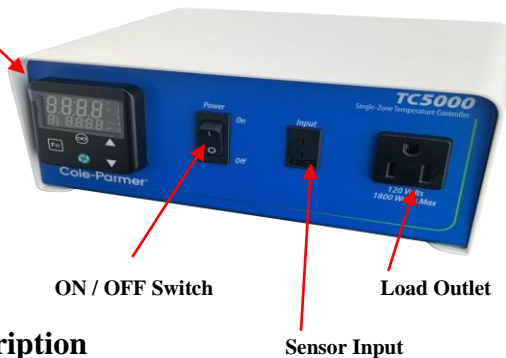

Load Fuse

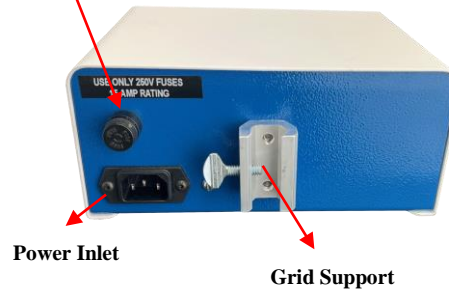


### General Description

This temperature control displays both the setpoint and process temperature and uses the most modern control technology for the best temperature stability. The Auto-tune feature minimizes setpoint overshoot and learns your process. Changing your setpoint is easily done with the interface keys. The controller can also perform ramp rate operations to allow the user to slowly raise the process temperature. There are several input types available. This unit comes with a 6' detachable power cord, grid support bracket, which is ideal for fume hood mounting to maximize bench space. This control is a microprocessor-based, digital indicating, automatic temperature control with a single input and a single output. It features an auto-tuning function that allows automatic setting of control parameters with a minimum of user input required. This family of control accepts a type "J", "K", "T" thermocouple, or RTD (model 3622564) input depending on the model ordered.

The control automatically sets the PID parameters through a "learning" sequence in the auto-tuning mode. PID parameters include proportional band, reset/integral and rate/derivative. User-friendly features include automatic LED indicators to aid in monitoring and setup, as well as dual LED displays for process temperature and set point indication. This control automatically stores all information in a non-volatile memory.

### Control Features

1. On/Off switch.
2. 15-amp output.
3. Control **PID with Auto-tune algorithm, solid state relay**
4. Circuit protection (fuse)
5. Low profile housing.
6. Safety feature- Output is de-energized if sensing device is unplugged or damaged.

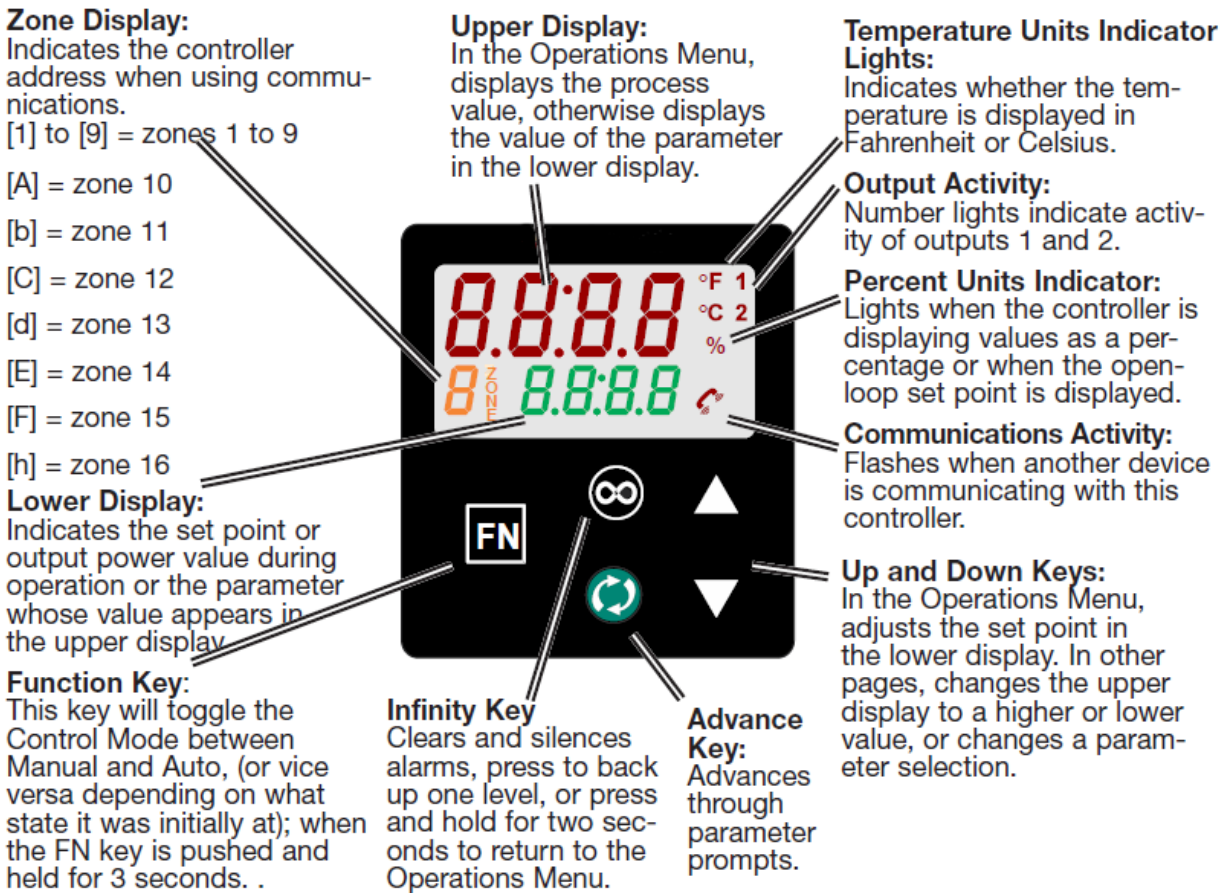

## How to Setup and Operate

First, connect the heater load and the sensor to the load and sensor receptacles on the front panel. Next, plug the control line cord into an appropriate 3-wire grounded power receptacle. Push the power switch to the "ON" position. Wait five seconds for the control unit to energize.

The control is shipped from the factory with the display reading in degrees C. If a display in degrees F is desirable, press the **UP/DOWN** arrow keys simultaneously for three seconds to access the Setup Page. Press the **Advance** key until the Celsius\_Fahrenheit parameter [C-F] is shown in the lower display. Press the **UP** arrow key to change from C to F.

After changing the display, press the **Infinity** key to display the set point temperature and the process temperature (temperature at the sensor).

To begin the heating process, use the **Up/Down** arrow keys to enter the desired set point temperature. This is shown in the lower display. Auto-tuning is recommended in all applications. Auto-tuning allows the control to set the PID parameters without those parameters being manually entered/calculated by the user. It also minimizes the amount of temperature overshoot of the set point and decreases the time required for the process to stabilize. See page 8 for instruction on performing the auto-tune function.

It is recommended that the auto-tune feature be used every time the process temperature set point has been changed. If the process is being repeated, the parameters learned in the initial setup are stored in the memory of the control and the auto-tune process isn't required to be repeated in most applications.

**Control Sensor**

Proper placement of the sensor can eliminate many problems associated in the total operation of the system. The probe should be placed so that it can detect any temperature change with little thermal lag. In a process that requires fairly constant heat output, the probe should be close to the heating source. In processes where heat demand is variable, the probe should be placed close to the work area. Experimenting with probe location should be tried to provide optimum results for your particular process.

**Specifications**

**Control Mode**

- Microprocessor-based, single input, single output.
- PID parameters
  - Proportional band: 1 to 999° F
  - Reset: 0.00 to 99.99 repeats per minute.
  - Integral: 0.00 to 99.99 minutes per repeat.
  - Rate or Derivative: 0.00 to 9.99 minutes.

**Operator Interface**

- Advance, Infinity, Up and Down keys, and ON/OFF switch.
- (2) 4 character LED display.
- Thermocouple receptacle and 3 wire load receptacle.

**Input**

- Type J, K , T thermocouple or (RTD): input grounded or ungrounded.
- Automatic cold junction compensation and break protection for sensor.
- Degrees F or degrees C display; user selectable (preset for degrees C)

<b>Range Type J:</b>	<b>-200°C to 1200°C</b>
<b>Range Type K:</b>	<b>-200°C to 1370°C</b>
<b>Range Type T:</b>	<b>-200°C to 400°C</b>
<b>Range Type RTD:</b>	<b>-200°C to 800°C</b>

**Primary Output (Heating or Cooling)**

- 15A @ 120VAC or 10A @ 240VAC depending on model

**Accuracy**

- Calibration accuracy: 0.1% of span.
- Temperature stability: 0.2 °F / °F rise in ambient maximum.
- Voltage stability: 0.01% of span / % of rated line voltage.

**Power**




- 50/60 Hz 5%
- Data retention upon power failure via nonvolatile memory.

**Operating Environment**


- 32 to 149 °F / 0 to 65 °C; 0 to 90% RH, non-condensing.

Operations Menu

Operations Menu

To enter the Operations Menu, press the Infinity Key  to return to the Home Page. Press the green Advance Key  to scroll through the various prompts found in the Operations Menu. Press the Infinity Key  at any point within the Operations Menu to return to the Home Page. Use the Up and Down Arrow Keys to increment or decrement settings or change selection.


Operations Menu

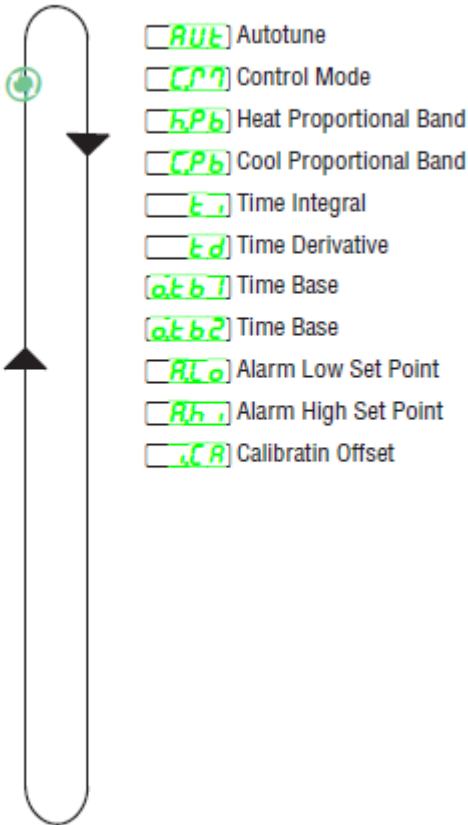

### OPERATION MENU

Upon power up of the control, using the advance key will scroll through the various prompts found in the Operations Menu. At any point within the Operations menu to return to the default display push the Infinity ^ key.

### Operations Menu

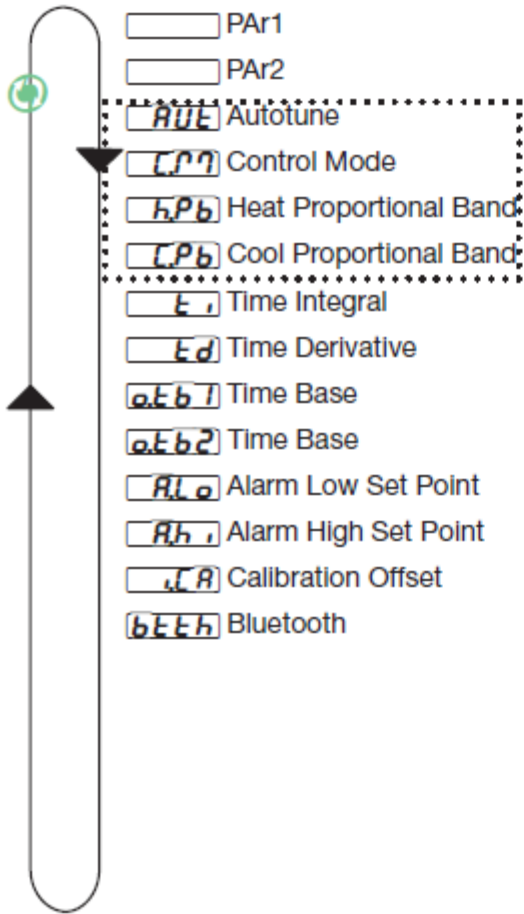


Display	Parameter Name Description
<input checked="" type="checkbox"/> <b>AUE</b> [AUT]	<p><b>Autotune</b> Start an autotune. While active the upper or left and lower or right display will flash [tUn1] and [Attn]. Appears if: Heat or cool algorithm set to PID Range: Defaults are shown below in bold <input type="checkbox"/> <b>no</b> No <input checked="" type="checkbox"/> <b>YES</b> Yes</p>
<input checked="" type="checkbox"/> <b>CPn</b> [C.M]	<p><b>Control Mode Active</b> View the current control mode. Appears if: Always Range: (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off <input checked="" type="checkbox"/> <b>AUEo</b> Auto <input checked="" type="checkbox"/> <b>CPn</b> Manual</p>
<input checked="" type="checkbox"/> <b>hPb</b> [h.Pb]	<p><b>Heat Proportional Band</b> Set the PID proportional band for the heat outputs. Appears if: Heat algorithm set to PID Range: (Defaults are shown in bold) 0 to 9,999.000°F or units 0 to 5,555.000°C <b>Units, 25.0°F or 14.0°C</b></p>
<input checked="" type="checkbox"/> <b>CPb</b> [C.Pb]	<p><b>Cool Proportional Band</b> Set the PID proportional band for the cool outputs. Appears if: Cool algorithm set to PID- Range (Defaults are shown in bold) 0 to 9,999.000°F or units 0 to 5,555.000°C <b>Units, 25.0°F or 14.0°C</b></p>

**OPERATION MENU (Continued)**

Upon power up of the control, using the advance key will scroll through the various prompts found in the Operations Menu. At any point within the Operations menu to return to the default display push the Infinity ^ key.

**Operations Menu**

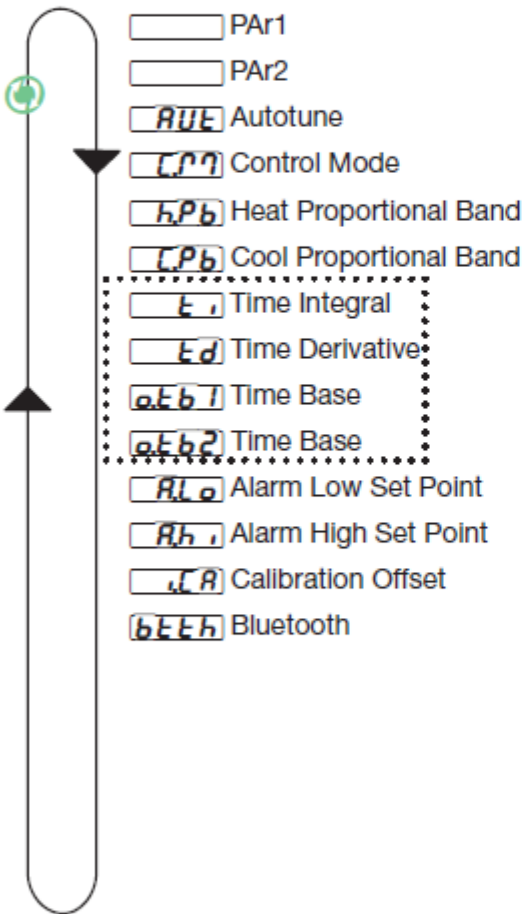


<b>ti</b> [ ti]	<b>Time Integral</b> Set the PID integral for the outputs. Appears if: Heat or cool algorithm set to PID Range (Defaults are shown in bold) 0 to 9,999 seconds per repeat <b>180</b>
<b>td</b> [ td]	<b>Time Derivative</b> Set the PID derivative time for the outputs. Appears if: Heat or cool algorithm set to PID Range (Defaults are shown in bold) 0 to 9,999 seconds <b>0 seconds</b>
<b>tb1</b> [o.tb1]	<b>Time Base Output 1</b> Set the time base for fixed-time-base control. Appears if: Output 1 set to heat or cool with control algorithm set to PID. Range (Defaults are shown in bold) 0.1 to 60.0 seconds (solid-state relay or switched dc) 5.0 to 60.0 seconds (mechanical relay & NO-ARC power control) <b>1 sec. [SSR &amp; sw dc], 20.0 sec. [mech. relay &amp; NO-ARC]</b>
<b>tb2</b> [o.tb2]	<b>Time Base Output 2</b> Set the time base for fixed-time-base control. Appears if: Output 2 set to heat or cool with control algorithm set to PID. Range (Defaults are shown in bold) 0.1 to 60.0 seconds (solid-state relay or switched dc) 5.0 to 60.0 seconds (mechanical relay & NO-ARC power control) <b>1 sec. [SSR &amp; sw dc], 20.0 sec. [mech. relay &amp; NO-ARC]</b>

**OPERATION MENU (Continued)**

Upon power up of the control, using the advance key will scroll through the various prompts found in the Operations Menu. At any point within the Operations menu to return to the default display push the Infinity ^ key.


**Operations Menu**

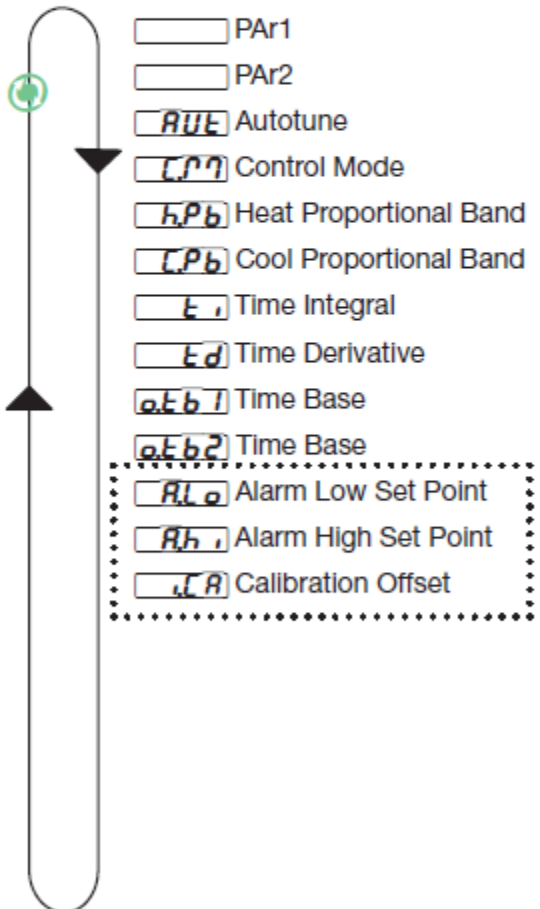

<p>[ALo] [A.Lo]</p>	<p><b>Alarm Low Set Point</b>                  Process - set the process value that will trigger a low alarm.                  Deviation - set the span of units (Must be negative number) from the closed loop set point that will trigger a low alarm.                  Appears if: If Alarm Type [A;ty] is set to Process or Deviation Alarm                  Range (Defaults are shown in bold)                  -1,999.000 to 9,999.000°F or units                  -1,128.000 to 5,537.000°C  <b>Units, 32.0°F or 0.0°C</b></p>
<p>[Ah] [A.hi]</p>	<p><b>Alarm High Set Point</b>                  Process - set the process value that will trigger a high alarm.                  Deviation - set the span of units from the closed loop set point that will trigger a high alarm.                  Appears if: If Alarm Type (A.ty) is set to Process or Deviation Alarm                  Range (Defaults are shown in bold)                  -1,999.000 to 9,999.000°F or units                  -1,128.000 to 5,537.000°C  <b>Units, 300.0°F or 150.0°C</b></p>
<p>[CA] [I.CA]</p>	<p><b>Calibration Offset</b>                  Set an offset value for a process output.                  Appears if: Always                  Range (Defaults are shown in bold)                  [-1,999.000 to 9,999.000°F or units,                  -1,110.555 to 5,555.000°C, <b>0.0</b></p>

## SET UP MENU

To enter the Setup Menu push and hold the up and down arrow keys for approximately 3 seconds. Push the green advance key to scroll through to the prompt of choice. Use the up and down arrow keys to change the range.

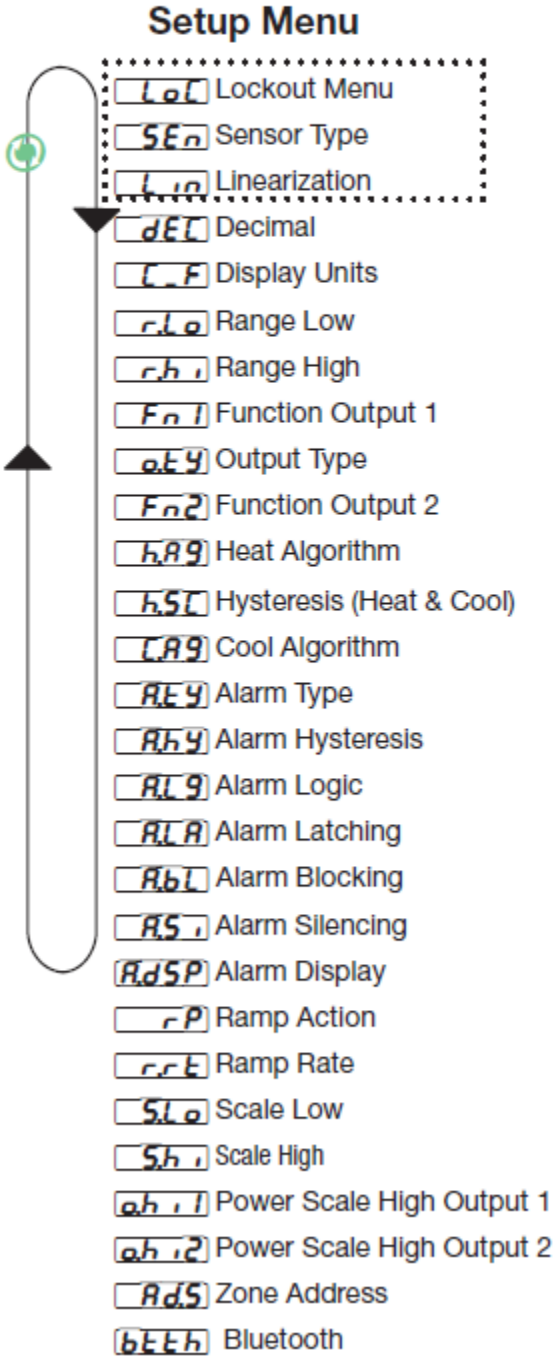

Setup Menu													
Display	Parameter Name Description												
<b>LoL</b> [LoC]	<p><b>Lockout Menu</b> Set the security clearance level. The user can access the selected level and all lower levels. Appears if: Always Range (Defaults are shown in bold) <b>1 to 5 Operations Menu and Setup Menu full access</b> You can change the security level at any level.</p> <ol style="list-style-type: none"> <li>1. Operations Menu (read only)*</li> <li>2. Operations Menu (Set point R/W)*</li> <li>3. Operations Menu (Set point R/W, Control Mode R/W)*</li> <li>4. Operations Menu R/W access)*</li> <li>5. Operations Menu and Setup Menu full R/W access</li> </ol>												
<b>SEn</b> [SEn]	<p><b>Sensor Type</b> Set the analog sensor type to match the device wired to this input. Appears if: Always Range (Defaults are shown in bold) <b>ET</b> Thermocouple, <b>uoLE</b> Volts dc, <b>PA</b> Milliamps dc <b>rQIH</b> RTD 100 Ω, <b>EhEr</b> Thermistor</p>												
<b>Lin</b> [Lin]	<p><b>Linearization</b> Set the linearization to match the thermocouple type wired to this input. For example, select [“H] for a type K Thermocouple. Appears if: Sensor Type is set to Thermocouple. Range (Defaults are shown in bold)</p> <table style="width: 100%; border: none;"> <tr> <td><b>b</b> B</td> <td><b>F</b> F</td> <td><b>r</b> R</td> </tr> <tr> <td><b>c</b> C</td> <td><b>J</b> J</td> <td><b>S</b> S</td> </tr> <tr> <td><b>d</b> D</td> <td><b>H</b> K</td> <td><b>E</b> T</td> </tr> <tr> <td><b>E</b> E</td> <td><b>n</b> N</td> <td></td> </tr> </table>	<b>b</b> B	<b>F</b> F	<b>r</b> R	<b>c</b> C	<b>J</b> J	<b>S</b> S	<b>d</b> D	<b>H</b> K	<b>E</b> T	<b>E</b> E	<b>n</b> N	
<b>b</b> B	<b>F</b> F	<b>r</b> R											
<b>c</b> C	<b>J</b> J	<b>S</b> S											
<b>d</b> D	<b>H</b> K	<b>E</b> T											
<b>E</b> E	<b>n</b> N												

### SET UP MENU (Continued)

To enter the Setup Menu, push and hold the up and down arrow keys for approximately 3 seconds. Push the green advance key to scroll through to the prompt of choice. Use the up and down arrow keys to change the range.


#### Setup Menu

- L<sub>o</sub>L** Lockout Menu
- S<sub>E</sub>n** Sensor Type
- L<sub>i</sub>n** Linearization
- d<sub>E</sub>L** Decimal
- C<sub>\_</sub>F** Display Units
- r<sub>L</sub>o** Range Low
- r<sub>h</sub>i** Range High
- F<sub>n</sub>1** Function Output 1
- o<sub>E</sub>y** Output Type
- F<sub>n</sub>2** Function Output 2
- H<sub>A</sub>g** Heat Algorithm
- h<sub>S</sub>T** Hysteresis (Heat & Cool)
- C<sub>A</sub>g** Cool Algorithm
- A<sub>E</sub>y** Alarm Type
- A<sub>H</sub>y** Alarm Hysteresis
- A<sub>L</sub>g** Alarm Logic
- A<sub>L</sub>A** Alarm Latching
- A<sub>b</sub>L** Alarm Blocking
- A<sub>S</sub>i** Alarm Silencing
- A<sub>d</sub>S<sub>P</sub>** Alarm Display
- r<sub>P</sub>** Ramp Action
- r<sub>r</sub>E** Ramp Rate
- S<sub>L</sub>o** Scale Low
- S<sub>h</sub>i** Scale High
- o<sub>h</sub>i1** Power Scale High Output 1
- o<sub>h</sub>i2** Power Scale High Output 2
- A<sub>d</sub>S** Zone Address
- b<sub>E</sub>E<sub>h</sub>** Bluetooth

Setup Menu (cont.)	
Display	Parameter Name Description
<input type="checkbox"/> <b>d<sub>E</sub>L</b> [ dEC]	<b>Decimal</b> Set the precision of the displayed value. Appears if: Always Range (Defaults are shown in bold) <input type="checkbox"/> <b>0</b> Whole, <input type="checkbox"/> <b>00</b> Tenths, <input type="checkbox"/> <b>000</b> Hundredths
<input type="checkbox"/> <b>C<sub>_</sub>F</b> [ C_F]	<b>Display Units</b> Select which units will be displayed. Appears if: Always Range (Defaults are shown in bold) <input type="checkbox"/> <b>F</b> °F <input type="checkbox"/> <b>C</b> °C
<input type="checkbox"/> <b>r<sub>L</sub>o</b> [ r.Lo]	<b>Range Low</b> For process signals, this value scales the units to minimum electrical units (0 volts or 4 mA). For TC, RTD or Thermistor, this value is the parameter for setting the minimum allowable setpoint . Appears if: Always Range (Defaults are shown in bold) -1,999.000 to 9,999.000 <b>0.0</b>
<input type="checkbox"/> <b>r<sub>h</sub>i</b> [ r.hi]	<b>Range High</b> For process signals, this value scales the units to maximum electrical units (10 volts or 20 mA). For TC, RTD or Thermistor, this value is the parameter for setting the maximum allowable setpoint . Appears if: Always Range (Defaults are shown in bold) -1,999.000 to 9,999.000
<input type="checkbox"/> <b>F<sub>n</sub>1</b> [ fn1]	<b>Function of Output 1</b> Select which function will drive this output. Appears if: If output 1 is ordered Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off, <input type="checkbox"/> <b>Cool</b> Cool, <input type="checkbox"/> <b>HEAT</b> Heat, <input type="checkbox"/> <b>ALP</b> Alarm
<input type="checkbox"/> <b>o<sub>E</sub>y</b> [ o.ty]	<b>Output Type</b> Select whether the process output will operate in volts or milliamps. Appears if: A process output (PM _ C _ F _ _ AAA _ _ ) Range (Defaults are shown in bold) <input type="checkbox"/> <b>vOL</b> Volts <input type="checkbox"/> <b>mA</b> Milliamps

### SET UP MENU (Continued)

To enter the Setup Menu push and hold the up and down arrow keys for approximately 3 seconds. Push the green advance key to scroll through to the prompt of choice. Use the up and down arrow keys to change the range.



#### Setup Menu

- LoL** Lockout Menu
- SEn** Sensor Type
- Lin** Linearization
- dEL** Decimal
- L.F** Display Units
- rLo** Range Low
- rHi** Range High
- Fn1** Function Output 1
- oEY** Output Type
- Fn2** Function Output 2
- hAg** Heat Algorithm
- hSC** Hysteresis (Heat & Cool)
- CAG** Cool Algorithm
- AEY** Alarm Type
- AhY** Alarm Hysteresis
- ALg** Alarm Logic
- ALA** Alarm Latching
- AbL** Alarm Blocking
- ASi** Alarm Silencing
- AdSP** Alarm Display
- rP** Ramp Action
- rrE** Ramp Rate
- SLo** Scale Low
- Shi** Scale High
- oHi1** Power Scale High Output 1
- oHi2** Power Scale High Output 2
- AdS** Zone Address
- bEEh** Bluetooth

Setup Menu (cont.)	
Display	Parameter Name Description
<input type="checkbox"/> <b>Fn2</b> [ fn2]	<b>Function of Output 2</b> Select which function will drive this output. Appears if: If output 2 is ordered Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off, <input type="checkbox"/> <b>CoOL</b> Cool, <input type="checkbox"/> <b>HEAT</b> Heat, <input type="checkbox"/> <b>ALPn</b> Alarm
<input type="checkbox"/> <b>hAg</b> [ h.Ag]	<b>Heat Algorithm</b> Set the heat control method. Appears if: Output 1 or 2 set to heat Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off, <input type="checkbox"/> <b>Pid</b> PID <input type="checkbox"/> <b>oNoF</b> On-Off
<input type="checkbox"/> <b>hSC</b> [ hSC]	<b>Hysteresis (Heat &amp; Cool)</b> Set the control switching hysteresis for on-off control. This determines how far into the "on" region the process value needs to move before the output turns on. Appears if: Heat or Cool Algorithm is set to On-Off. Range (Defaults are shown in bold) 0 to 9,999.000°F or units 0 to 5,555.000°C <b>Units, 3.0°F or 2.0°C</b>
<input type="checkbox"/> <b>CAG</b> [ C.Ag]	<b>Cool Algorithm</b> Set the cool control method. Appears if: If Output 1 or 2 is set to cool Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off, <input type="checkbox"/> <b>Pid</b> PID <input type="checkbox"/> <b>oNoF</b> On-Off
<input type="checkbox"/> <b>AEY</b> [ A.ty]	<b>Alarm Type</b> Selects whether the Alarm will be a Process Alarm/Absolute Temperature; or a Deviation Alarm/Temperature with respect to Setpoint. Appears if: Always Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off, <input type="checkbox"/> <b>PrAL</b> Process Alarm <input type="checkbox"/> <b>dEARL</b> Deviation Alarm

### SET UP MENU (Continued)

To enter the Setup Menu push and hold the up and down arrow keys for approximately 3 seconds. Once there, push the green advance key to scroll through to the prompt of choice and then use the up and down arrow keys to change the range.

#### Setup Menu

- L<sub>o</sub>L** Lockout Menu
- S<sub>E</sub>n** Sensor Type
- L<sub>i</sub>n** Linearization
- d<sub>E</sub>L** Decimal
- L<sub>U</sub>F** Display Units
- r<sub>L</sub>o** Range Low
- r<sub>h</sub>i** Range High
- F<sub>n</sub>1** Function Output 1
- o<sub>E</sub>Y** Output Type
- F<sub>n</sub>2** Function Output 2
- h<sub>R</sub>g** Heat Algorithm
- h<sub>S</sub>L** Hysteresis (Heat & Cool)
- C<sub>R</sub>g** Cool Algorithm
- A<sub>E</sub>Y** Alarm Type
- A<sub>H</sub>Y** Alarm Hysteresis
- A<sub>L</sub>g** Alarm Logic
- A<sub>L</sub>A** Alarm Latching
- A<sub>b</sub>L** Alarm Blocking
- A<sub>S</sub>i** Alarm Silencing
- A<sub>d</sub>S<sub>P</sub>** Alarm Display
- r<sub>P</sub>** Ramp Action
- r<sub>r</sub>E** Ramp Rate
- S<sub>L</sub>o** Scale Low
- S<sub>h</sub>i** Scale High
- o<sub>h</sub>i1** Power Scale High Output 1
- o<sub>h</sub>i2** Power Scale High Output 2
- A<sub>d</sub>S** Zone Address
- b<sub>E</sub>E<sub>H</sub>** Bluetooth

Setup Menu (cont.)	
Display	Parameter Name Description
<input type="checkbox"/> <b>A<sub>H</sub>Y</b> [ A.hy]	<p><b>Alarm Hysteresis</b> Set the hysteresis for an alarm. This determines how far into the safe region the process value needs to move before the alarm can be cleared.</p> <p>Appears if: When alarm type is set to process or deviation alarm Range (Defaults are shown in bold) 0.001 to 9,999.000°F or units 0.001 to 5,555.000°C <b>Units, 1.0°F or 1.0°C</b></p>
<input type="checkbox"/> <b>A<sub>L</sub>g</b> [A.Lg]	<p><b>Alarm Logic</b> Select what the output condition will be during the alarm state.</p> <p>Appears if: Always Range (Defaults are shown in bold)</p> <p><input type="checkbox"/> <b>A<sub>L</sub>L</b> Close on Alarm <input type="checkbox"/> <b>A<sub>L</sub>o</b> Open on alarm</p>
<input type="checkbox"/> <b>A<sub>L</sub>A</b> [ A.LA]	<p><b>Alarm Latching</b> Turn alarm latching on or off. A latched alarm has to be turned off by the user.</p> <p>Appears if: When alarm type is set to process or deviation alarm Range (Defaults are shown in bold)</p> <p><input type="checkbox"/> <b>n<sub>L</sub>A<sub>E</sub></b> Non-Latching <input type="checkbox"/> <b>L<sub>A</sub>E</b> Latching</p>
<input type="checkbox"/> <b>A<sub>b</sub>L</b> [ A.bL]	<p><b>Alarm Blocking</b> Select when an alarm will be blocked. After startup and/or after the set point changes, the alarm will be blocked until the process value enters the normal range.</p> <p>Appears if: When alarm type is set to process or deviation alarm Range (Defaults are shown in bold)</p> <p><input type="checkbox"/> <b>o<sub>F</sub>F</b> Off <input type="checkbox"/> <b>S<sub>E</sub>r</b> Startup <input type="checkbox"/> <b>S<sub>E</sub>P<sub>E</sub></b> Set Point <input type="checkbox"/> <b>b<sub>o</sub>E<sub>H</sub></b> Both</p>

### SET UP MENU (Continued)

To enter the Setup Menu push and hold the up and down arrow keys for approximately 3 seconds. Once there, push the green advance key to scroll through to the prompt of choice and then use the up and down arrow keys to change the range.

### Setup Menu


- 
- LoL** Lockout Menu
  - SEn** Sensor Type
  - Lin** Linearization
  - dEL** Decimal
  - L.F** Display Units
  - rLo** Range Low
  - rhi** Range High
  - Fn1** Function Output 1
  - oEY** Output Type
  - Fn2** Function Output 2
  - hA9** Heat Algorithm
  - hST** Hysteresis (Heat & Cool)
  - LA9** Cool Algorithm
  - AEY** Alarm Type
  - AHY** Alarm Hysteresis
  - AL9** Alarm Logic
  - ALR** Alarm Latching
  - AbL** Alarm Blocking
  - ASi** Alarm Silencing
  - AdSP** Alarm Display
  - rP** Ramp Action
  - rrE** Ramp Rate
  - SLo** Scale Low
  - Shi** Scale High
  - oHi1** Power Scale High Output 1
  - oHi2** Power Scale High Output 2
  - AdS** Zone Address
  - bEEh** Bluetooth

Setup Menu (cont.)	
Display	Parameter Name Description
<input type="checkbox"/> <b>ASi</b> [ A.Si]	<b>Alarm Silencing</b> Turn alarm silencing on to allow the user to disable the output tied (configured) to this alarm Appears if: When alarm type is set to process or deviation alarm Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off <input type="checkbox"/> <b>oN</b> On
<input type="checkbox"/> <b>AdSP</b> [ A.dSP]	<b>Alarm Display</b> Display an alarm message when an alarm is active. Appears if: When alarm type is set to process or deviation alarm Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off <input type="checkbox"/> <b>oN</b> On
<input type="checkbox"/> <b>rP</b> [ rP]	<b>Ramp Action</b> Select when the controller's set point will ramp to the defined end set point. Appears if: Always Range (Defaults are shown in bold) <input type="checkbox"/> <b>oFF</b> Off, <input type="checkbox"/> <b>SEr</b> Startup <input type="checkbox"/> <b>SEPE</b> Set Point Change, <input type="checkbox"/> <b>boEH</b> Both
<input type="checkbox"/> <b>rrE</b> [ r.rt]	<b>Ramp Rate</b> Set the rate for the set point ramp. Set the time units for the rate with the Ramp Scale parameter. Appears if: Ramp Action is set to Startup, Set Point or Both Range Range is 0.0 to 9,999 deg F, or 0.0 to 5.555 deg. C per hour. (Defaults are shown in bold) <b>1.0°F degrees or units per hr.</b>
<input type="checkbox"/> <b>SLo</b> [ S.Lo]	<b>Scale Low Output 1</b> Set the scale low for process output in electrical units. This value; in volts or milliamps, will correspond to 0% PID power output. Appears if: Output 1 is a Process set to heat or cool Range (Defaults are shown in bold) -100.0 to 100.0, <b>0.0</b>

### SET UP MENU (Continued)

To enter the Setup Menu, push and hold the up and down arrow keys for approximately 3 seconds. Once there, push the green advance key to scroll through to the prompt of choice and then use the up and down arrow keys to change the range.

### Setup Menu

- 
- [LoL] Lockout Menu
  - [SEn] Sensor Type
  - [Lin] Linearization
  - [dEL] Decimal
  - [L.F] Display Units
  - [rLo] Range Low
  - [rh,] Range High
  - [Fn1] Function Output 1
  - [oEY] Output Type
  - [Fn2] Function Output 2
  - [hA9] Heat Algorithm
  - [hST] Hysteresis (Heat & Cool)
  - [CA9] Cool Algorithm
  - [AEY] Alarm Type
  - [AHY] Alarm Hysteresis
  - [AL9] Alarm Logic
  - [ALR] Alarm Latching
  - [AbL] Alarm Blocking
  - [AS,] Alarm Silencing
  - [AdSP] Alarm Display
  - [rP] Ramp Action
  - [rRE] Ramp Rate
  - [SLo] Scale Low
  - [Sh,] Scale High
  - [oh,1] Power Scale High Output 1
  - [oh,2] Power Scale High Output 2
  - [AdS] Zone Address
  - [bEEh] Bluetooth

Setup Menu (cont.)	
Display	Parameter Name Description
[Sh,] [S.hi]	<b>Scale High Output 1</b> Set the scale high for process output in electrical units. This value; in volts or milliamps, will correspond to 100% PID power output. Appears if: Output 1 is a Process set to heat or cool Range (Defaults are shown in bold) -100.0 to 100.0, <b>10.0</b>
[oh,1] [o.hi1]	<b>Power Scale High Output 1</b> Set maximum value of output 1 range. Appears if: Output 1 is Switched and set to heat or cool Range (Defaults are shown in bold) 0.0 to 100%, <b>100.0</b>
[oh,2] [o.hi2]	<b>Power Scale High Output 2</b> Set maximum value of output 2 range. Appears if: Output 2 is Switched and set to heat or cool Range (Defaults are shown in bold) 0.0 to 100% <b>100.0</b>
[PAR1] [PAR1]	<b>Upper or Left Display</b> Select parameter to display. Appears if: Always Range (Defaults are shown in bold) <b>[ALPU] Active Process Value,</b> [Rh,] Alarm High Set Point [ALo] Alarm Low Set Point, <b>[nonE]</b> None
[PAR2] [PAR2]	<b>Lower or Right Display</b> Select parameter to display. Appears if: Always Range (Defaults are shown in bold) <b>[ALSP] Active Set Point,</b> [Rh,] Alarm High Set Point [ALo] Alarm Low Set Point, <b>[nonE]</b> None
[AdS] [Ad.S]	<b>Zone Address - Standard Bus Communication</b> Set zone address from 1-16. Appears if: Always Range (Defaults are shown in bold) 1-16, <b>1</b>
[bEEh] [btth]	<b>Bluetooth</b> Sets the power to the Bluetooth® module to on or off. Appears if: Always. Range (Defaults are shown bold) <b>On</b> , Off The Bluetooth radio can only be turned on via the front panel once turned off.

### How to Tune and Operate

Manual operation provides direct (time proportioned % time) control from -100% to 100%. A negative value is allowed only when Ot 1=Cool. Automatic operation provides sensory feedback ON/OFF or PID control. When the operation transfers from automatic to manual operation, the power level from automatic operation is retained and restored to the previous set point.

The % LED indicates manual operation. The LED is on when in Manual operation and off when in AUTO operation. When the LED is flashing, press the **Infinity** key again within 5 seconds to complete the change in operation. If the sensor is open and LOC=0, 1 or 2, the control switches to Manual operation if the output was stable before the break occurred.


### **Tuning-Automatic**

It is recommended that auto-tuning be used each time a new process is initiated. If the process is being duplicated, the parameters learned in the initial auto tune setup are stored in the memory of the control and the auto-tune procedure isn't required in most applications.

Auto-tuning: This control can automatically tune the PID parameters to fit the characteristics of your particular thermal system.

Auto-tune can only be used when Output 1 setting is hEAAt and heat algorithm (h.Ag, page 12) is set for PID mode. The heat algorithm setting is set for PID from the factory. When the auto-tune sequence has started, the lower display will flash between tun1-attn and the main operation screen.

Changing the set point, while in auto-tune mode, will not change the set point value. The displayed value will change but the control has locked in the original set point value till the auto tune is complete or has been turned off by the user. Once the control has finished "learning" the system, it returns to standard PID control with the values automatically set. Auto-tuning is complete within 80 minutes. In order for the control to successfully complete auto-tune, the process must cross the 90% set point value four times within 80 minutes after auto-tune has started. If this does not happen within the 80 minute time limit, Pb1 remains at 0 and the control will operate with the factory default PID values.

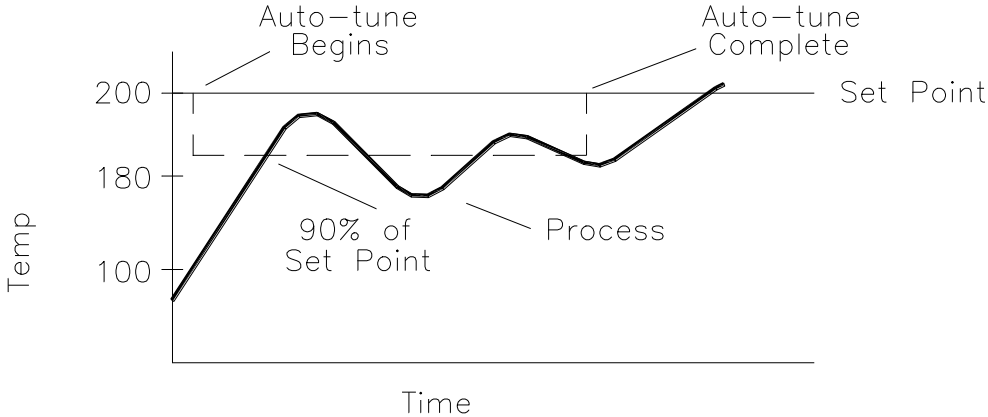

To start auto-tuning:

1. Set the set point value with the arrow keys to the desired temperature of operation.
2. Press the **Advance** key until the Aut prompt appears in the lower display.
3. Change the parameter to **YES** with the arrow keys.

4. Press the **Infinity** key to display the set point and sensor temperatures. While the control is in the auto-tuning mode, the display will alternate between the set point and the "Aut" prompt. When tuning is complete, the lower display indicates only the set point.
5. When tuning is complete, the lower display indicates setpoint only and Aut reverts to OFF. The control installs appropriate PID tuning parameters and retains them in the non-volatile memory.

\*To abort auto-tuning, the operator must reset Aut to OFF. Auto-tuning may also be aborted by cycling power off and on. In all cases, aborting auto-tuning restores all original factory default values.

### Temperature overshoot

In most cases this is due to a small sample size being heated with a high wattage heater. The control is set at the factory with the Power Scale High Output 1 (o.hil) with a value of 100. This value can be adjusted in the setup menu, reference page 13 for instructions on how to change value. If excessive temperature overshoot is occurring, adjust the value to 50 and start the heating process from ambient room temperature. If excessive temperature overshoot is still occurring, decrease to a lower value and repeat heating process from ambient room temperature. If the process temperature never makes it to the set point temperature value, increase the power scale high output 1 value. This process can take several value changes to obtain proper temperature control with your particular application setup. It is recommended to initiate the Auto-Tune process again after the proper power scale high output 1 value has been determined for your application.

**Error Code Definitions and Actions:**

Indication	Description	Possible Causes	Corrective Action
<b>EC11</b> Error Input	Sensor does not provide a valid signal to controller	Sensor improperly wired or open, Incorrect setting or sensor type, or Calibration corrupt	Correct wiring or replace sensor, Match setting to sensor, or Check calibration of controller
<b>EUN1</b> Autotuning 1	Controller is autotuning the control loop	User started the autotune function	Wait until autotune complets or disable autotune feature
<b>rP1</b> Ramping 1	Controller is ramping to new set point	Ramping feature is activated	Disable ramping feature if not required
<b>ALh1</b> Alarm High <b>ALh2</b> <b>ALh3</b> <b>ALh4</b>	Sensor input above high alarm set point	Temperature is greater than alarm set point	Check casue of over temperature, or Set alarm source to proper setting

## **GLOSSARY**

**Automatic prompts:** Data entry points where a microprocessor-based control "prompts" or asks the operator for information input.

**Auto-tune:** Automatically tunes the parameters to fit the characteristics of your particular thermal system.

**Cold junction compensation:** Electronic means to compensate for the effect temperature at the cold junction.

**Cycle time:** The time necessary to complete a full On-through-Off period in a time proportioning control system.

**Derivative/Rate:** Anticipatory action that senses the rate of change of the process, and compensates to minimize overshoot and undershoot.

**Default parameters:** The parameters (programmed instructions) permanently stored in microprocessor software to provide the data base.

**Droop:** Difference in temperature between set point and stabilized process temperature.

**Hysteresis:** In On/Off control, the temperature change necessary to change the output from On to full Off.

**Input (sensor):** Process variable information being supplied to the instrument.

**Integral/Reset:** Control action that automatically eliminates offset, or "droop", between set point and actual process temperature.

**Offset:** Adjustment to actual input temperature and to the temperature valves the control uses for display and control.

**ON/OFF control:** Control of temperature about a set point by turning the output full On below set point and full Off above set point.

**Output:** Action in response to difference between set point and process variable.

**Overshoot:** Condition where temperature exceeds stepping due to initial power up or process changes.

**Parameter:** a physical property whose value determines the response of a electronic control to given inputs.

**PID:** Proportioning control with auto-reset and rate.

**Process variable:** Thermal system element to be regulated, such as time, temperature, relative humidity, etc.

**Proportional band:** Span of temperature about the set point where time proportional control action takes place.

**Set point:** Intended value of the process variable.

**Thermal system:** A regulated environment consisting of a heat source, heat transfer medium, sensing device and a process variable control.

**Thermocouple:** Temperature sensing device that is constructed of two dissimilar metals wherein a measurable, predicated voltage is generated corresponding to temperature.

**Thermocouple break protection:** Fail-safe operation that assures output shutdown upon an open thermocouple condition.

Time Proportioning Control: Action which varies the amount of ON and OFF time when "close" to the set point (within the proportional band). This variance is proportional to the difference between the set point and the actual process temperature.

**Maintenance**

Simple preventative maintenance steps include keeping the controller clean. Protect it from overload, excessive dirt, oil and corrosion.

**Warranty**

Cole-Parmer, for itself does hereby offer a warranty for products from the date of receipt by the user, under normal and proper usage, against defects in workmanship and materials for 12 months, and will repair or replace any defective part(s) without charge when same is shipped Prepaid to Cole-Parmer from which the product was originally purchased.

Should the nature of any defect require that the product, or any constituent portion thereof, be returned to Cole-Parmer, Vernon Hill, Illinois, prepaid for service, a condition precedent to any return shall be the procurement of authorization from Cole-Parmer assigning a **Return Goods Number** to the product or part requiring service.

Parts and accessories manufactured by others are warranted only to the extent of the regular warranty of the manufacturer or supplier of such materials and only insofar as Cole-Parmer is able to transfer the benefits of warranty coverage, if any, to the user. Any adequately warranted defective part or accessory manufactured or supplied by others may be exchanged through Cole-Parmer for a replacement part is shipped prepaid and received at Cole-Parmer within 30 days from the date any replacement part is obtained by the user.

This warranty supersedes and is given in lieu of all implied warranties, and is void if the user causes damages from improper usage of product under normal operating conditions.

12 MONTHS LIMITED WARRANTY ON ALL PARTS AND LABOR IS GIVEN BY COLE-PARMER.

CATALOG NUMBER 3622561, 3622562, 3622563, or 3622564

SERIAL NUMBER \_\_\_\_\_

DATE OF PURCHASE \_\_\_\_\_

**Cole-Parmer®**  
625 East Bunker Court, MS 18  
Vernon Hills, Illinois 60061-1844, U.S.A  
Phone: 847-549-7600 Fax: 847-549-7676  
( 800) 323-4340