

Cole-Parmer®

User Manual

Cole-Parmer® TC9600 Advanced Multiparameter Temperature Controller with Thermocouple, Thermistor, and RTD Inputs

Models 89800-13 and 89800-14



THE STANDARD IN PRECISION MEASUREMENT



Table of Contents

Introduction	3
Unpacking	3
Controller Description—Front Panel	4
Controller Description—Back Panel	5
Controller Description—Active Display Screen	6
Setup and Operation	7
Initial Setup	7
Basic Operation Setup.....	7
Auto Tune Setup	7
PID Control Mode Setup (Default Mode)	8–15
ON/OFF Control Mode Setup	16–22
Ramp & Soak / Ramp Rate Control Mode Setup	23–29
Data Acquisition Software	30–33
Safety	34
Specifications for Controller Models	35–36
Specifications for Sensor Inputs	37
Screen Flow Charts	38–39
General Setup Flow Chart	38
Advanced Menu Setup Flow Chart	39
RTD Wiring Instructions	40
Ferrite Clip Installation	40
Alternative Power Cords—for Various Countries	41
Output Cord Adapters—for Various Countries	42
Approvals	43
Maintenance	43

Introduction

The Cole-Parmer TC9600 Advanced Multiparameter Temperature Controller (Models 89800-13 and 89800-14) is designed to regulate a user-defined set point temperature. The Ramp and Soak feature allow up to a 40-step profile using; time, hold, soak and end. Which allows you to create the ideal profile for your process. The microprocessor control uses a PID algorithm that provides precise heating/cooling tolerances for demanding applications in the laboratory or industrial settings. Advanced features include a USB communication port, audible alarm, external alarm relay output, detachable power cord. and grid support bracket that is ideal for fume hood mounting to maximize bench space.

The TC9600 is a microprocessor-based, digital indicating, single-loop temperature controller. It has a single output located on the back of the controller. It features an auto-tuning function that allows automatic configuration of the PID parameters with minimum user configuration required. The TC9600 has multiple inputs that accept thermocouple, RTD, and thermistor sensors.


The controller automatically sets the PID parameters through a "learning" sequence in the Tune mode. PID parameters include proportional band, reset/integral, and rate/derivative. The user-friendly LED display is to aid in monitoring and setup of the controller. The controller automatically store all user parameters in nonvolatile memory.

Unpacking

Check individual parts against the list of items below. If anything is missing or damaged, please contact your instrument supplier immediately.

1. Temperature controller
2. Grid support bracket (attached to the back of the controller)
3. 6-ft (3-m) detachable IEC power cord. Model 89800-13 comes standard with a US 120 VAC plug; Model 89800-14 comes standard with a US 220 VAC plug.
See page 41 for additional cords.
4. Electrical noise canceling ferrite clip (connects to sensor input; see page 40 for reference)
5. User manual
6. Quick-start guide

Controller Description — Front Panel

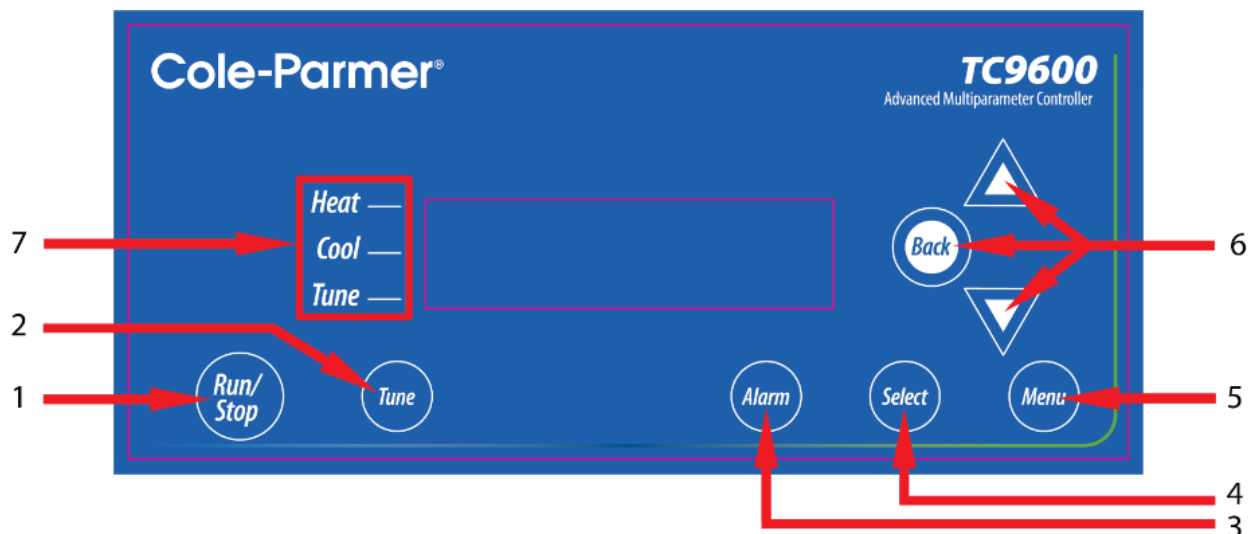

1. RUN/STOP

Pressing RUN/STOP once will start the control process if the Temperature Controller is stopped, or stop the control process if the Temperature Controller is running. If the Temperature Controller is running, “Heat” and “Cool” on-screen indicators will illuminate appropriately in the “Alarm/Action Display”.

2. TUNE Button

Pressing TUNE once will start the AUTO TUNE cycle. AUTO TUNING must be enabled in the setup mode for this key to function. (See page 14, Screen 16)

3. ALARM Button

In an ALARM situation, the screen will display either “WARNING” or “ERROR” with the corresponding message. A “WARNING” will not stop the control process. An “ERROR” will stop the control process.

- A. *MANUAL RESET* mode: Pressing ALARM once will silence the audible alarm and clear the on-screen alarm message. If the alarm situation is still present the “Alarm/Action Display” will remain illuminated. The alarm and on-screen message will not clear automatically, even if the system is no longer in an alarm situation.
- B. *AUTO RESET* mode: Pressing ALARM once will silence the audible alarm and clear the on-screen alarm message. If the alarm situation is still present, the “Alarm/Action Display” will remain illuminated. If the system leaves an alarm situation, the system will automatically silence the alarm and clear the on-screen alarm message.

4. SELECT Button

Pressing SELECT once will cycle through user-configurable control set points. All user-configurable set points will be underlined with a greyed outline. The selected Set Point will be underlined. Change the set point with the UP and DOWN arrow buttons.

5. MENU Button

The MENU button provides access to all user-configurable setup parameters of the controller. Pressing this key once will scroll through parameter options. Pressing and holding this key will exit to the home screen, saving any changes made up to that point.

6. UP, DOWN, BACK Buttons

The UP and DOWN arrow buttons will increase or decrease the value of the set point selected (underlined). Pressing the UP or DOWN arrow keys will increase or decrease *numerical* entries by the least significant digit. The rate of acceleration will increase as shown in the table below, starting from the least significant digit. Pressing and holding the UP or DOWN arrow key will increase or decrease *text* entries without an acceleration factor. The back button moves backwards through the General and Advanced Setup Menus. To exit to the main screen from either menu, press and hold the MENU button.


Numbers...	Increase/Decrease by...
0.0 – 0.9	0.1
0 – 9	1
10 – 100	10
100 +	50

Table 1. Acceleration Factor Table

7. HEAT, COOL, AND TUNE Indicators

When any of these modes are active they will have a block indicator on the display showing they are active. **Example:** When the control is in the heat mode and it is applying power to the heater output there will be an indicator block on the display to show the heater output is active.

Controller Description — Back Panel





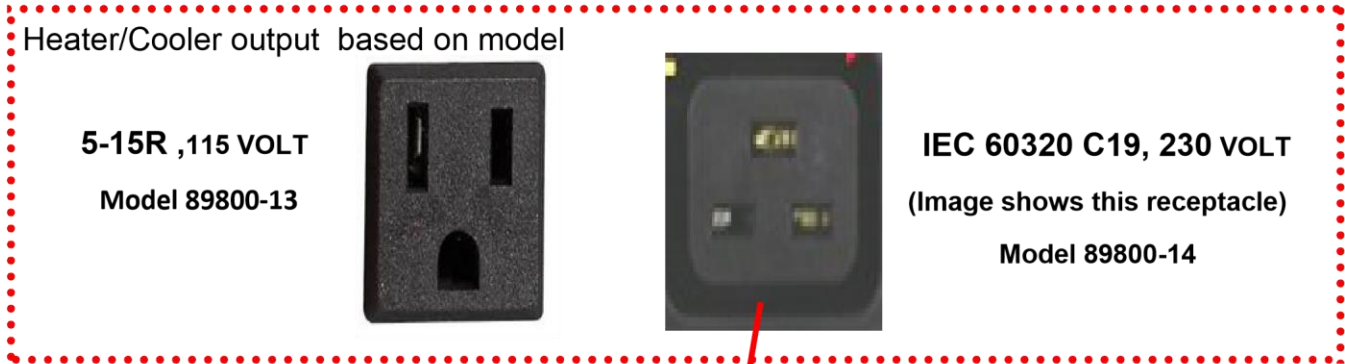


1. IEC power cord connection (see page 41 for additional cords)
2. Heater/cooler output (see pages 42 for optional cord adapters dependent on country)
3. Power switch
4. Grid support bracket
5. External alarm output
6. USB port
7. Thermistor input: accepts type 400 and 700 with Cole-Parmer phone plug connector
8. Thermocouple input: accepts type J, K, N, R, S, T, B, and E with mini-connector
9. RTD input: accepts type Alpha 0.003850, Alpha 0.003850 manufactured with MgO and Alpha 0.003916 with ANSI 3 blade connector. (RTD probes with ANSI connectors purchased from Cole-Parmer or Davis Instruments before May of 2013 will need the wiring in the probe connector modified. See page 40 for wiring instructions.)
10. RTD input: accepts Cole-Parmer 3-pin round connector
11. Fuse

Controller Description — Active Display Screen

MAIN SCREENS


RAMPING/SOAKING
INDICATOR
(R): Ramping
(S): Soaking

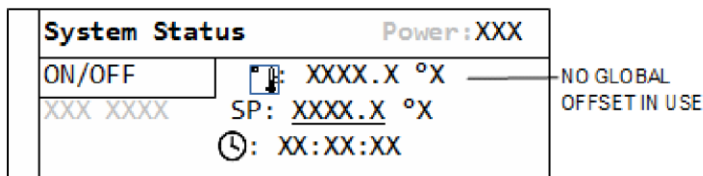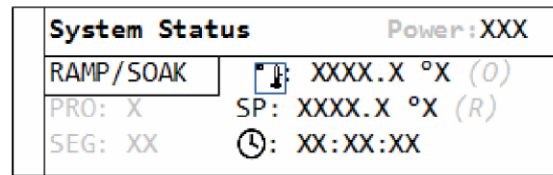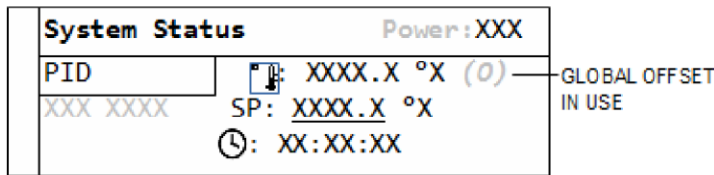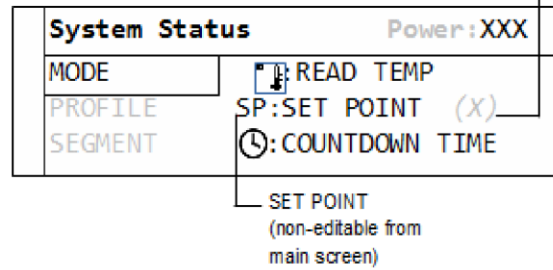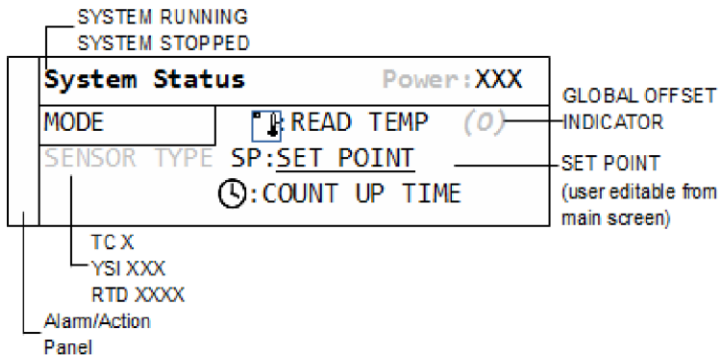

RAMP/SOAK MODE

PID and ON/OFF MODE

ALARM/ACTION PANEL

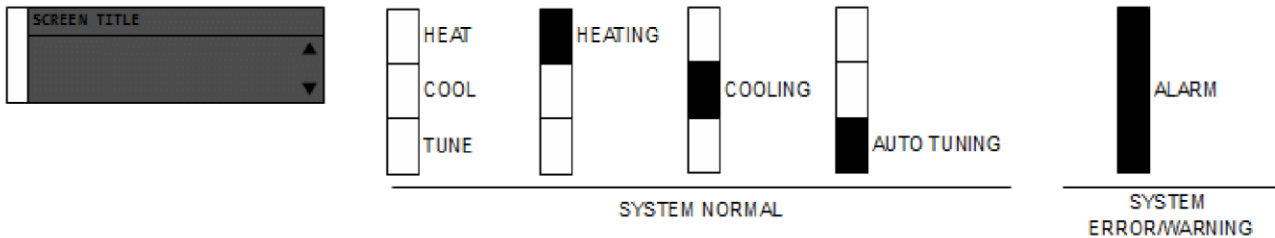

NAVIGATION



Setup and Operation

Initial Setup

- Install controller in a safe operating area.
- Plug the heating or cooling device (sold separately) into the output connector, located on the back of the controller.
- Connect input sensor into the proper connection input on the back of the controller.
- Place the ferrite clip over lead wire of sensor (see page 40 for reference picture of installation).
- Connect external alarm output to the controller (optional).
- Connect USB cable to the USB connector, located on the back of the controller (optional).
- Plug supplied AC cord into the IEC input connector, located on the back of the controller.

Basic Operation Setup

- Turn power switch ON, located on the back of controller.
- Follow the instructions on the “welcome” screen.
 - Press SELECT key to read a brief description of each key on the front of controller.
 - Press the MENU key to skip the instructions and enter into the main operation screen.
- Press the SELECT key to make a user-editable field active. A line will appear under the field when the field is active for editing.
- Use the UP/DOWN arrow keys to adjust the value that is active in a user-editable field.
- Enter the user-configurable setup by pressing the MENU key from the system status screen.
- Use the MENU key to advance through each menu setting.
- Hold UP and press DOWN arrow keys to jump directly to Output Power adjustment screen.
- Depending on process operation, follow the directions to set up the controller for the proper operation of your process:
 - PID (factory default) on pages 8-15
 - ON/OFF on pages 16-22
 - Ramp & Soak on pages 23-29
- All changed settings will be retained in memory when returning to the System Status screen.
- Use the screen flow charts (see pages 38-39) to have a visual of the controls menu layout.

Auto Tune Operation

- Set up your process as noted in the initial setup.
- Verify that the Auto Tune feature is enabled in the menu settings.
- From the main operation screen set the set point temperature.
- Press the TUNE button and the indicators showing Heat and Tune will be illuminated on display.
- The Set Point value cannot be altered after the Tune process has started. The value is locked until the Tune process is complete or aborted by the user.
- Please note that the controller will tune at 90% of the set point. Don't be alarmed if the controller does not reach your entered set point in the tune mode.
- Stopping the Tune operation prior to it finishing will cause the PID settings to be returned to factory default values.
- PID values will be saved in nonvolatile memory.

PID Control Mode Setup:

(FACTORY DEFAULT MODE)

Screen # 1 - Selecting Sensor Type

- Press the MENU key one time to select sensor input type.
- Use UP/DOWN arrow keys to change sensor type.
- Press MENU key when complete.

Sensor Type
<u>THERMOCOUPLE</u>

Screen # 2 - Selecting Sensor Type

- If **Thermocouple** is selected on screen # 1:
- Use the UP/DOWN arrow keys to select the desired thermocouple type: J, K, N, R, S, T, B, or E.
- Press MENU key to advance to screen #3 after selection has been made.

Thermocouple Type
Type <u>J</u>
Range: ±XXX to XXXX °X

- If **Thermistor** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired thermistor type: 400 or 700.
- Press MENU key to advance to screen #3 after selection has been made.

Thermistor Type
Type <u>400</u>
Range: ±XXX to XXXX °X

- If **RTD** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired RTD type: Alpha 003850, 003850 w/ MgO, or 003916.
- Press MENU key to advance to screen #3 after selection has been made.

RTD Type
Al pha <u>003850</u>
Range: ±XXX to XXXX °X

Screen # 3 - Selecting Temperature Scale


- Use the UP/DOWN keys to select the desired temperature scale: Celsius °C, Fahrenheit °F, Kelvin K, Reaumur °Ré, or Rankine °Ra.
- Press MENU key to advance to screen # 4 after selection has been made.

Temperature Scale
<u>CELSIUS °C</u>

PID Control Mode Setup:


Screen # 4 - Selecting Alarm Setting

- Use the UP/DOWN arrow keys to change between Auto Reset, Manual Reset and Off.
- Press MENU key to advance to next screen.
 - If Auto or Manual is selected, the controller will advance to screen # 5.
 - If selection is OFF, the controller will advance to screen # 9.

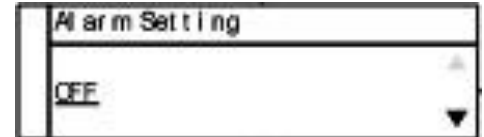

Screen # 5 - Selecting Alarm Mode

- Use the UP / DOWN arrow keys to select alarm mode.
- Selectable alarm modes:
 - **Process High:** The alarm will be activated if the temperature rises above the user set alarm value.
 - **Process Low:** The alarm will be activated if the temperature goes below the user set alarm value.
 - **Process High/Low:** The alarm will be activated if the temperature goes above or below the high and low temperature levels set by user.
 - **Deviation High:** Alarm activation point that changes with the change of the process set point.
Example: The user sets deviation high alarm at 30°C and has a process set point of 100° C. If the temperature reaches 130°C, the alarm will be activated. The user changes process set point to 60°C. If process temperature reaches 90°C, the alarm will be activated.
 - **Deviation Low:** Alarm activation point that changes with the change of the process set point.
Example: The user sets the deviation low alarm at 30°C and the process set point is 100° C. If the process temperature is 70°C or less, the alarm will be activated. If the process set point temperature is set at 70°C, then if the process temperature is 40°C or less, it will activate the alarm.
 - **Deviation High/Low:** Same as Deviation High and Deviation Low alarms. It has the capability of having an alarm activate when the set point temperature goes above or below a set point temperature.
 - **Deviation Band:** Allows the user to set the alarm to be activated within a specified temperature range.
- Press the MENU key to advance to screen # 6 after selecting alarm mode.

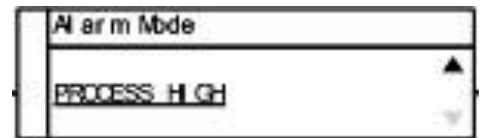

PID Control Mode Setup:

Screen # 6 - Setting Alarm Mode Variables

- This screen reflects the variable settings for the selected alarm mode from screen # 5.

-Set Point—Process High

- Press the UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Process High	▲
HIGH <u>XXXX.X</u> °X	▼

-Set Point—Process Low

- Press the UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Process Low	▲
LOW <u>XXXX.X</u> °X	▼

-Set Point—Process High/Low

- Press the UP/DOWN arrow keys to adjust the temperature value.
- Press the SELECT key to toggle between High and Low. There will be a line below the active field.
- When complete, press MENU key to advance to screen # 7.

Process High/Low	▲
High: <u>XXXX.X</u> °X	▼
Low: <u>XXXX.X</u> °X	

-Set Point—Deviation High

- Press the UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation High	▲
HIGH <u>XXXX.X</u> °X	▼

-Set Point—Deviation Low

- Press the UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation Low	▲
LOW <u>XXXX.X</u> °X	▼

PID Control Mode Setup:

-Set Point—Deviation High/Low

- Press the UP/DOWN arrow keys to adjust the temperature value.
- Press the SELECT key to toggle between the High and Low. There will be a line below the active field.
- When complete, press MENU key to advance to screen # 7.

Deviation High/Low	
High: XXXX.X °X	▲
Low: XXXX.X °X	▼

-Set Point—Deviation Band

- Press the UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation Band	
BAND: XXXX.X °X	▲
	▼

Screen # 7 - Setting Hysteresis Alarm

- Use the UP/DOWN arrow keys to adjust value.
- The user-defined value when the alarm will exit an alarm condition. **Example:** User has a Process High alarm programmed with a value of 150°C and has set the Alarm Hysteresis of 1°C. The alarm has been activated after the temperature reached 160°C. When temperature reaches 149°C, the alarm condition will not be active.
- When complete, press MENU key to advance to screen # 8.

Alarm Hysteresis	
X.XXXX °X	▲
	▼

Screen # 8 - Setting Audible Alarm


- Use UP/DOWN arrow keys to change value to either ON audible alarm or OFF audible alarm.
- When complete, press MENU key to advance to screen # 9.

Audible Alarm	
ON	▲
	▼

PID Control Mode Setup:

Screen # 9 - Advanced Menu Gateway

- Use the UP/DOWN arrow keys to change selection to ENTER.
- When complete, press MENU key to advance to screen # 10.
- The following screens are in the advance portion of the setup menu:

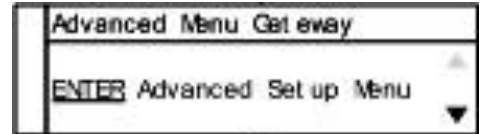

Screen # 10 - Calibration Gateway

Screen # 11 - Global Sensor Offset

Screen # 12 - Over Temperature Stop

Screen # 13 - Loop Break Stop

Screen # 14 - Control Action

Screen # 15 - Control Mode

Screen # 16 - Auto Tune

Screen # 17 - Proportional Gain Setup

Screen # 18 - Integral Gain Setup

Screen # 19 - Derivative Rate Gain Setup

Screen # 20 - Power Rate Control


Screen # 21 - Run Time

Screen # 22 - Power Failure Control

PID Control Mode Setup:

Screen # 10 - Calibration Gateway

- Screen will default to SKIP Calibration Setup.
- **WARNING:** Do not enter the calibration gateway unless you have the proper calibration equipment to calibrate the sensor inputs associated with this controller. Changing calibration settings in the controller can cause errors in sensor temperature readings and operation of the control. For instructions on how to use the calibration functions, contact your instrument supplier.
- With SKIP selected, press the MENU key to advance to screen # 11.

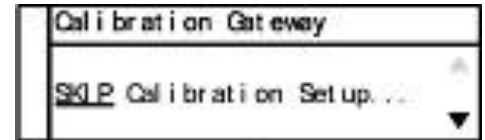


Screen # 11 - Global Sensor Offset Calibration

- Use the UP/DOWN arrow keys to change the offset value.
- The use of another temperature device is required to determine the correct offset needed.
- Adjusting this value is offsetting the sensor input temperature being displayed.
- When complete, press MENU key to advance to screen # 12.




Screen # 12 - Over Temperature Stop

- Use UP/DOWN arrow keys to change the value.
- When sensor temperature goes above this value, the output device will be turned off.
- When complete, press MENU key to advance to screen # 13.

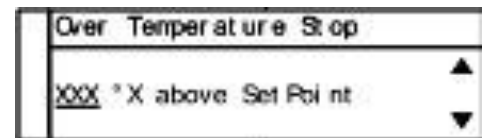

Screen # 13 - Loop Break Stop


- Default value is ENABLED. Use UP/DOWN arrow keys to change the value.
- Loop Break Stop is a feature that will stop the process output while the controller is in the run mode. If the controller senses there is no change in temperature in a user set amount of time, the output will be turned off and error message will be displayed on the screen along with an audible alarm.
- Set to DISABLED if this function is not required.
- When complete, press MENU key to advance to screen # 14.

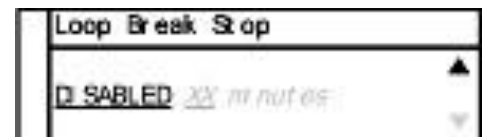

PID Control Mode Setup:


Screen # 14 - Control Action

- Use UP/DOWN arrow keys to change the value.
- Set for the process that is being performed: Heat or Cool
- When complete, press MENU key to advance to screen # 15.

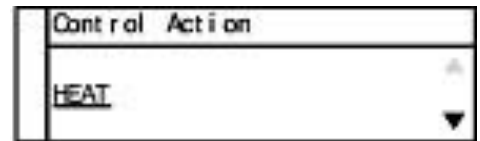

Screen # 15 - Control Mode

- Use UP/DOWN arrow keys to change the value.
- Default value is PID mode. (If different, select PID mode)
- When complete, press MENU key to advance to screen # 16.

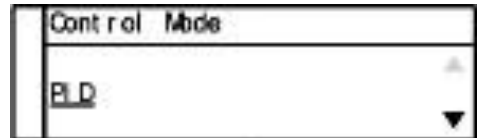


Screen # 16 - Auto Tune

- Default value is ENABLED. Use UP/DOWN arrow keys to toggle between ENABLED and DISABLED.
- Select ENABLED to use the Tune key on the front panel to activate the Auto Tune function. This function needs to remain ENABLED until the controller has been properly tuned to the user's particular application.
- NOTE: the control auto tunes to 90% of the set point.
- When complete, press MENU key to advance to screen # 17.




Screen # 17 - Proportional Gain Setup

- Use UP/DOWN arrow keys to change the value.
- **WARNING:** Do not change these values unless you have experience in manual setup of PID (Proportional, Integral, and Derivative) of this controller. These values are derived from running the auto tune function. Altering the value after auto tune has been completed will affect the performance of the system.
- When complete, press MENU key to advance to screen # 18.

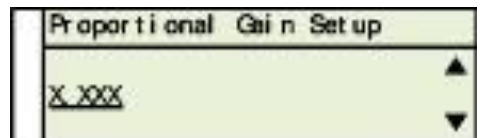

Screen # 18 - Integral Gain Setup


- Use UP/DOWN arrow keys to change the value.
- **WARNING:** Do not change these values unless you have experience in manual setup of PID (Proportional, Integral, and Derivative) of this controller. These values are derived from running the auto tune function. Altering the values after auto tune has been completed will affect the performance of the system.
- When complete, press MENU key to advance to screen # 19.

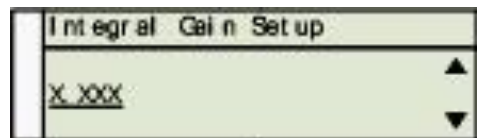

PID Control Mode Setup:


Screen # 19 - Derivative Rate Gain Setup

- Use UP/DOWN arrow keys to change the value.
- **WARNING:** Do not change these values unless you have experience in manual setup of PID (Proportional, Integral, and Derivative) of this controller. These values are derived from running the auto tune function. Altering the value after auto tune has been completed will affect the performance of the system.
- When complete, press MENU key to advance to screen # 20.

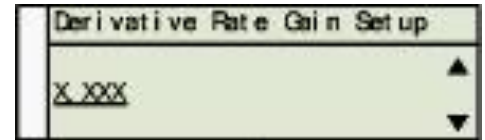


Screen # 20 - Power Output Control

- Use UP/DOWN arrow keys to change the value.
- Default value is 100 percent.
- This feature allows the user to reduce the output power to the output heating device—useful in applications where temperature overshoot is occurring after auto tune has been completed.
- Experimenting with this value will be necessary because every process setup is different.
- When complete, press MENU key to advance to screen # 21.

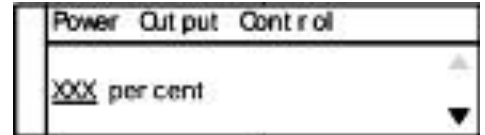


Screen # 21 - Run Time

- Use UP/DOWN arrow keys to change the value.
- Use SELECT key to toggle between the HR and MIN fields.
- Run Time is a safety feature while the control is in the ON/OFF or PID mode. The user can set a timer to have the output turned OFF if the time expires.
- When complete, press MENU key to advance to screen # 22.

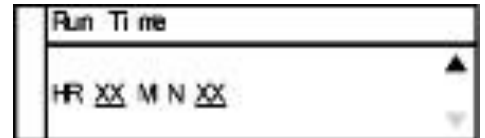

Screen # 22 - Power Failure Control

- Use UP/DOWN arrow keys to change the value.
- Default value is STOP:
 - STOP* — If the controller loses power while in the run mode, it will not resume the process once the power has been restored.
 - RESUME* — If the controller loses power while in the run mode, it will resume the operating process once the power has been restored.
- When complete, press MENU key to advance to screen # 23.

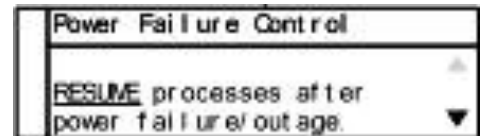


Screen # 23 - Advanced Menu Exit

ON/OFF Control Mode Setup


Screen # 1 - Selecting Sensor Type

- Press the MENU key one time to select sensor input type.
- Use UP/DOWN arrow keys to change sensor type.
- When complete, press MENU key to advance to next screen.

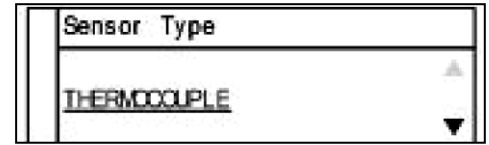


Screen # 2 - Selecting Sensor Type


- If **Thermocouple** is selected on screen # 1:
- Use the UP/DOWN arrow keys to select the desired thermocouple type: J, K, N, R, S, T, B, or E.
- Press MENU key to advance to screen #3 after selection has been made.

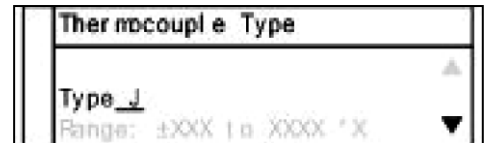

- If **Thermistor** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired thermistor type: 400 or 700.
- Press MENU key to advance to screen #3 after selection has been made.

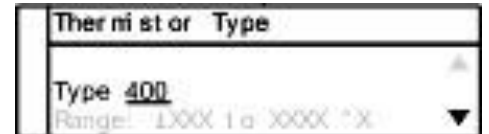


- If **RTD** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired RTD type: Alpha 003850, 003850 w/ MgO, or 003916.
- Press MENU key to advance to screen #3 after selection has been made.

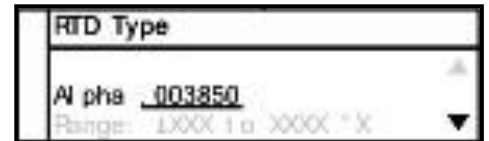


Screen # 3 - Selecting Temperature Scale

- Use the UP/DOWN keys to select the desired temperature scale: Celsius °C, Fahrenheit °F, Kelvin K, Reaumur °Ré, or Rankine °Ra.
- Press MENU key to advance to screen # 4 after selection has been made.

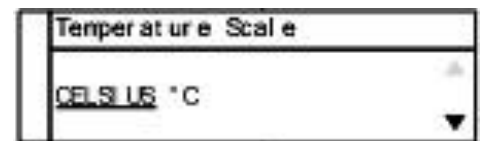

Screen # 4 - Selecting Alarm Setting


- Use the UP/DOWN arrow keys to change between Auto Reset, Manual Reset, and Off.
- Press MENU key to advance to next screen:
 - If Auto or Manual is selected, the controller will advance to screen # 5.
 - If selection is OFF, the controller will advance to screen # 9.

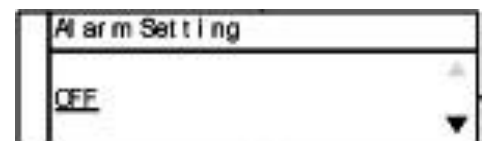

ON/OFF Control Mode Setup

Screen # 5 - Selecting Alarm Mode

- Use the UP/DOWN arrow keys to select alarm mode.
- Selectable alarm modes:

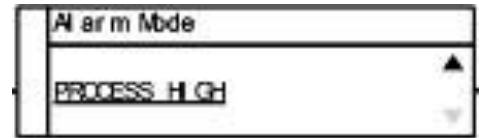

- **Process High:** The alarm will be activated if the temperature rises above the user set alarm value.
 - **Process Low:** The alarm will be activated if the temperature goes below the user set alarm value.
 - **Process High/Low:** The alarm will be activated if the temperature goes above or below the high and low temperature levels set by user.
 - **Deviation High:** Alarm activation point that changes with the change of the process set point. **Example:** The user sets deviation high alarm at 30°C and has a process set point of 100°C. If the temperature reaches 130°C, the alarm will be activated. The user changes process set point to 60°C. If process temperature reaches 90°C, the alarm will be activated.
 - **Deviation Low:** Alarm activation point that changes with the change of the process set point. **Example:** User sets the deviation low alarm 30°C and the process set point is 100°C. If the process temperature is 70°C or less, the alarm will be activated. If the process set point temperature is set at 70°C, then if the process temperature is 40°C or less, it will activate the alarm.
 - **Deviation High/Low:** Same as Deviation High and Deviation Low alarms. It has the capability of having an alarm activate when the set point temperature goes above or below a set point temperature.
 - **Deviation Band:** Allows the user to set the alarm to be activated within a specified temperature range.
- Press the MENU key to advance to screen # 6 after selecting alarm mode.

ON/OFF Control Mode Setup

Screen # 6 - Setting Alarm Mode Variables

- This screen reflects the variable settings for the selected alarm mode from screen # 5.

-Set Point—Process High

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Process High
HIGH <u>XXXX.X</u> °X

- Set Point—Process Low

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Process Low
LOW <u>XXXX.X</u> °X

- Set Point—Process High/Low

- Use UP/DOWN arrow keys to adjust the temperature value.
- Press the SELECT key to toggle between High and Low. There will be a line below the active field.
- When complete, press MENU key to advance to screen # 7.

Process High/Low
High: <u>XXXX.X</u> °X
Low: <u>XXXX.X</u> °X

- Set Point—Deviation High

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation High
HIGH <u>XXXX.X</u> °X

- Set Point—Deviation Low

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation Low
LOW <u>XXXX.X</u> °X

-Set Point—Deviation High/Low


- Use UP/DOWN arrow keys to adjust the temperature value.
- Press the SELECT key to toggle between High and Low. There will be a line below the active field.
- When complete, press MENU key to advance to screen # 7.

Deviation High/Low
High: <u>XXXX.X</u> °X
Low: <u>XXXX.X</u> °X

ON/OFF Control Mode Setup


- Set Point—Deviation Band

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press the key to advance to screen # 7.

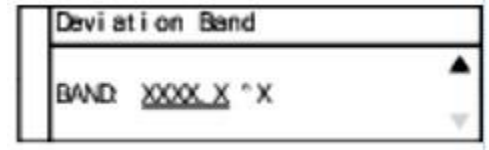


Screen # 7 - Setting Hysteresis Alarm

- Use the UP/DOWN arrow keys to adjust value.
- The user-defined value when the alarm will exit an alarm condition. **Example:** The user has a Process High alarm programmed with a value of 150°C and has set the Alarm Hysteresis of 1°C. The alarm has been activated after the temperature reached 160°C. When temperature reaches 149°C, the alarm condition will not be active.
- When complete, press MENU key to advance to screen # 8.

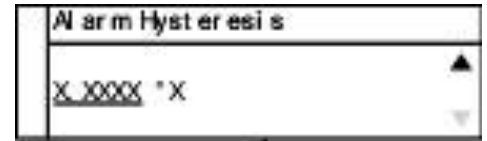


Screen # 8 - Setting Audible Alarm

- Use UP/DOWN arrow keys to change value to either ON audible alarm or OFF audible alarm.
- When complete, press MENU key to advance to screen # 9.

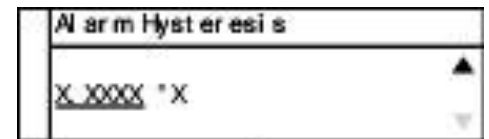

Screen # 9 - Advanced Menu Gateway

- Use UP/DOWN arrow keys to change selection to ENTER.
- When complete, press MENU key to advance to screen # 10.

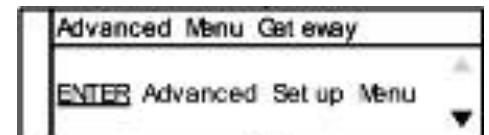

Screen # 10 - Calibration Gateway

Screen # 11 - Global Sensor Offset

Screen # 12 - Over Temperature Stop

Screen # 13 - Loop Break Stop

Screen # 14 - Control Action

Screen # 15 - Control Mode

Screen # 16 - On/Off Control


Screen # 17 - Power Output Control

Screen # 18 - Run Time

ON/OFF Control Mode Setup:


Screen # 10 Calibration Gateway

- Screen will default to SKIP Calibration Setup.
- **WARNING:** Do not enter the calibration gateway unless you have the proper calibration equipment to calibrate the sensor inputs associated with this controller. Changing calibration settings in the control can cause errors in sensor temperature readings and operation of the control. For instructions on how to use the calibration functions, contact your instrument supplier.
- With SKIP selected, press the MENU key to advance to screen # 11.

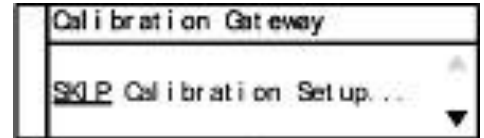


Screen # 11 - Global Sensor Offset Calibration

- Use the UP/DOWN arrow keys to change the offset value.
- The use of another temperature device is required to determine the correct offset needed.
- Adjusting this value is offsetting the sensor input temperature being displayed.
- When complete, press MENU key to advance to screen # 12.

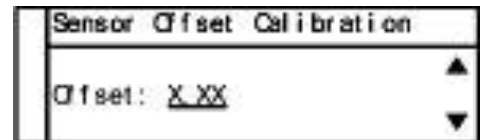


Screen # 12 - Over Temperature Stop

- Use the UP/DOWN arrow keys to change the value.
- When sensor temperature goes above this value, the output device will be turned off.
- When complete, press MENU key to advance to screen # 13.

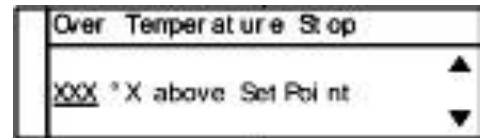

Screen # 13 - Loop Break Stop


- Default value is ENABLED. Use the UP/DOWN arrow keys to change the value.
- Loop Break Stop is a feature that will stop the process output while the control is in the run mode. If the control senses there is no change in temperature in a user set amount of time the output will be turned off and error will be displayed on the screen along with a audible alarm.
- Set to DISABLED if this function is not required.
- When complete, press MENU key to advance to screen # 14.

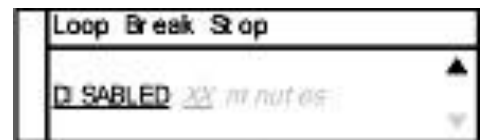

ON/OFF Control Mode Setup:


Screen # 14 Control Action

- Use the UP/DOWN arrow keys to change the value from Heat to Cool.
- Set for the particular process that is being performed.
- When complete, press MENU key to advance to screen # 15.

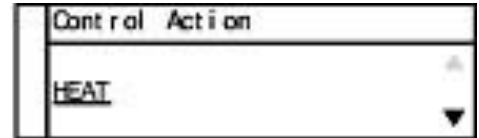


Screen # 15 - Control Mode

- Use UP/DOWN arrow keys to change the value from the default mode of PID to On/Off.
- When complete, press MENU key to advance to screen # 16.

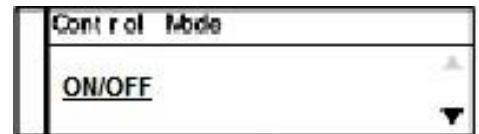


Screen # 16 - ON/OFF Control

- Use the UP/DOWN arrow keys to adjust value.
- Cycle time is the rate that the output cycles on and off. The recommended time is 1 second.
- If a mechanical relay is being used from the output, it is recommended that the time be increased to more than 1 second. Mechanical relays are slower than solid-state relays and require more time to operate correctly. Damage to the relay and control can occur if not set correctly.
- When complete, press MENU key to advance to screen # 17.

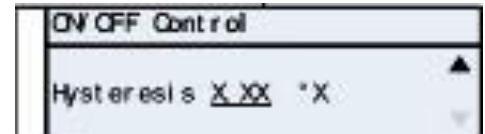


Screen # 17 - Power Output Control

- Use UP/DOWN arrow keys to change the value.
- Default value is 100 percent.
- This feature allows the user to reduce the output power to the output heating device—useful in applications where temperature overshoot is occurring after Auto Tune has been completed.
- Experimenting with this value will be necessary because every process setup is different.
- When complete, press MENU key to advance to screen # 18.

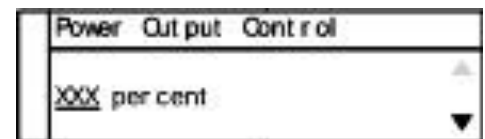

Screen # 18 Run Time

- Use the UP/DOWN arrow keys to change the value.
- Use the SELECT key to toggle between HR and MIN fields.

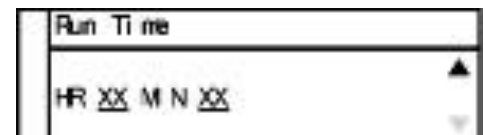


ON/OFF Control Mode Setup:

- Run Time is a safety feature while the control is in the On/Off or PID mode. The user can set a timer to have the output turned OFF if the time expires.
- When complete, press MENU key to advance to screen # 19.

Screen # 19 - Power Failure Control

- Use the UP/DOWN arrow keys to change the value.
- Default value is STOP:

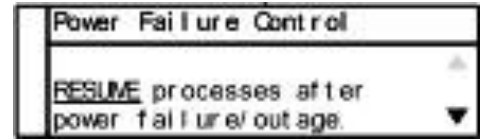

STOP — If the controller loses power while in the run mode, it will not resume the process once the power has been restored.

RESUME — If the controller loses power while in the run mode, it will resume the operating process once the power has been restored.

- When complete, press MENU key to advance to screen # 20.


Screen # 20 - Advanced Menu Exit

- Setup is complete.
- Automatically will exit to the main operation screen.

Ramp & Soak Mode Setup:

Screen # 1 - Selecting Sensor Type

- Press the MENU key one time to select sensor input type.
- Use UP/DOWN arrow keys to change sensor type.
- Press MENU key when complete.

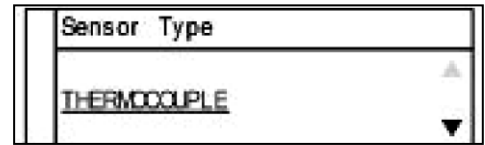


Screen # 2 - Selecting Sensor Type


- If **thermocouple** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired thermocouple type: J, K, N, R, S, T, B, or E.
- Press MENU key to advance to screen # 3 after selection has been made.



- If **thermistor** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired thermistor type: 400 or 700.
- Press MENU key to advance to screen # 3 after selection has been made.

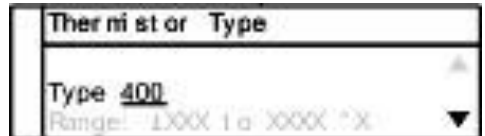


- If **RTD** is selected on screen # 1:
- Use UP/DOWN arrow keys to select the desired RTD type: Alpha 003850, 003850 w/ MgO, or 003916.
- Press menu to advance to screen # 3 after selection has been made.

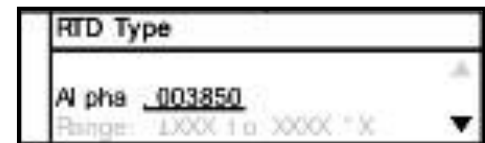

Screen # 3 - Selecting Temperature Scale


- Use the UP/DOWN keys to select the desired temperature scale: Celsius °C, Fahrenheit °F, Kelvin K, Reaumur °Ré, or Rankine °Ra.
- Press MENU key to advance to screen # 4 after selection has been made.

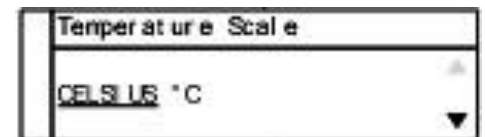

Ramp & Soak Mode Setup:


Screen # 4 - Selecting Alarm Setting

- Use the UP/DOWN arrow keys to change between Auto Reset and Manual Reset.
- Press MENU key to advance to next screen:
 - If Auto Reset or Manual Reset was selected, continue to screen # 5.
 - If selection is OFF, skip to screen # 9.

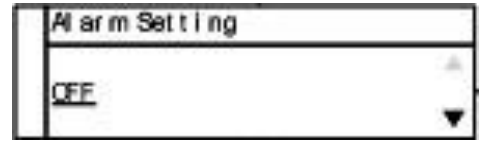

Screen # 5 - Alarm Mode

- Use the UP/DOWN arrow keys to select alarm mode.
- Selectable alarm modes:
 - **Process High:** The alarm will be activated if the temperature rises above the user set alarm value.
 - **Process Low:** The alarm will be activated if the process value goes below the user set alarm value.
 - **Process High/Low:** The alarm will be activated if the temperature goes above or below a high and low temperature set by user.
 - **Deviation High:** Alarm activation point that changes with the change of the process set point. Example: The user sets deviation high alarm at 30°C and has a process set point of 100°C. If the temperature reaches 130°C the alarm will be activated. The user changes process set point to 60°C. If process temperature reaches 90°C the alarm will be activated.
 - **Deviation Low:** Alarm activation point that changes with the change of the process set point. Example: The user sets the deviation low alarm 30°C and the process set point is 100°C. If the process temperature is 70°C or less, the alarm will be activated. If the process set point temperature is set at 70°C, then if the process temperature is 40°C or less it will activate the alarm.
 - **Deviation High/Low:** Same as Deviation High and Deviation Low alarms. It has the capability of having an alarm activate when the set point temperature goes above or below a set point temperature.
 - **Deviation Band:** Allows the user to set the alarm to be activated within a specified temperature range.
- Press the menu key to advance to screen # 6 after selecting alarm mode.

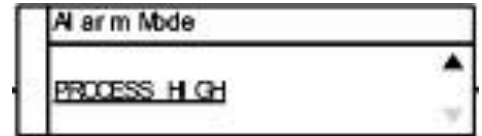

Ramp & Soak Mode Setup:

Screen # 6 - Setting Alarm Mode Variables

- This screen reflects the variable settings for the selected alarm mode from screen # 5.

-Set Point—Process High

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Process High
HIGH <u>XXXX.X</u> °X

- Set Point—Process Low

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Process Low
LOW <u>XXXX.X</u> °X

- Set Point—Process High/Low

- Use UP/DOWN arrow keys to adjust the temperature value.
- Press the SELECT key to toggle between High and Low. There will be a line below the active field.
- When complete, press MENU key to advance to screen # 7.

Process High/Low
High: <u>XXXX.X</u> °X
Low: <u>XXXX.X</u> °X

- Set Point—Deviation High

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation High
HIGH <u>XXXX.X</u> °X

- Set Point—Deviation Low

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press MENU key to advance to screen # 7.

Deviation Low
LOW <u>XXXX.X</u> °X

Set Point—Deviation High/Low


- Use UP/DOWN arrow keys to adjust the temperature value.
- Press the SELECT key to toggle between High and Low. There will be a line below the active field.
- When complete, press MENU key to advance to screen # 7.

Deviation High/Low
High: <u>XXXX.X</u> °X
Low: <u>XXXX.X</u> °X

Ramp & Soak Mode Setup:


- Set Point—Deviation Band

- Use UP/DOWN arrow keys to adjust the temperature value.
- When complete, press the key to advance to screen # 7.

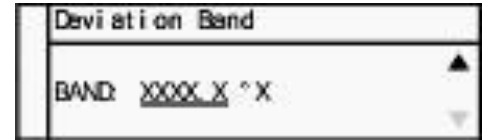


Screen # 7 - Setting Hysteresis Alarm

- Use the UP/DOWN arrow keys to adjust value.
- The user-defined value when the alarm will exit an alarm condition. **Example:** The user has a Process High alarm programmed with a value of 150°C and has set the Alarm Hysteresis of 1°C. The alarm has been activated after the temperature reached 160°C. When temperature reaches 149°C, the alarm condition will not be active.
- When complete, press MENU key to advance to screen # 8.

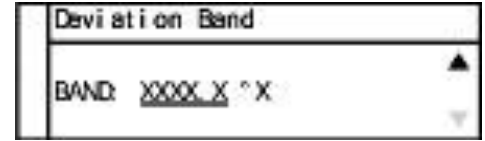


Screen # 8 - Setting Audible Alarm

- Use UP/DOWN arrow keys to change value to either ON audible alarm or OFF audible alarm.
- When complete, press MENU key to advance to screen # 9.

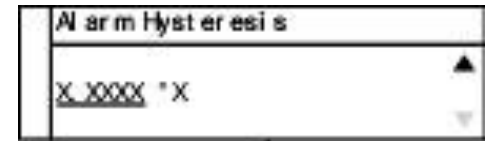

Screen # 9 - Advanced Menu Gateway

- Use UP/DOWN arrow keys to change selection to ENTER.
- When complete, press MENU key to advance to screen # 10.
- The following screens are in the advanced portion of the setup menu:

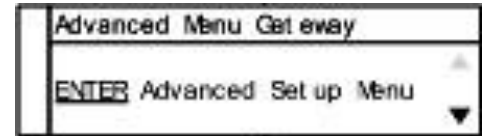

Screen # 10 - Calibration Gateway

Screen # 19– Run Time

Screen # 11 - Global Sensor Offset

Screen # 20 - Power Failure Control

Screen # 12 - Over Temperature Stop

Screen # 21 - Advanced Menu Exit

Screen # 13 - Loop Break Stop

Screen # 14 - Control Action

Screen # 15 - Control Mode


Screen # 16 - Ramp & Soak Profile

Screen # 17 - Assured Soak

Ramp & Soak Mode Setup:


Screen # 10 - Calibration Gateway

- Screen will default to SKIP Calibration Setup.
- **WARNING:** Do not enter the calibration gateway unless you have the proper calibration equipment to calibrate the sensor inputs associated with this controller. Changing calibration settings in the control can cause errors in sensor temperature readings and operation of the control. For instructions on how to use the calibration functions, contact your instrument supplier.
- With SKIP selected, press the MENU key to advance to screen # 11.

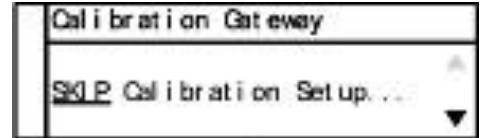


Screen # 11 - Global Sensor Offset Calibration

- Use the UP/DOWN arrow keys to change the offset value.
- The use of another temperature device is required to determine the correct offset needed.
- Adjusting this value is offsetting the sensor input temperature being displayed.
- When complete, press MENU key to advance to screen # 12.

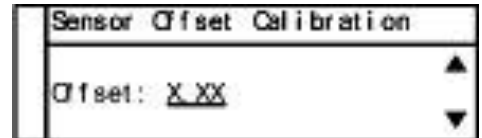


Screen # 12 - Over Temperature Stop

- Use the UP/DOWN arrow keys to change the value.
- When sensor temperature goes above this value, the output device will be turned off.
- When complete, press MENU key to advance to screen # 13.

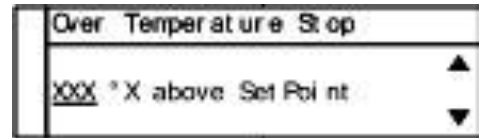

Screen # 13 - Loop Break Stop


- Default value is ENABLED. Use the UP/DOWN arrow keys to change the value.
- Loop Break Stop is a feature that will stop the process output while the control is in the run mode. If the control senses there is no change in temperature in a user set amount of time the output will be turned off and error will be displayed on the screen along with a audible alarm.
- Set to DISABLED if this function is not required.
- When complete, press MENU key to advance to screen # 14.

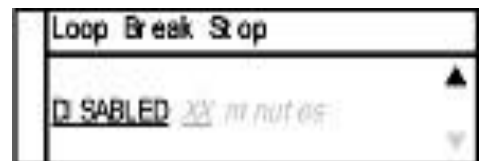

Ramp & Soak Mode Setup:


Screen # 14 - Control Action

- Use the UP/DOWN arrow keys to change the value from Heat to Cool.
- Set for the particular process that is being performed.
- When complete, press MENU key to advance to screen # 15.

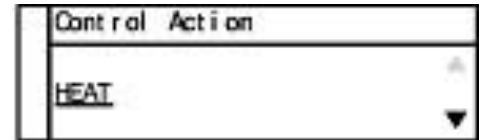


Screen # 15 - Control Mode

- Use UP/DOWN arrow keys to change the value from the default mode of PID to On/Off.
- When complete, press MENU key to advance to screen # 16.

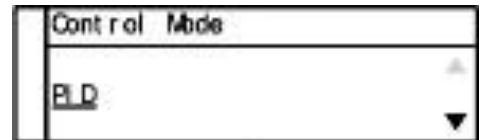


Screen # 16 - Ramp & Soak Profile

- Use UP/DOWN arrow keys to change the profile number.
- Profiles # 1 through # 9 are available to set up for different processes.
- When complete, press MENU key to advance to screen # 17.

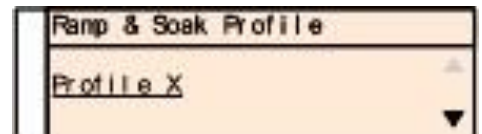

Screen # 17 - Assured Soak

- Use UP/DOWN arrow keys to change value.
- If Assured Soak is set to YES, the profile will not advance to the next step until it reaches a certain temperature set in the profile.
If set to NO, the profile will continue on even if the soak temperature has not been met.
- When complete, press MENU key to advance to screen # 18.

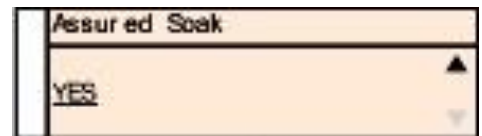

Ramp & Soak Mode Setup:

Screen # 18 - Segment Definition Setup

- Use UP/DOWN arrow keys to change the value.
- Use the SELECT key to change between user-defined sections:

Segment Set up	Power: XXX
SEG XX	XXXX.X °X ▲
End Time: HR XX MN XX	
Rate: XXX.X	▼

- **SEG (Segment):** There are 1 to 16 programmable segments per profile.
- **Temp (Temperature):** Controller is expected to meet this temperature within a user defined time. Two user-editable fields: HR (Hour) and MIN (Minutes).
- **End Time:** Time control is expected to complete segment.
- **Rate:** User defined rate per segment, degrees per minute.
- **Power Rate:** Sets the output power duty cycle for the selected segment.

Press Menu key to advance to next screen.

Screen # 19 - Run Time

- Use the UP/DOWN arrow keys to change the value.
- Use the SELECT key to toggle between HR and MIN fields.
- Run Time is a safety feature while the control is in the On/Off or PID mode. The user can set a timer to have the output turned OFF if the time expires.
- When complete, press MENU key to advance to screen # 20.

Run Time
HR XX MN XX ▲
▼

Screen # 20 - Power Failure Control

- Use the UP/DOWN arrow keys to change the value.
- Default value is STOP:

STOP — If the controller loses power while in the run mode, it will not resume the process once the power has been restored.

RESUME — If the controller loses power while in the run mode,

it will resume the operating process once the power has been restored.

- When complete, press MENU key to advance to screen # 21.

Power Failure Control
RESUME processes after power failure/outage. ▲
▼

Screen # 21 - Advanced Menu Exit


Data Acquisition Software Installation

1. Download software from Cole-Parmer website. Search for product 89800-13 or 89800-14.
2. Data Acquisition software is located in Technical Resources section of the product page.
3. Extract the downloaded file after download is complete.
To install the software execute the setup.exe file.

4. Press the Next button to continue with the installation of software.



5. Press the Next button or make any changes needed for your installation.

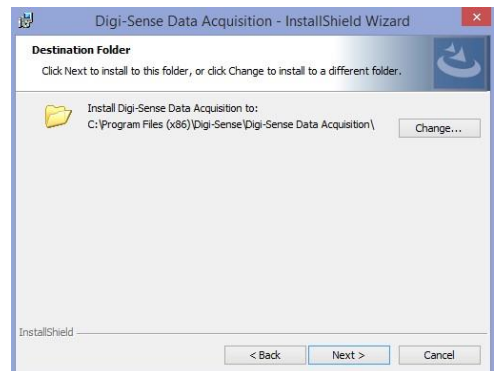

6. Press the Next button to confirm the installation of the software.




7. Installation is complete. Press the Finish button to close Installation window.
8. Software will be located in Cole-Parmer folder

Data Acquisition Software COM Port Setup

1. Connect USB cable between the Cole-Parmer TC9600 controller and your computer.
2. Turn the controller ON.
3. Start the Data Acquisition software.
4. Click the COM Port Settings button in the software and select the appropriate COM port.

COM Port Settings...

5. If you do not know the COM port the controller is using, follow these steps to select correct port:
 - Go into Device Manager. To enter device manager, go to the START button and type the word device manager.
 - Click on Device Manager
 - In Device Manager, find the option POTS (COM & LPT), then click on the plus symbol.
 - Find the associated port that has been assigned by the computer to the controller.

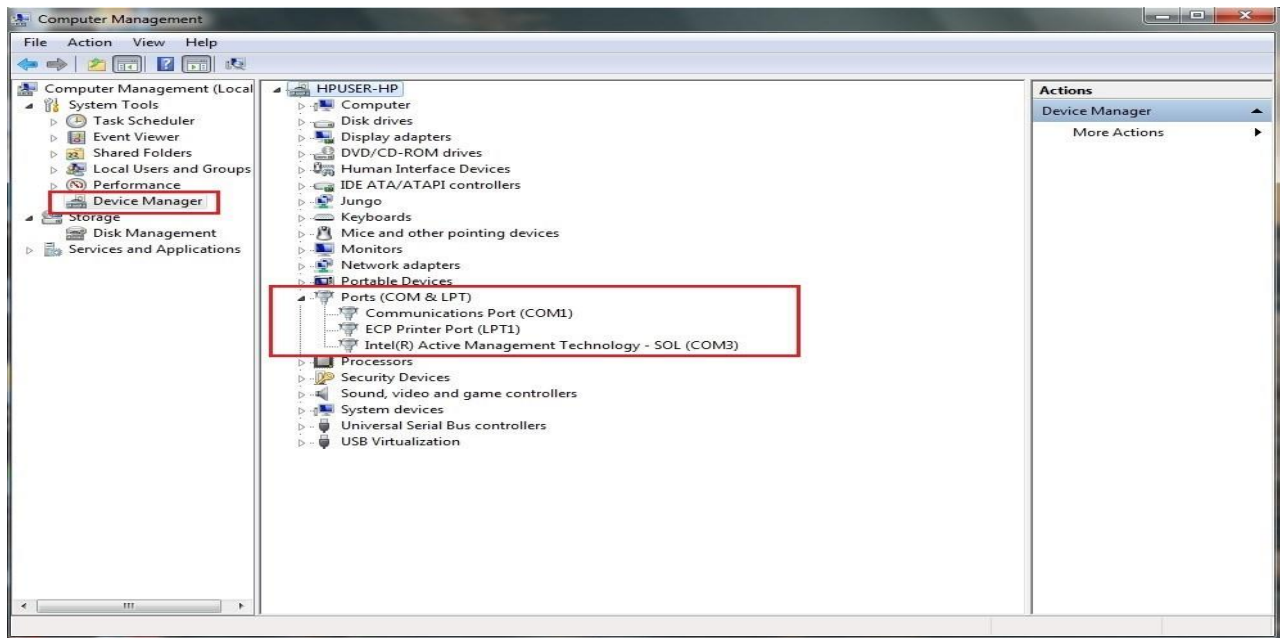

6. After COM port has been determined, press the Connect button.



7. If connection is successful, the sensor temperature will appear in the current temp location and status will change to connected.

Data Acquisition Software Setup

Main operation screen:

The screenshot shows the 'Digi-Sense Data Acquisition Utility' window. The interface is divided into several sections:

- Device Section:** Contains 'Current Temp:' and 'Set Temp:' input fields, both with a red box around them and a label 'Sensor Temperature' pointing to the 'Current Temp' field. Next to each field are 'Up' and 'Down' buttons. Below these are radio buttons for 'F' and 'C', with 'C' selected.
- Connection Section:** Shows 'Status: Not connected' with a red box and label 'Status of USB connection to control'. Below it is a 'COM Port Settings...' button with label 'Press button to change COM Port'. At the bottom are 'Connect' and 'Disconnect' buttons with label 'Use these buttons to Connect or Disconnect USB communications to the computer'.
- Logging Section:** Includes 'Current logging file:' with a text box and a browse button ('...') with label 'Enter filename and location for data log file'. Below is 'Logging Interval:' set to '1.00' with label 'Rate the data is saved. Defaults to 1 second intervals.'. A checked 'Enable graph' checkbox has label 'Live graph function'. At the bottom are 'Start', 'Stop', 'Read File', and 'Clear Graph' buttons with label 'Data Logging control buttons'.
- Graph Area:** A large plot area on the right with a y-axis from 0.0 to 1.2 and an x-axis with '01' markers. It is labeled 'Live Graph Area'.

The status bar at the bottom left of the window displays 'Not connected'.

Data Acquisition Software Setup

Setting Temperature Set Point

- Use the UP and DOWN buttons beside Set Temp to adjust temperature set point.
- The user can also manually type in the temperature by placing the mouse pointer in Set Temp box. Then type in set point value and press the Set button.
- This will update the temperature on the Cole-Parmer controller.


Device _____
Current Temp: _____ °C Up
Set Temp: _____ °C Set Down
 F C

Data Logging Setup

Logging _____
Current logging file: _____ ...

1. Enter file name into Current logging file. Press the button to change file directory path.

2. Data logging interval defaults at 1.00 seconds. Change the value by left-clicking with pointer located inside the box.

3.  Logging Interval: _____ Type in new interval time and left-click with mouse in the graph area.

4. Enable graph Enable Graph will have a check in the box as the default setting. If the live graph is not wanted, left-click with mouse pointer over the box to turn off the feature.

5. Start/stop data logging and graphing function:

Start button — Select this button to start graphing and data logging.

Stop button — Stops the data logging and graphing when the Start button is active.

Start	Stop
Read File	Clear Graph

Read File button — Loads a previous data file for viewing on graph.

Clear Graph button — Clears all data on the graph.

Safety Precautions



DANGER: DO NOT REMOVE COVER! HIGH VOLTAGE IS PRESENT IN THE CONTROL. Contact vendor for service.



DANGER: If high voltage is present on external temperature sensor from outside source, high voltage will be present at the control.



DANGER: Fire protection and control damage: Replace all fuses with the correct fuse replacement. Reference page 35 for model 89800-13 and page 36 for model 89800-14.



WARNING: Specifications for the power cord: see page 35 for model 89800-13 and page 36 for model 89800-14 for proper replacement cord. Additional input power cords for various countries are listed on page 41.



WARNING: Use of separate temperature limit control is recommended were a fault condition could occur and result in a fire or other hazardous condition.

Specifications for TC9600 Controller — Model 89800-13

Power input: 120 VAC $\pm 10\%$, 50/60 Hz $\pm 3\%$, 15 amp, 1800 watts max load

Operating environment: 32 to 77°F (0 to 25°C); 90% RH, noncondensing

Maximum altitude: 2187 yd (2000 m)

Pollution degree: 2 (normally only nonconductivity pollution occurs)

Installation category II: local level (connect to branch circuit and not directly to a main circuit, such as a fuse panel)

Storage: 32 to 140°F (0 to 60°C); 5 to 80% RH, noncondensing

Fuse: 250 volt, 15 amp rated (fast-acting)

AC line cord: SJT-14-3 14 AWG, 15 amp, 125 VAC, less than 9 ft (3 m) in length

Heater/cooler output: max voltage: 120 VAC $\pm 10\%$, 15 amp, 50/60 Hz $\pm 3\%$, 1800 watts max load

External alarm output: SPDT relay — 24 VAC/VDC, 1 amp

Process memory: data retention upon power failure via nonvolatile memory

Dimensions (W x H x D): 8" x 3.75" x 9" (20.3 x 9.5 x 22.9 cm)

Specifications for TC9600 Controller — Model 89800-14

Power input: 230 VAC $\pm 10\%$, 50/60 Hz $\pm 3\%$, 10 amp, 2300 watts max load

Operating environment: 32 to 77°F (0 to 25°C); 90% RH, noncondensing

Maximum altitude: 2187 yd (2000 m)

Pollution degree: 2 (normally only nonconductivity pollution occurs)

Installation category II: local level (connect to branch circuit and not directly to a main circuit, such as a fuse panel)

Storage: 32 to 140°F (0 to 60°C); 5 to 80% RH, noncondensing

Fuse: 250 volt, 10 amp rated (fast-acting)

AC line cord: SJT-14-3 14 AWG, 15 amp, 240 VAC, less than 9 ft (3 m) in length

Heater/cooler output: max voltage: 240 VAC $\pm 10\%$, 15 amp, 50/60 Hz $\pm 3\%$, 3600 watts max load

External alarm output: SPDT relay — 24 VAC/VDC, 1 amp

Process memory: data retention upon power failure via nonvolatile memory

Dimensions (W x H x D): 8" x 3.75" x 9" (20.3 x 9.5 x 22.9 cm)

Specifications for Sensor Input

- Thermocouple (grounded or nongrounded), RTD (Type Alpha 0.003850 and Alpha 0.003916), and Thermistor (Type 400 and 700)
- Automatic cold junction compensation and break protection for sensor

Ranges

Thermocouples	
Type J	-310 to 1832°F (-190 to 1000°C)
Type K	-328 to 2502°F (-200 to 1372°C)
Type N	-328 to 2372°F (-200 to 1300°C)
Type R	32 to 3214°F (0 to 1768°C)
Type S	32 to 3214°F (0 to 1768°C)
Type T	-328 to 752°F (-200 to 400°C)
Type B	392 to 3272°F (200 to 1800°C)
Type E	-328 to 1832°F (-200 to 1000°C)
RTDs	
Type 0.003850 100Ω	-328 to 1472°F (-200 to 800°C)
Type 0.003850 100Ω w/ MgO	-328 to 1472°F (-200 to 800°C)
Type 0.003916 100Ω	-328 to 680°F (-200 to 630°C)
Thermistors	
Type 400	-22 to 212°F (-30 to 100°C)
Type 700	5 to 212°F (-15 to 100°C)

Sensor accuracy for thermocouples and RTDs


Calibration accuracy
 Thermocouple types J, K, T, E, N, $\pm 0.1\%$ of span or $\pm 1^\circ\text{C}$
 Thermocouple types B, R, S, $\pm 0.2\%$ of span or $\pm 2^\circ\text{C}$
 RTD; $\pm 0.1\%$ of span
 Accuracy span is 1000°F (540°C) minimum

Sensor accuracy for thermistors


Calibration accuracy
 Thermistors: $\pm 0.4\%$ of span or $\pm 0.5^\circ\text{C}$
 Accuracy span is 266°F (130°C) minimum

Screen Flow Charts

-ADVANCED MODEL-

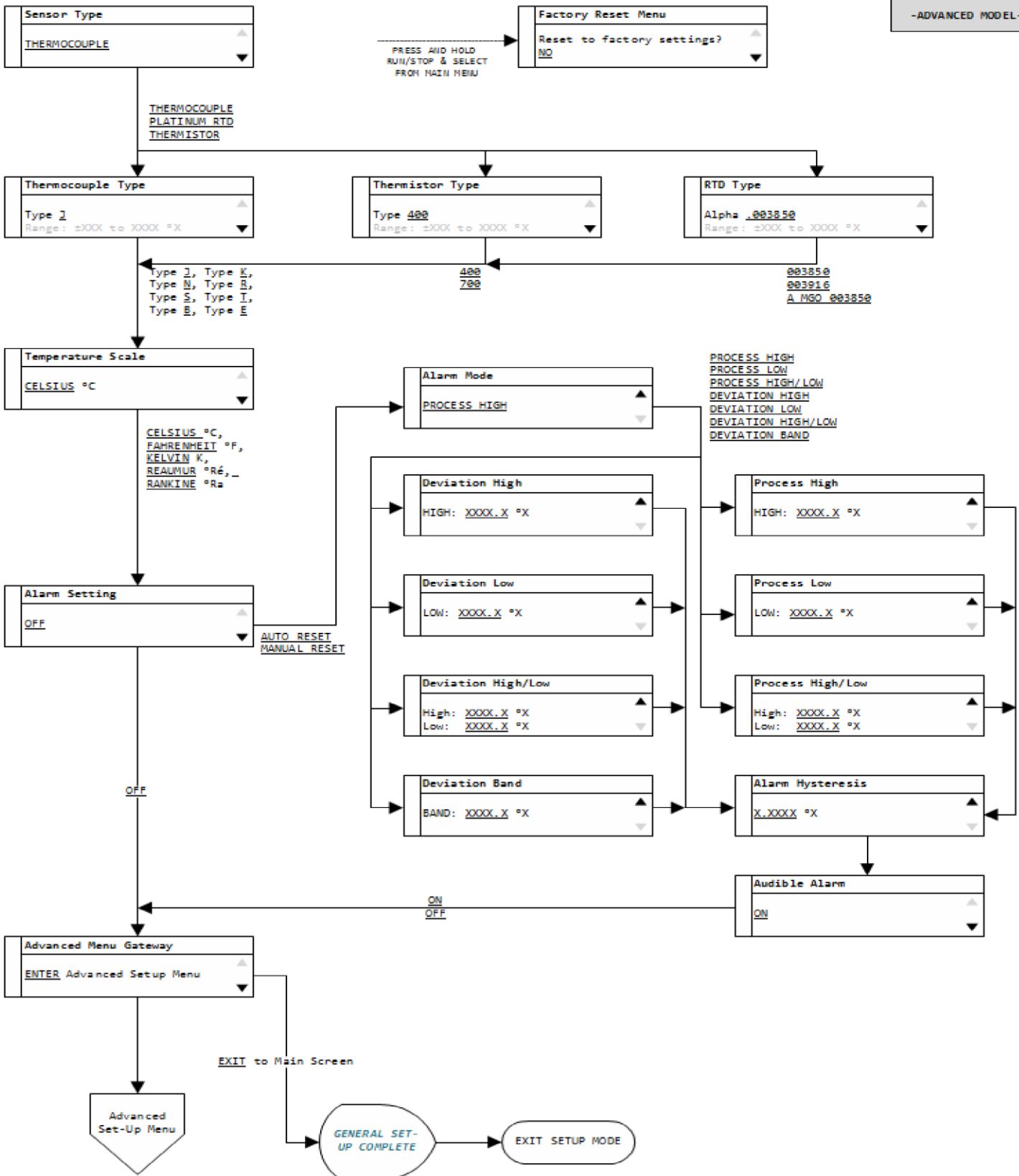

Screen Flow Charts

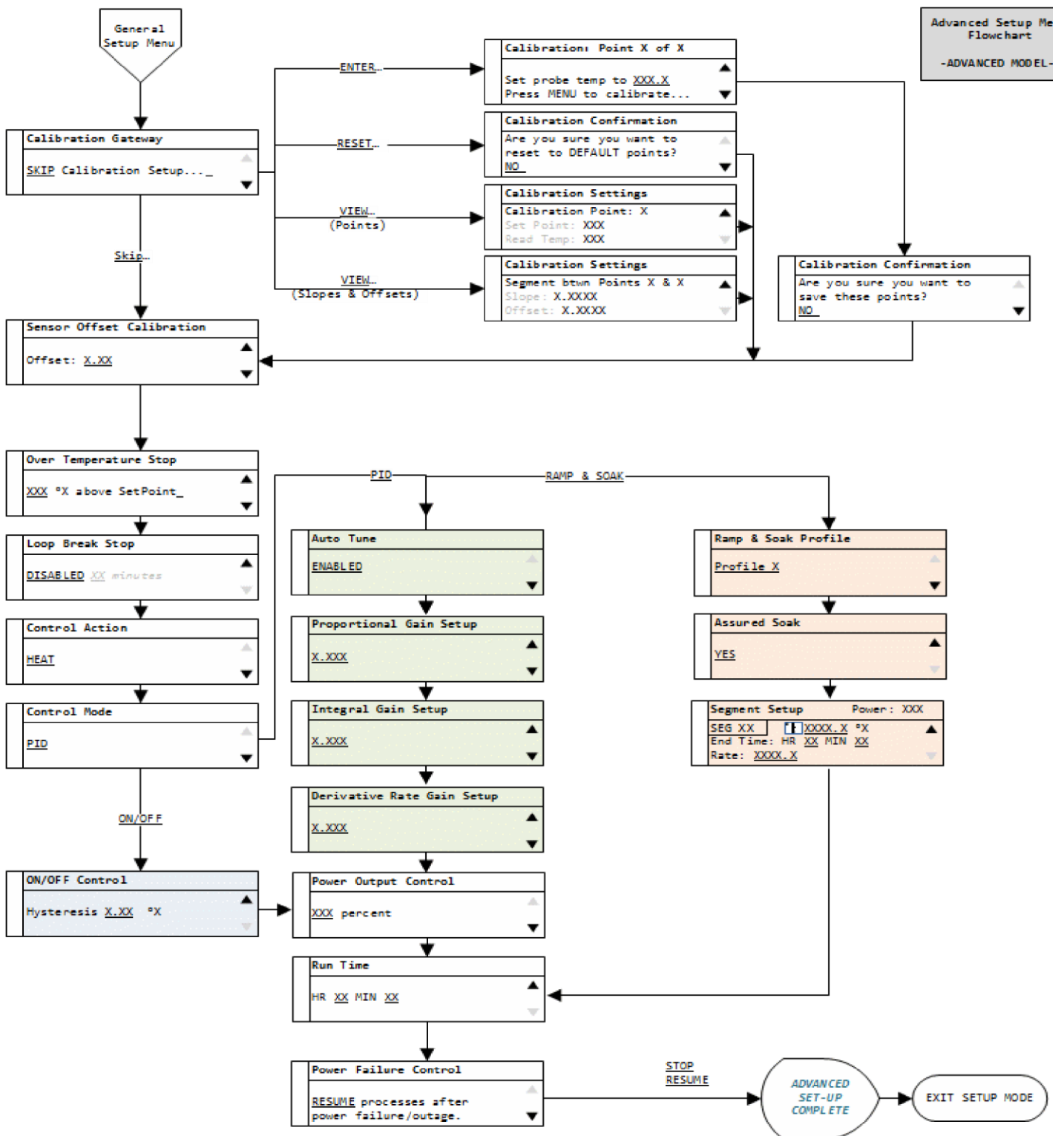

Correct Wiring of ANSI (3-Blade) RTD Probe

- If the controller shows an “open” sensor alarm when using an RTD sensor with an ANSI Style 3-blade connector:
 - Please verify that the probe is wired correctly to industry standards.
 - Open the connector of the probe and confirm the wiring is per the picture below.
 - If the wirings does not match, please rewire as shown.



Ferrite Clip Installation




Alternative Power Cords — for Various Countries

A detachable cord/plug set is automatically included with both models of the temperature controller:

- Model 89800-13 includes a US 120 VAC plug
- Model 89800-14 includes a US 220 VAC plug

Below is a ordering table for the available cord/plug sets. Cord/plug sets feature a country-specific male plug on one end and an IEC 320 female plug on the other end. Order a cord/plug set to replace a lost or damaged set or to use your temperature controller in another country.

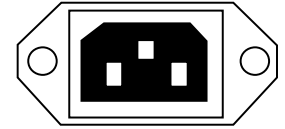

IEC 320 socket

Illustration	Country	Catalog number
	US Standard	50001-68
	Australia, Japan	50001-60
	Denmark	50001-62
	India	50001-64
	Israel	50001-69
	Europe	50001-70
	England	50001-72
	Switzerland	50001-74
	Italy	50001-76
	US (NEMA)	50001-78

Output Cord Adapters – for Various Countries

A detachable cord set is not included with the temperature controller:

- Model 89800-13 includes a US 120 VAC female plug
- Model 89800-14 includes a IEC 60320 C19, 230 VAC female plug

Below is an ordering table for the available cord sets. Cord sets feature a country specific female plug on one end and an IEC 360320 C19 male plug on the other end that will plug into the controller. Order a cord set to use your heating or cooling devices with your 230 VAC temperature controller in IEC 60320 C19 another country. Each cord set is 12 in (30.5 cm) in length.



Illustration	Country	Catalog number
	Australia, Japan	80800-24
	Denmark	80800-21
	India	80800-23
	Israel	80800-28
	Europe	89800-19
	England	89800-22
	Switzerland	80800-26
	Italy	80800-27
	US (NEMA)	80800-29

CE Approval

Conforms to the following Product Specifications:

EMC: EC Directive 204/108/EC
Using the following: EN 61326-1-2006
EN 55011: 2009 ISM RADIO FREQUENCY EQUIPMENT, CLASS B
EN 61000-3-2:2006 (IEC 61000-3-2: 2005) HARMONICS
En 61000-3-3:2008 (IEC 6100-3-3:2008) FLICKER
Immunity Test:
EN 61000-4-2: 2001-4 ESD
EN 61000-4-3: 2006-02 RADIATED RF SUSCEPTIBILITY
EN 61000-4-4: 2004-07 ELECTRICAL FAST TRANSIENT / BURST
EN 61000-4-5: 2001-04 SURGE
EN 61000-4-6: 2004-11 RF CONDUCTED SUSCEPTIBILITY, COMMON MODE
EN 61000-4-11: 2004-03 VOLTAGE DIPS AND INTERRUPTIONS

Safety: 2006/95/EC using EN61010-1

UL Approval—File E207546, Vol.

Standard: UL 61010-1, 3rd Edition, 2012-04-17 (ELECTRICAL EQUIPMENT FOR MEASUREMENT, CONTROL, AND LABORATORY USE - Part 1: General Requirements)
CAN/CSA-C22.2 No. 61010-1, 3rd Edition, 2012-04, (ELECTRICAL EQUIPMENT FOR MEASUREMENT, CONTROL, AND LABORATORY USE - Part 1: General Requirements)

Certification Type: Listing

CCN: OGTK, OGTK7 (Laboratory Use Electrical Equipment)

Maintenance

- Simple preventive maintenance steps include keeping the controller clean. Protect it from overload, excessive dirt, oil and corrosion.
- Cleaning: If cleaning is necessary, use only a damp cloth with water only. Wipe only the exterior of the control chassis.

CATALOG NUMBERS **89800-13 and 89800-14**

SERIAL NUMBER _____

DATE OF PURCHASE _____

For Product and Ordering Information, Contact:

Cole-Parmer[®]
essentials

Toll-Free: 1-800-323-4340

Phone: 1-847-549-7600

Fax: 1-847-247-2929

ColeParmer.com